

Jul 21, 2016

eDNA Fish Metabarcoding PCR; 12S & cytB

DOI

dx.doi.org/10.17504/protocols.io.fckbiuw

Harriet Johnson

EvoHull



Harriet Johnson

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.fckbiuw

External link: https://github.com/SchlossLab/MiSeq_WetLab_SOP/blob/master/MiSeq_WetLab_SOP_v4.md

Protocol Citation: Harriet Johnson 2016. eDNA Fish Metabarcoding PCR; 12S & cytB. [protocols.io](https://dx.doi.org/10.17504/protocols.io.fckbiuw)
<https://dx.doi.org/10.17504/protocols.io.fckbiuw>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](https://creativecommons.org/licenses/by/4.0/), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

Created: July 20, 2016

Last Modified: March 21, 2018

Protocol Integer ID: 3180

Abstract

One-step metabarcoding PCR. Used to amplify either mitochondrial 12S or cytochrome B regions from environmental DNA sampled from freshwater to identify fish species. PCR product contains the sequencing primers ready for sequencing on the Illumina MiSeq (custom oligos required for sequencing).

Guidelines

References to primer sequences and their use with this method:

Kocher, Thomas D., et al. "Dynamics of mitochondrial DNA evolution in animals: amplification and sequencing with conserved primers." *Proceedings of the National Academy of Sciences* 86.16 (1989): 6196-6200.

Riaz, Tiayyba, et al. "ecoPrimers: inference of new DNA barcode markers from whole genome sequence analysis." *Nucleic Acids Research* (2011): gkr732.

Kelly, Ryan P., et al. "Using environmental DNA to census marine fishes in a large mesocosm." *PloS one* 9.1 (2014): e86175.

Hänfling, Bernd, et al. "Environmental DNA metabarcoding of lake fish communities reflects long-term data from established survey methods." *Molecular ecology* (2016).

Safety warnings

- ⚠ Custom oligos are required for Illumina sequencing of this PCR product. In order to also sequence PhiX on the run, spike in the custom oligos to the supplied standard Illumina sequencing primers.

The sequencing library preparation protocol will also be added here. I have put a link to the Schloss library prep protocol for now (see links).

1 **Dual indexed primers:**

16 forward primers, 24 reverse primers: FA, FB, RA, RB. Can mix and match to create 384 unique combinations to identify multiplexed samples. See 'guidelines' for references to primer sequences.

Dilute primers to 10 uM concentration with water.

Plate primers in matrix on 96 well plate or strip tubes to create unique combinations:
eg.

RA701 - —→ RA712
FA501
|
FA508

Note

Multiple dilutions are recommended for making working concentration primers to minimise handling of stock primers and to avoid contamination.

Expected amplicon size:

Cytochrome B: 460 bp region (F primer: 74, R primer: 70), total: 604 bp
12S: 106 bp region (F primer: 67, R primer: 62), total: 235 bp

Note

Positive control: *Rhamphochromis esox*, a pelagic species in its natural habitat and isn't common in UK aquariums (according to Alan Smith). Info from Lynsey Harper.

SEPA Jan/Feb 2016: 2 µl 1/100 dilution DNA.

EA, April/May 2016 LFC project: 0.5 ul 1/500 dilution DNA.

Negative control: Molecular grade water

Note

No. of cycles were reduced from 40x to 35x, told of 'updated' protocol when I started, do not know when this was changed.

All rxns I have performed have been 35x.

2 **PCR reaction setup**

20 ul rxn: Windermere samples:

10 µl Q5 2x mastermix



1 μ l F primer
1 μ l R primer
2 μ l DNA
6 μ l H₂O

20 μ l rxn: Scottish Loch samples:

10 μ l Q5 2x mastermix
1 μ l F primer
1 μ l R primer
4 μ l DNA
4 μ l H₂O

35 μ l rxn: LFC samples:

17.5 μ l Q5 2x mastermix
1.75 μ l F primer
1.75 μ l R primer
3.5 μ l DNA
10.5 μ l H₂O

PCR strip tubes with individual lids, instead of open plates.

Note

Multiple dilutions are recommended for making working concentration primers to minimise handling of stock primers and to avoid contamination.

Expected amplicon size:

Cytochrome B: 460 bp region (F primer: 74, R primer: 70), total: 604 bp
12S: 106 bp region (F primer: 67, R primer: 62), total: 235 bp

Note

Positive control: *Rhamphochromis esox*, a pelagic species in its natural habitat and isn't common in UK aquariums (according to Alan Smith). Info from Lynsey Harper.

SEPA Jan/Feb 2016: 2 μ l 1/100 dilution DNA.

EA, April/May 2016 LFC project: 0.5 μ l 1/500 dilution DNA.

Negative control: Molecular grade water

Note

No. of cycles were reduced from 40x to 35x, told of 'updated' protocol when I started, do not know when this was changed.

All rxns I have performed have been 35x.

3 Thermal cycling parameters:

Initial denaturation: 98°C 5 min;

35X of

1. 98°C 10 sec
2. 58 °C 15 sec
3. 72°C 20 sec;

Final extension 72°C 5 min

Store at 10°C

Note

Multiple dilutions are recommended for making working concentration primers to minimise handling of stock primers and to avoid contamination.

Expected amplicon size:

Cytochrome B: 460 bp region (F primer: 74, R primer: 70), total: 604 bp

12S: 106 bp region (F primer: 67, R primer: 62), total: 235 bp

Note

Positive control: *Rhamphochromis esox*, a pelagic species in its natural habitat and isn't common in UK aquariums (according to Alan Smith). Info from Lynsey Harper.

SEPA Jan/Feb 2016: 2 µl 1/100 dilution DNA.

EA, April/May 2016 LFC project: 0.5 ul 1/500 dilution DNA.

Negative control: Molecular grade water

Note

No. of cycles were reduced from 40x to 35x, told of 'updated' protocol when I started, do not know when this was changed.

All rxns I have performed have been 35x.

4 Visualise PCR product via gel electrophoresis:

large gel:

1.5g agarose
100 ml 0.5x TBE
10 ul GelRed

Run at 90V 40 mins.

2-3 ul product
1.5ul Easy Ladder

5 Triplicate repeat

SEPA Jan/Feb 2016:
20 ul rxns, performed in triplicate repeat and then pooled.