

Dec 06, 2024

Version 1

EdU detection on tissue sections V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.e6nvwbdrzvmk/v1>

Ashley Seifert¹, Sarah Calve¹

¹University of Kentucky



Ashley Seifert

University of Kentucky

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.e6nvwbdrzvmk/v1>

Protocol Citation: Ashley Seifert, Sarah Calve 2024. EdU detection on tissue sections. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.e6nvwbdrzvmk/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: December 06, 2024



Last Modified: December 06, 2024

Protocol Integer ID: 114538

Keywords: ASAPCRN, seifert lab protocol for fluorescent immunohistochemistry, detection on tissue section, seifert lab protocol, seifert lab, sensitive detection of dna synthesis, procedure for edu detection, fluorescent immunohistochemistry, immunohistochemistry, edu detection, sensitive detection, tissue section, embedded tissue section, detection, dna, vivo, lab, dna synthesis, chemical method

Abstract

This protocol describes the procedure for EdU detection used by the Seifert Lab. It assumes a starting sample of a paraffin-embedded tissue section and refers to the Seifert Lab protocol for fluorescent immunohistochemistry [add link when published].

It is adapted from the procedure described in A. Salic, T.J. Mitchison, A chemical method for fast and sensitive detection of DNA synthesis *in vivo*, *Proc. Natl. Acad. Sci. U.S.A.* 105 (7) 2415-2420, <https://doi.org/10.1073/pnas.0712168105> (2008).

Protocol materials

⊗ Tris-buffered saline (TBS), 1x solution **Fisher Scientific Catalog #BP24721**

⊗ Hoechst 33342 **Catalog #H3570**

⊗ Prolong Gold **Thermo Fisher Scientific Catalog #P36930**



- 1 Deparaffinize and rehydrate slides per normal IHC protocol
- 2 If using subsequent antibodies for IHC, perform antigen retrieval



- 3 Rinse slides two times in

Tris-buffered saline (TBS), 1x solution **Fisher Scientific Catalog #BP24721**

- 4 *OPTIONAL*

Block for 00:05:00 to 00:10:00 in appropriate blocking solution



Note

This is not necessary for rodent ear or skin tissue

- 5 Leave slides in TBS until reaction cocktail is ready
- 6 Prepare EdU reaction cocktail:

	A	B	C
	Component	Volume (μL)	Notes
	2 M Tris, pH 8.5	5	100 mL = 24.2 g
	50 mM CuSO ₄	2	1 mL = 7.98 mg
	0.5 M ascorbic acid	20	1 mL = 88 mg *UNSTABLE AT RT, MAKE FRESH
	AlexaFluor Azide (0.25 mg/mL)	0.2	0.5 mg in 2 mL DMSO
	ddH ₂ O	73	



	A	B	C
	Total Volume:	100	

Note

- **IMPORTANT:** use the reaction cocktail within  00:15:00 of preparing
- Use half the indicated AlexaFluor Azide to reduce background staining (0.1 μL per 100 μL)
- 10 mM EdU = 2.55 mg/mL
- too much ascorbic acid messes up phalloidin staining
- 250 nM AlexaFluor azide (AF488) for staining of live cells is 0.45 $\mu\text{L}/\text{mL}$
- Use half the indicated AlexaFluor Azide to reduce background staining (0.1 μL per 100 μL)

7 Remove TBS and blot slides

8 Add reaction cocktail to each slide

Note

 100 μL to  250 μL usually works well

9 Incubate for  00:30:00 at  Room temperature protected from light

10 Wash slides two times for  00:05:00 each in TBS

11 **EITHER:**

Continue with normal fluorescent IHC protocol from blocking step





Note

IMPORTANT: keep slides protected from light for all steps to reduce bleaching of EdU signal

12 **OR:**

Counterstain nuclei by incubating with [M] 1 µg/mL

 Hoechst 33342 **Catalog #H3570** for  00:05:00

12.1 Rinse 3 times with ddH₂O

12.2 Mount with  Prolong Gold **Thermo Fisher Scientific Catalog #P36930**

*