

Jun 05, 2025

Version 1

Efficient Preparation of Competent Cells and Recombinant Protein Expression in E. coli BL21 (DE3) Using IPTG Induction V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.kqdg3wzzqv25/v1>

Durvesh Smita. Burhade¹

¹Parul University

Durvesh Smita. Burhade: SCIENCE IS ELEGANT

Durvesh Burhade Lab



Durvesh Burhade

Parul University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kqdg3wzzqv25/v1>

Protocol Citation: Durvesh Smita. Burhade 2025. Efficient Preparation of Competent Cells and Recombinant Protein Expression in E. coli BL21 (DE3) Using IPTG Induction. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.kqdg3wzzqv25/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 03, 2025

Last Modified: June 05, 2025

Protocol Integer ID: 219368

Keywords: BL21 (DE3), Competent Cells, Transformation, Induction & Protein Expression Screening, recombinant protein expression, inducing protein expression, protein expression, efficient preparation of competent cell, using iptg induction reproducible method, iptg induction reproducible method, using iptg, preparing tss, plasmid, competent cell, bl21

Abstract

Reproducible method for preparing TSS-competent *E. coli* BL21 (DE3) cells, transforming them with plasmids, inducing protein expression using IPTG and expression analysis by SDS-PAGE.





Competent Cells Preparation by TSS Method (DAY 1)

- 1 Selection of a healthy strain of BL21 (DE3) *E. coli* Culture from an Agar Plate
- 2 Inoculate the isolated colony and incubate
 200 rpm, 37°C, 24:00:00 Overnight


Competent Cells Preparation by TSS Method (DAY 2)

40m

- 3 Prepare 50 mL of L.B Agar and autoclaved it at 121 °C for 15 Minutes
 2 g of L.B Agar in 50 mL Water
- 4 Prepare 100 mL of L.B Broth
 2.5 g in 100 mL of Water
- 5 Dilute the overnight culture in a 1:100 ratio
 1 mL of overnight culture in 99 mL Autoclaved L.B Broth
- 6 Incubate at 180 rpm, 37°C Till it reaches 0.5 - 0.6 OD
- 7 Prepare the TSS Solution (10 mL Total Volume)
10 % of PEG for a 10 mL solution is 1 g
5 % of DMSO for a 10 mL solution is 500 µL
50 mM MgCl₂ for a 10 mL solution is 47 mg
Add all the components in 8.45 g of Autoclaved L.B Broth to make total volume of 10 mL
- 8 After reaching 0.5 OD, cells should be arrested by incubating the culture on ice for 00:30:00 (30 Minutes) 30m
- 9 Cells were harvested by centrifuging at 4000 rpm, 4°C, 00:10:00 , (10 Minutes) 10m




- 10 Discard the supernatant. Suspend the cells by adding  1 mL of TSS Solution by tapping them to ice
- 11 After resuspension of cells, Aliquote  50 µL in each 1.5 mL Eppendorf tubes (Pre-Chilled)

Checking & Validation (DAY 2)









- 12 To check the experiment, add  50 µL of competent cells in 25 ml Agar Plates + Kanamycin Antibiotics
In this step, if we can't see any colonies, it is likely expected that competent cells are ready.

Transformation (DAY 3)

1d 1h 30m







- 13 Make 50 mL L.B Agar and pour it into the plates (25 mL each)
In the upcoming steps, this L.B Media will be used
- 14 Make 25 mL L.B Broth and store in 50 mL Falcon tube
In the upcoming steps, this L.B Media will be used
- 15 Thawed the cells by keeping the eppendorf containing competent cells at  0 °C on ice
- 16 Add  100 ng Plasmid to the Eppendorf
100ng= 0.1 ug
- 17 Incubate the Eppendorf for 15 Minutes on Ice
- 18 Heat-shock the Eppendorf containing competent cells  42 °C for 45 Seconds
- 19 Put Eppendorf on ice for 2-4 Minutes





- 20 Add  800 μL of L.B Broth in Eppendorf and incubate at  180 rpm, 37°C, 01:00:00 1h
This step is known as regeneration
- 21 After regeneration, harvest the cell by centrifugation  4000 rpm, 37°C, 00:30:00 30m
- 22 Discard the supernatant and add  200 μL of L.B Broth , resuspend it till you see a clear solution 
(Don't resuspend the cells vigorously, it may damage transformant cells.)
- 23 Pour  100 μL of Transformant cells into an agar plate (25 mL plate) and add  10 μL Kanamycin
- 24 Incubate it overnight at  180 rpm, 37°C, 24:00:00 1d
- 25 Colonies are likely to appear if they have an uptake of plasmid


Induction & Expression Screening (DAY 4-5)

1d 20h 15m

- 26 Make 40 mL L.B Broth and aliquot 10 mL each in 15 mL Falcon
- 27 Pick a healthy, medium-sized, isolated colony from an agar plate and inoculate it in  10 mL of L.B Broth (Primary Culture) ; Incubate at  180 rpm, 37°C, 24:00:00 1d
- 28 Take 1%  0.1 μL of the Transformant cells from an overnight culture and inoculate them into 10 mL Falcon (All 4 of them). Incubate them at  180 rpm, 37°C Till it reaches 0.5 OD (Secondary culture)
- 29 Add 0.5 mM of IPTG into two Falcon tubes and 0.1 mM of IPTG into two Falcon tubes.
- 30 Incubate two Falcon tubes containing 0.1 & 0.5 mM concentration of IPTG at  25 °C 20h
 16:00:00

Incubate two Falcon tubes containing 0.1 & 0.5 mM concentration of IPTG at  37 °C

 04:00:00

31 Harvest all 4 Falcon Tubes at  4000 rpm, 37°C, 00:15:00
(At their respective timing)

15m

32 Run SDS-PAGE and check the protein expression