


Jun 14, 2019

Electroporation of *Caecitellus sp.* with FITC-dextran

DOI

dx.doi.org/10.17504/protocols.io.35kgq4w

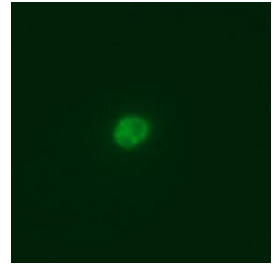

Nick Irwin¹, Elisabeth Hehenberger¹, Patrick Keeling¹

¹University of British Columbia

Protist Research to Opti...



Nick Irwin

University of British Columbia

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.35kgq4w

Protocol Citation: Nick Irwin, Elisabeth Hehenberger, Patrick Keeling 2019. Electroporation of *Caecitellus sp.* with FITC-dextran. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.35kgq4w>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: In development

We are still developing and optimizing this protocol

Created: June 13, 2019

Last Modified: June 14, 2019

Protocol Integer ID: 24460



Abstract

Caecitellus sp. is a small phagotroph belonging to the biocosoecid order of the stramenopiles. *Caecitellus* is found across global oceans but its role in marine ecosystems, as with many heterotrophic flagellates, is unclear.

We isolated a strain of *Caecitellus* (small subunit rRNA gene was 99.53% identical to *Caecitellus parvulus* and *Caecitellus paraparvulus*) from a contaminated culture of *Oxyrrhis marina* and assessed its amenability to transformation.

Fluorescein isothiocyanate conjugated to dextran was successfully introduced into *Caecitellus* sp. using a high voltage exponential decay pulse (1000 V, 10 μ F, ∞ Ω) from a BioRad GenePulser Xcell (cuvette width 0.2 cm). Cell viability was confirmed by observing motility and fluorescence 24 h after electroporation (Figure 1). Lower voltage pulses did not result in FITC uptake. Estimates of transformation efficiency were difficult to obtain due to the small size of the cells.

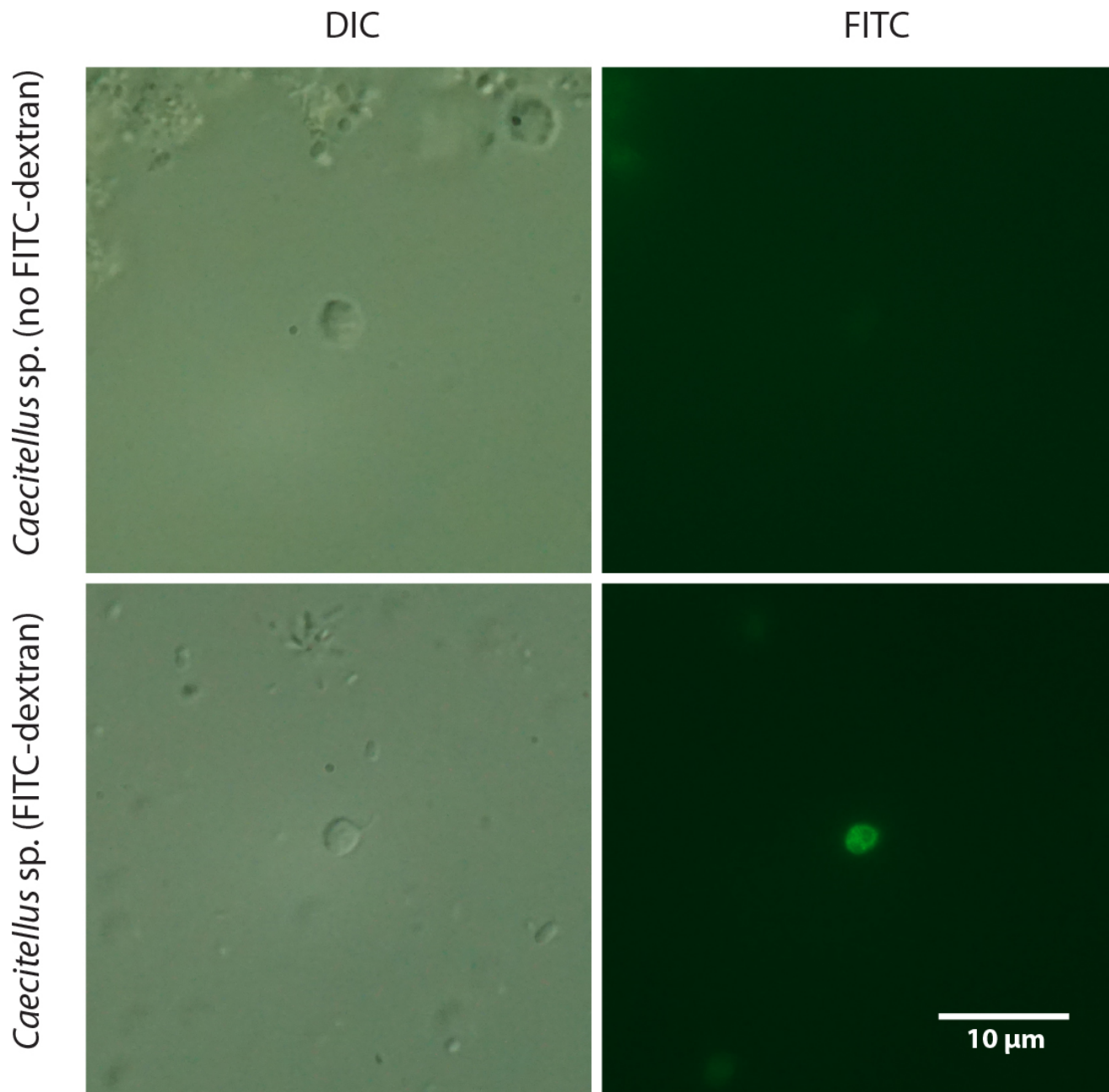


Figure 1. Electroporation of *Caecitellus* sp. with FITC-dextran.

Unfortunately, the isolated strain of *Caecitellus* perished in an incubator malfunction. However, these results may be transferable to *Caecitellus parvulus*, a model for heterotrophic flagellates.



Culturing

- 1 Cells were grown in f/2 media in a 20 C incubator on a 12:12 light:dark cycle.

Cells fed on natural populations of bacteria in the culture and did not need supplemental food.

Electroporation

- 2 Spin down cells at 3000 x g for 10 minutes.

Resuspend cells in 100 uL of BioRad Gene Pulser electroporation buffer with 2 mg/mL FITC-dextran.

Electroporate the cells using a high voltage exponential decay pulse (1000 V, 10 μ F, ∞ Ω) from a BioRad GenePulser Xcell (cuvette width 0.2 cm).

After electroporation, add 900 uL of f/2 and transfer to a 1.5 mL centrifuge tube.

Wash the cells once by spinning at 7000 xg for 3 minutes, removing the supernatant, and resuspending in 1 mL fresh f/2

Place cells in a 24-well plate and allow them to recover at room temperature in the dark for 24 hours

Observations

- 3 Observe the cells 24 hours after electroporation.

Successful transformants will appear strongly fluorescent green and will be swimming.