

Oct 26, 2018

Electroporation protocol for *Vibrio natriegens*

DOI

dx.doi.org/10.17504/protocols.io.u3geyiw

Josef Hoff¹

¹Phillips University



Josef Hoff

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.u3geyiw>

Protocol Citation: Josef Hoff 2018. Electroporation protocol for *Vibrio natriegens*. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.u3geyiw>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 26, 2018



Last Modified: October 26, 2018

Protocol Integer ID: 17224

Keywords: electroporation protocol for vibrio natriegens electroporation protocol, vibrio natriegens electroporation protocol, electroporation protocol, electroporation, sterile filtration, chilled cell, chilled cells from the cold, dna suspension, chilled electroporation cuvette, electrocompetent cell, vibrio natriegen, electroportor, microcentrifuge tube, filtration, v2 salt,



Abstract

Electroporation protocol

Weinstock paper:

Matthew T Weinstock, Eric D Heseck, Christopher M Wilson, Daniel G Gibson

Vibrio natriegens as a fast-growing host for molecular biology

Nature Methods volume 13, pages 849–851 (2016)

To prepare before:

recovery medium

(BHI + v2 salts (204 mM NaCl, 4.2 mM KCl, 23.14mM MgCl₂), and 680 mM sucrose) sterile filtration

* i have a box of aliquots in the freezer and pre heat them before use

-A vial of competent cells is retrieved from storage at –80 °C and allowed to thaw on ice.

-Plasmid DNA and electrocompetent cells are combined and gently mixed in a chilled 1.5-mL microcentrifuge tube.

-The cell–DNA suspension is transferred to a **chilled** electroporation cuvette with a 0.1-cm gap size.

-Cells are electroporated with 900V (in our Electroportor we cant set other parameters)

* in the Weinstock paper they recommend depending on the strain 700–900V, 25 µF and 200 Ω

-Cells are **immediately** recovered in 500 µL **preheated** (50°C) recovery medium and transferred to a 1,5-mL tube.

* we preheat the media to 50°C because the recovery media is cooled down by pipetting and up taking the chilled cells from the cold cuvettes

-The cells are recovered by incubating at 37 °C for 1.5h.

(also put the agar plates for preheating in the incubator at 37°C)

-The cells are centrifuged down for one minute at 3000g. The supernatant is then discarded.

- the pellet is resuspended in the leftover of the media and plated out on

warm agar plates containing appropriate antibiotic.

* for Chloramphenicol 2µg/mL ; for Kanamycine 200µg/mL; for Carbenicillin 200µg/mL

-The plates are incubated for several hours or overnight at 37 °C for colonies to appear.

Attachments



Electroporation prot...

48KB

Troubleshooting

