

Apr 08, 2019 Version 2

Electrotransformation of Clostridium species V.2

DOI

dx.doi.org/10.17504/protocols.io.zvsf66e

Chin Yee Tan¹

¹Duke University



Chin Yee Tan

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.zvsf66e

Protocol Citation: Chin Yee Tan 2019. Electrotransformation of Clostridium species. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.zvsf66e>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: In development

We are still developing and optimizing this protocol

Created: April 08, 2019













Last Modified: April 08, 2019

Protocol Integer ID: 22162

Abstract

A brief protocol for electrotransformation of Clostridium species



- 1 Inoculate  10 μ L BHI-supplemented broth with  100 μ L stock culture overnight.
- 2 Use the overnight culture to inoculate  100 mL BHI-supplemented broth to a starting density of OD 0.02.
- 3 Harvest early-exponential phase culture (OD 0.2 to 0.25) by centrifugation by 12,000g x  00:15:00 at  25 °C .
- 4 Wash once in  10 mL SML electroporation buffer .
- 5 Resuspend in  10 mL SMP electroporation buffer .
- 6 Take  400 μ L of aliquots and and mix with  500 ng of DNA, transfer to prechilled cuvettes with 0.2cm gap.
- 7 Incubate on ice for  00:10:00 .
- 8 Electroporate at the following parameters: 25 μ F, resistance 200 Ω , voltage 1.8kV.
- 9 Immediately transfer the cells into  10 mL of BHI broth and incubate for  03:00:00 .
- 10 Plate cells in dilutions on solid selective and non-selective BHI agar.