

May 30, 2019

 Elisa

DOI

<https://dx.doi.org/10.17504/protocols.io.3iagkae>

Alex Andre¹

¹paris 6



Alex Andre

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.3iagkae>

Protocol Citation: Alex Andre 2019. Elisa. protocols.io <https://dx.doi.org/10.17504/protocols.io.3iagkae>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 30, 2019

Last Modified: May 30, 2019



Protocol Integer ID: 23842

Keywords: elisa titration, antibodies in the serum, antibody, serum, patient

Abstract

Titration of antibodies in the serum of patients.

Guidelines

prepare the plate the day before

Materials

MATERIALS

⊗ PBS

⊗ AggreWell™ 400 24-well plate, 1 pack 1 Plate **STEMCELL Technologies Inc. Catalog #34411**

⊗ anti-G3BP1 and anti-PABP1 primary antibodies **Abcam**

test

Safety warnings

! inactivate the serum

Before start

put the reagents at RT



coating

1

add ab at 1M 1 micromolar (μ M)

2

wash and keep at 4 °C

Blocking

3

add 100ul BSA

4

incubate 01:00:00

20 °C

Adding of sample

5

add the serum in the plate

6

incubate 20 °C

01:00:00

7

Wash

Detection



8 Add Ab at [M] 10 micromolar (μM)

9 incubate 🌡 Room temperature ⌚ 00:00:00