



Aug 04, 2024

🌐 Endocytosis and internalization assay in primary neuronal culture

📖 [Science Advances](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.4r3l22j5ql1y/v1>

Arpine Sokratian¹, andrew.west west²


¹Duke Univeristy; ²Duke University

West lab protocols



Arpine Sokratian

Duke Univeristy

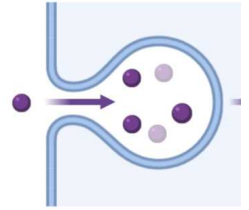

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l22j5ql1y/v1>

Protocol Citation: Arpine Sokratian, andrew.west west 2024. Endocytosis and internalization assay in primary neuronal culture. protocols.io <https://dx.doi.org/10.17504/protocols.io.4r3l22j5ql1y/v1>

**Manuscript citation:****DOI:** [10.1126/sciadv.adi8716](https://doi.org/10.1126/sciadv.adi8716)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working**We use this protocol and it's working****Created:** January 22, 2024**Last Modified:** August 05, 2024**Protocol Integer ID:** 93919

Keywords: ASAPCRN, endocytosis in primary neuronal culture, synuclein fibril structure, internalization assay in primary neuronal culture, endocytosis pathway, synuclein, quantifying endocytosis, primary neuronal culture, endocytosis, experimental protein, fibril, internalization assay, protein

Funders Acknowledgements:**ASAP**

Grant ID: ASAP-020527

Abstract

This protocol outlines a method for quantifying endocytosis in primary neuronal cultures, specifically focusing on different α -synuclein fibril structures. By utilizing pH-sensitive dye conjugated to experimental proteins/fibrils, we can identify endocytosis pathways and analyze intensity changes in the pHrodo-specific channel from real-time acquired images at various time points.

Materials

 EIPA **caymanchem Catalog #14406** Dyngo **Abcam Catalog #ab120689** Methyl- β -cyclodextrin **Merck MilliporeSigma (Sigma-Aldrich) Catalog #C4555** Pitstop® 2 **Abcam Catalog #ab120687** Wortmannin **Merck MilliporeSigma (Sigma-Aldrich) Catalog #W1628-1MG**



Protocol materials

⊗ EIPA caymanchem Catalog #14406

⊗ Dyngo Abcam Catalog #ab120689

⊗ Methyl- β -cyclodextrin Merck MilliporeSigma (Sigma-Aldrich) Catalog #C4555

⊗ Pitstop® 2 Abcam Catalog #ab120687

⊗ Wortmannin Merck MilliporeSigma (Sigma-Aldrich) Catalog #W1628-1MG

⊗ Hoechst 33342 Catalog #H3570

⊗ LysoTracker®; Green DND-26 - Special Packaging Thermo Fisher Catalog #L7526

⊗ pHrodo®; Red Dextran, 10,000 MW, for Endocytosis Thermo Fisher Catalog #P10361



Preparation of endocytosis inhibitors

- 1 For the assay we use DIV7 primary hippocampal neuron culture plated in 48-well plates
- 2 Prepare stock solution of the endocytosis inhibitors in DMSO:

A	B
2.3 mM	Wortmannin
50 mM	Dyngo
10 mM	Pitstop 2
384 mM	methyl- β -cyclodextrin (M β CD)
50 mM	ethyl-isopropyl amiloride

Recommended concentrations for stock solutions (stocks can be freezed at -80C for couple of days)

- 3 At the day of the experiment, thaw down the stock solutions in a water bath and dilute the drugs to reach concentrations:

A	B	C	D	E	F	G
Ethyl-isopropyl amiloride (EIPA)	50 uM	500uM: 10 ul of stock + 990ul PBS				
Dyngo	10 uM	10 mM: 20ul of the drug + 80 ul (50% DMSO)	for 100 uM of Dyngo: 5 ul of the 10mM dyngo + 495 ul of PBS			
Wortman nin	0.2 uM	10 uM: 4ul of the stock +	1uM: 200 ul of 10 uM + 800 ul of PBS			

	A	B	C	D	E	F	G
			916 ul PBS				
	Pitstop 2	15 uM	150 uM: 15ul of P + 984 ul PBS				
	Methyl- β- cyclodext rin (MβCD)	2 mM	20 mM: 25 ul of stock + 475 ul PBS				

Drugs are prepared to be diluted to final concentration in the cell culture

4 Before adding protein/drug of interest, add endocytosis inhibitors 30 minutes before the treatment. As calculated for 48-well plates with 3 mL of the media, 30 ul of diluted drugs are sufficient. ** Note: Be careful adding the diluted drugs (should be RT). Place the plate back to the cell culture incubator.

5 After 30 minutes, add the protein/drug of interest supplement with

 Hoechst 33342 **Catalog #H3570**

As for pHrodo-conjugated fibrils, please, see the protocol describing the conjugation process (step 5):

As a control for endocytosis add lysotracker

 LysoTracker[®]; Green DND-26 - Special Packaging **Thermo Fisher Catalog #L7526**

and phrodo-10kDA dextran

 pHrodo[®]; Red Dextran, 10,000 MW, for Endocytosis **Thermo Fisher Catalog #P10361**

Protocol

NAME



Preparation of mouse and human α -synuclein fibrils

CREATED BY

Arpine Sokratian

Preview



- 6 After two hours of incubation, start taking images of each well at the high-speed (high sensitive mode)
Continue acquiring images every two to four hours up to 48 hours of incubation.
- 7 Calculate the images using cell count (Hoechst) for normalization and calculate the signal for each time-point as a proportion to the maxima of the signal

Calculate the signal from LysoTracker after 30 minutes of incubation and compare the control (no endocytosis inhibitors) to the experimental reactions.