

Feb 10, 2020

Endometrium dissociation with trypsin

DOI

dx.doi.org/10.17504/protocols.io.72dhqa6

Regina Hoo¹, Roser Vento-Tormo¹

¹Wellcome Sanger Institute

Vento-Tormo

Tech. support email: rv4@sanger.ac.uk



Regina Hoo

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.72dhqa6

Protocol Citation: Regina Hoo, Roser Vento-Tormo 2020. Endometrium dissociation with trypsin . **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.72dhqa6>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 08, 2019

Last Modified: February 10, 2020

Protocol Integer ID: 28453

Abstract

This protocol is for enrichment of epithelial glands on endometrium; following the step from "Endometrium dissociation with collagenase".

Guidelines

Human samples including tissue, blood and bodily fluids have the potential to harbour HG2 and Hazard Group 3 (HG3) organisms, specifically Blood Borne Viruses (BBVs,); and for brain tissue, CNS tissue and CSF, prions. In the UK we can work with such samples at CL2 on the condition that we do not intend to culture any of the organisms that might be contained in the samples and that the samples haven't already been identified by tests or diagnosis as containing HG3 organisms.

Materials

MATERIALS

 10% FBS RPMI

 FBS **Invitrogen - Thermo Fisher**


 RPMI 1640 Medium **Thermo Fisher Scientific Catalog #11875093**

 Trypsin-EDTA (0.25%) phenol red **Thermo Fisher Scientific Catalog #25200072**

 PBS **Invitrogen - Thermo Fisher**

 DNase I **Sigma Catalog #4716728001**

Safety warnings

-  Samples are unscreened human tissues, please adhere to Biological Safety at Containment Level 2 work procedures.



1 Prepare Trypsin-EDTA-DNaseI mix:

Product	Stock	Final volume (10 ml)	Concentration
Trypsin-EDTA (0.25%) phenol red		9.9 ml	
DNaseI	10mg/ml	100 ul	0.1 mg/ml

2 Following step 10 from "Endometrium dissociation with collagenase" protocol, where pieces of tissue are retained on the cell strainer.

Collect pieces retained on the filter by inverting the filter into a new 50 mL tub and adding 45 ml of PBS with a 1 mL pipette.

3 Centrifuge at 450 g for 00:05:00 together with collagenased sample. Discard supernatants.

4 Resuspend the pieces with 10 mL Trypsin/EDTA 0.25% and incubate for 00:20:00 at 37 °C while shaking.

5 Add 20 mL of RPMI with 10% FBS.



6 Filter material slowly and carefully (very dens liquid) through a 100 µM strainer. Discard retained tissue.




7 Centrifuge at 500 rcf for 00:05:00 at 4 °C . Discard supernatant.


Same step as 14 in "Endometrium dissociation with collagenase" protocol

8 {optional} Resuspend sample with 5 mL of RLB mix and incubate for 00:10:00

*RLB preparation: Dilute 10x RLB stock with water

After RBC lysis, add  10 mL of PBS and centrifuge at 500 rcf for  00:05:00

- 9 Resuspend cells in  1 mL of PBS in an Eppendorf.
- 10 Centrifuge at 500 rcf for  00:05:00 at  4 °C .

Discard supernatant.
- 11 Resuspend with  250 µL of PBS (final volume depending on cell number)
- 12 Filter cells through a cell strainer snap cap tube.
- 13 Count live cells using trypan blue and haematocytometer.
- 14 {Optional} Resuspend cells in freezing medium to a concentration of 1×10^7 cells and aliquot into cryogenic storage vials.