

Oct 17, 2019

Energy solution preparation for OnePot PURE cell-free system

DOI

dx.doi.org/10.17504/protocols.io.8ayhsfw



Konstantinos Ragios¹

¹EPFL - EPF Lausanne

iGEM EPFL



Konstantinos Ragios

EPFL - EPF Lausanne

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.8ayhsfw>

External link: <https://pubs.acs.org/doi/10.1021/acssynbio.8b00427>

Protocol Citation: Konstantinos Ragios 2019. Energy solution preparation for OnePot PURE cell-free system. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.8ayhsfw>



License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 16, 2019

Last Modified: October 17, 2019

Protocol Integer ID: 28728

Keywords: energy, PURE, cell-free, iGEM, expression, tRNA, NTPs, energy solution preparation for onepot pure cell, protein expression in onepot pure cell, onepot pure cell, energy solution preparation, pure cell, protein expression, energy solution, free system in this protocol, cell, protein, free system

Abstract

In this protocol we explain the procedure to create the Energy solution used for protein expression in OnePot PURE cell-free system.

Materials

Material/Consumables:

- Amino Acids (Sigma-Aldrich: LAA21-1KT)
- Magnesium acetate (Sigma-Aldrich: M0631)
- Potassium glutamate (Sigma-Aldrich: 49601)
- DTT (SantaCruz Biotech: sc-29089B)
- ATP (ThermoFisher: R0481)
- GTP (ThermoFisher: R0481)
- CTP (ThermoFisher: R0481)
- UTP (ThermoFisher: R0481)
- tRNA (Roche:10109541001)
- Creatine phosphate (Sigma-Aldrich: 27920)
- Folinic acid (Sigma-Aldrich: PHR1541)
- Spermidine (Sigma-Aldrich: S2626)
- HEPES (Sigma-Aldrich: H0887-100ML)
- biDistilled water

Equipment:

- Vortex
- Nanodrop Spectrophotometer

Troubleshooting

Before start

Keep all the components in ice while preparing the Energy solution.

- 1 For a 2.5x Energy Solution add the materials needed to a tube. The final concentration of the components is presented in Table 1.

Note

Before adding each component make sure it is totally melted and then vortex for a few seconds

2

	Compound	Concentration	Units
	Amino Acids	0.75	<i>mM</i>
	Magnesium acetate	29.5	<i>mM</i>
	Potassium glutamate	250	<i>mM</i>
	DTT	2.5	<i>mM</i>
	ATP	5	<i>mM</i>
	GTP	5	<i>mM</i>
	CTP	2.5	<i>mM</i>
	UTP	2.5	<i>mM</i>
	tRNA	130	<i>UA260/mL</i>
	Creatine phosphate	50	<i>mM</i>



	Folin ic acid	0.05	<i>mM</i>
	Sper midi ne	5	<i>mM</i>
	HEP ES	125	<i>mM</i>

Note

Because tRNA degrades over time before adding it to the solution you need to check the A_{260} units.

- 3 To complete the solution add biDistilled water until you reach the desirable volume, and then vortex for a few seconds
- 4 The Energy solution is stored at -80°C .