



Jul 13, 2023

Enriching and isolating phages in liquid culture

DOI

<https://dx.doi.org/10.17504/protocols.io.ewov1oeeylr2/v1>

Adair Borges¹

¹Arcadia Science

Arcadia Science



Arcadia Science

Arcadia Science

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.ewov1oeeylr2/v1>

Protocol Citation: Adair Borges 2023. Enriching and isolating phages in liquid culture. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.ewov1oeeylr2/v1>

Manuscript citation:

Borges AL, Dutton RJ, McDaniel EA, Radkov A, Reiter T, Weiss ECP. (2023). Isolation of a phage with an arabinosylated genome from a cheese microbial community. <https://doi.org/10.57844/arcadia-743p-ty94>



License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 05, 2023

Last Modified: January 29, 2024

Protocol Integer ID: 82921

Keywords: phage from microbial community, phages in liquid culture, isolating phage, using liquid enrichment, liquid enrichment, microbial community, phage, liquid culture

Abstract

This protocol explains how we isolate phage from microbial communities using liquid enrichment. For a protocol on enriching and isolating on solid agar plates, check out our **companion protocol**.

Liquid growth enrichment

- 1 Set up cultures of your target host strains. We use this protocol for cheese bacterial isolates, which grow best in LB at 25 °C with shaking at 200 rpm, and reach high density after two days.
- 2 After cells have grown to high density, prepare your infection media. We used LB supplemented with 1 mM MgSO₄ and 1 mM CaCl₂ (final concentrations). Add your phage extracts to this media, either individually for many parallel enrichments, or pooled together.

Note

We find that in liquid, enrichments are generally dominated by a single lytic phage. If you are screening multiple phage extracts, and expect that each extract might have a different phage targeting your host of interest, we recommend separate parallel enrichments for each host and phage source. If you instead find that phage × host matches are rare in your system, it might make sense to pool all your phage extracts to reduce the total number of enrichments you need to do.

- 3 Aliquot your infection media into 3 mL glass culture tubes, or deep-well 96-well plates. We generally use 0.25–1 mL of media for these enrichments.
- 4 Subculture your hosts into the infection media at a 1:100 dilution. For instance, you would add 10 µL of host culture to 1 mL of infection media.
- 5 Incubate your enrichments at 25 °C with shaking at 200 rpm overnight.
- 6 Harvest infected subcultures: Move infected cultures into 1.5 mL tube + 50 µL CHCl₃. Spin down for 5 min at max speed, then transfer the supernatant to fresh 1.5 mL tube + 50 µL CHCl₃.
- 7 Use uninfected parent cultures to pour phage plates using the **double agar overlay method**. Add 200 µL of culture to 3 mL of molten (but not overly hot) top agar. Pour top agar onto a room-temperature LB plate, swirling to evenly distribute the agar across the plate before it cools.

**Note****Top agar recipe**

100 mL LB

0.25 g agarose (0.25%)

100 μ L 1 M MgSO_4 (1 mM)

Autoclave to sterilize. Reheat to melt in the microwave, then make sure you let it cool to a point where it is warm to the touch but not scalding, or you will kill your bacterial hosts.

- 8 Once the plate is cooled, spot the phage-enriched supernatant onto the phage plate. If you only have one enrichment per host, you can use a larger volume (like 10 μ L) in one large drop that you can gently swirl with the pipette tip to distribute. If you have many enrichments per host, you should pipette smaller, 2–3 μ L spots onto the host in a grid pattern.
- 9 Incubate plates at room temperature. Plaques will show up in 24–72 hours.
- 10 Pick plaques using a pipette tip, and transfer into 300 μ L of SM buffer with 50 μ L of CHCl_3 .
- 11 Purify these phages by diluting the picked plaque down to a concentrate that will give single plaques, then infecting 200 μ L of host culture with 10 μ L of diluted phage, and finally plating out using the double agar overlay method. Repeat 2 $\times$.

Protocol references

Kropinski AM, Mazzocco A, Waddell TE, Lingohr E, Johnson RP. (2009). Enumeration of Bacteriophages by Double Agar Overlay Plaque Assay. https://doi.org/10.1007/978-1-60327-164-6_7