

Nov 13, 2025

Enzolution™ POLQ Helicase ATPase Assay System Technical Manual

DOI

<https://dx.doi.org/10.17504/protocols.io.e6nvw4n17lmk/v1>

info¹

¹BellBrook Labs LLC

BellBrook Labs

Tech. support phone: +1 (608) 443-2400 email: techsupport@bellbrooklabs.com



info info

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.e6nvw4n17lmk/v1>

Protocol Citation: info 2025. Enzolution™ POLQ Helicase ATPase Assay System Technical Manual. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.e6nvw4n17lmk/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 30, 2025

Last Modified: November 13, 2025

Protocol Integer ID: 231162

Keywords: dependent atpase activity of the polq helicase domain, dna polymerase theta, purified human polq helicase domain, dependent dna polymerase, polq inhibitor, dependent atpase activity, atpase activity, human polq helicase domain, dependent helicase, important roles in double strand break repair, polq helicase domain, formation by the atpase activity, double strand break repair, atp, dependent endonuclease activity, multifunctional enzyme with template, assay, multifunctional enzyme, inhibitor dose response measurement, dna

Abstract

The Enzolution™ POLQ Helicase ATPase Assay System is intended for use with the Transcreener® ADP² Assay Kits to measure the DNA-dependent ATPase activity of the POLQ Helicase domain. POLQ (also known as DNA Polymerase Theta), is a multifunctional enzyme with template-dependent DNA polymerase, ATP-dependent helicase, and dNTP-dependent endonuclease activities that play important roles in double strand break repair. ADP formation by the ATPase activity of the POLQ Helicase domain can be easily measured by the Transcreener® ADP² assay, a far-red, competitive fluorescence assay that enables single addition, mix-and-read detection in a continuous or endpoint format. The assay has been optimized and extensively validated for high throughput screening (HTS) and inhibitor dose response measurements using most multimode plate readers.

The Enzolution™ POLQ Helicase ATPase Assay System provides all reagents required to screen and profile POLQ inhibitors, including the purified human POLQ Helicase domain (amino acids 67-894, N-terminal 6xHis) and ssDNA, when used with the Transcreener® ADP² Assay Kits, which are available with FP, FI and TR-FRET readouts. The protocol is configured for 384-well plates; use of different multi-well plate formats will require adjustment of reagents concentrations utilized in the assay.

Introduction

- 1 The Enzolution™ POLQ Helicase ATPase Assay System is intended for use with the Transcreener® ADP² Assay Kits to measure the DNA-dependent ATPase activity of the POLQ Helicase domain. POLQ (also known as DNA Polymerase Theta), is a multifunctional enzyme with template-dependent DNA polymerase, ATP-dependent helicase, and dNTP-dependent endonuclease activities that play important roles in double strand break repair. ADP formation by the ATPase activity of the POLQ Helicase domain can be easily measured by the Transcreener® ADP² assay, a far-red, competitive fluorescence assay that enables single addition, mix-and-read detection in a continuous or endpoint format. The assay has been optimized and extensively validated for high throughput screening (HTS) and inhibitor dose response measurements using most multimode plate readers.

The Enzolution™ POLQ Helicase ATPase Assay System provides all reagents required to screen and profile POLQ inhibitors, including the purified human POLQ Helicase domain (amino acids 67-894, N-terminal 6xHis) and ssDNA, when used with the Transcreener® ADP² Assay Kits, which are available with FP, FI and TR-FRET readouts. The protocol is configured for 384-well plates; use of different multi-well plate formats will require adjustment of reagents concentrations utilized in the assay.

Key Applications:

- Screening for POLQ Helicase inhibitors (or activators)
- Generating dose response curves and IC₅₀ values for POLQ Helicase inhibitors
- Kinetic and mechanistic analyses

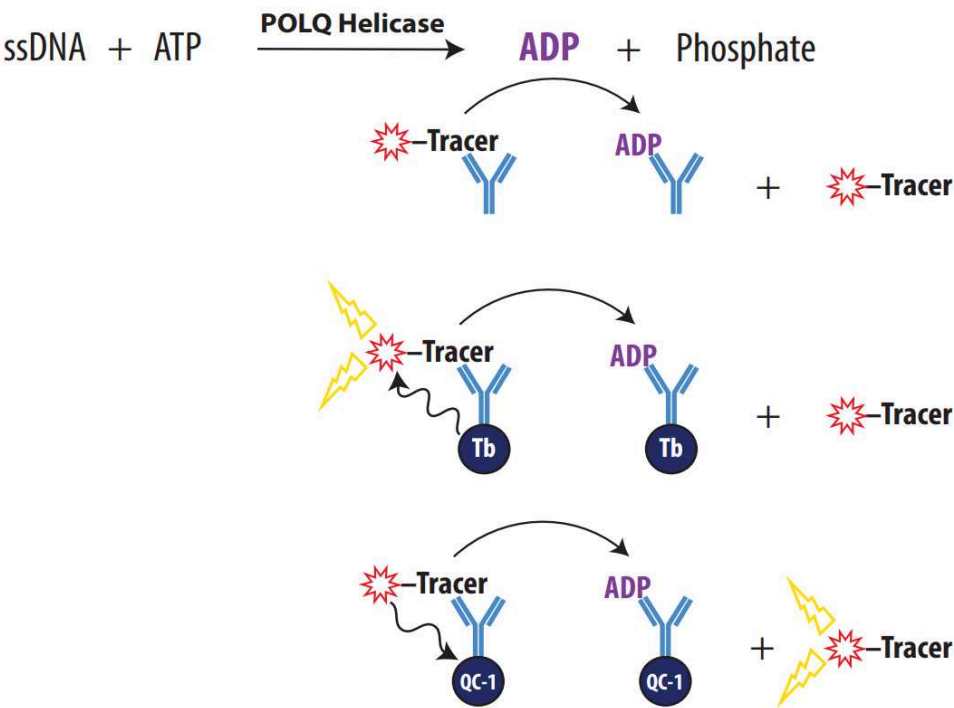


Figure 1. Schematic Overview of the Enzolution™ POLQ Helicase ATPase Assay System with the Transcreener® ADP² Assays.

For FP readout (A): ADP produced by the ATPase activity of POLQ Helicase displaces an Alexa Fluor® 633 Tracer from the ADP² antibody, resulting in decreased fluorescence polarization.

For TR-FRET readout (B): ADP produced by the ATPase activity of POLQ Helicase displaces a HiLyte647 Tracer from the ADP² Antibody conjugated to terbium (Tb), resulting in a decrease in TR-FRET.

For FI readout (C): ADP produced by the ATPase activity of POLQ Helicase displaces an Alexa Fluor® 594 Tracer from the ADP² Antibody conjugated to an IRDye® QC-1 quencher, resulting in an increase in fluorescence intensity

Product Specifications

2

Product	Quantity	Part #	Part #
Enzolution™ POLQ Helicase ATPase Assay System	1,000 assays*	FP	3034-1K-FP
		FI	3034-1K-FI
		TR-FRET	3034-1K-TR

Product	Quantity	Part #	Part #
	10,000 assays*	FP	3034-10K-FP
		FI	3034-10K-FI
		TR-FRET	3034-10K-TR

*The exact number of assays depends on the enzyme reaction conditions. The kits are designed for use with 384-well plates, using a 10 μ L Enzyme Reaction and a 20 μ L Complete Assay volume.

Storage

Enzymes should be stored at -80°C ; other reagents can be stored at -20°C . Though we have confirmed that the POLQ Helicase Enzyme is stable to 6 freeze-thaw cycles, we recommend aliquoting the enzyme and snap-freezing for multiple uses to minimize loss of activity.

Use the reagents provided in this kit within 6 months from date of receipt.

Materials Provided

2.1

Component	Composition	Notes
POLQ Helicase Enzyme	0.1 mg/mL (1.1 μ M)* in 20 mM HEPES, 500 mM NaCl, 5% glycerol, 0.5 mM TCEP, pH 7.5	Amino acids 67-894, N-terminal 6xHis, 93.1 kDa. Sufficient enzyme is included in the kit to complete at least 1,000 assays (Part # 3034-1K) or 10,000 assays (Part # 3034-10K).
POLQ ssDNA	40 μ M in H_2O	Supplied as a 30-bp ssDNA oligomer
Enzyme Assay Buffer A, 10X	500 mM TRIS (pH 7.5), 10 mM MgCl_2 , 0.1% Triton	Use the Enzyme Assay Buffer A in the Enzyme Reaction and for preincubation with inhibitors. Changes to the assay buffer could affect enzyme activity and/or detection of ADP.
384-Well Low Volume Assay Plates	Corning #4514 - FP and FI Only Corning #4513 - TR-FRET Only	Polystyrene non-binding surface assay plates in either a 3-pack (1,000+ Assays) or a 30-pack (10,000+ Assays). We strongly recommend the use of these plates as inconsistent results have been observed with other plates.

*The exact concentration may vary from batch to batch. Please refer to the Certificate of Analysis for an accurate concentration.

Materials Required But Not Provided

2.2

Component	Notes
Ultrapure Nuclease Free Water	Some deionized water systems are contaminated with nucleases that can degrade both nucleotide substrates and products, reducing assay performance. Use nuclease free water such as: Invitrogen Part # AM9930
Plate Reader	A multimode microplate reader configured to measure FP, FI, or TR-FRET is required. Transcreener® Assays have been validated on the following instruments: BioTek Synergy™2 and Synergy™4; BMG Labtech PHERAstar® Plus and CLARIOstar® Plus; Molecular Devices SpectraMax™ Paradigm; Perkin Elmer EnVision® and ViewLux; and Tecan Infinite® F500, Safire2™, and M1000. Full list of compatible plate readers and settings.
Liquid Handling Devices	Use liquid handling devices that can accurately dispense submicroliter volumes into 384-well plates.
Laboratory Incubator	An incubator model that is capable of maintaining temperature stability at 30°C is required.

Transcreener® ADP² FP Assay - SOLD SEPARATELY

Component	Composition	Notes
ADP ² Antibody	3.4 mg/mL solution in PBS with 10% glycerol*	1,000 assays (Part # 3010-1K) or 10,000 assays (Part # 3010-10K).
ADP ² Alexa Fluor® 633 Tracer	400 nM solution in 2 mM HEPES (pH 7.5) containing 0.01% Brij-35	1,000 assays (Part # 3010-1K) or 10,000 assays (Part # 3010-10K).
Stop & Detect Buffer B, 10X	200 mM HEPES (pH 7.5), 400 mM EDTA, and 0.2% Brij-35	The EDTA in the Stop & Detect Buffer B quenches the POLQ Enzyme Reaction by chelating Mg ²⁺ . The final concentrations of Mg ²⁺ and EDTA in the Complete Assay are 1 mM and 20 mM, respectively.
ATP	5 mM	The ATP supplied in this kit can be used for the POLQ Helicase ATPase



Component	Composition	Notes
		Enzyme Reaction and the ATP/ADP standard curve.
ADP	5 mM	The ADP is used for the ATP/ADP standard curve.

Transcreener® ADP² FI Assay - SOLD SEPARATELY

Component	Composition	Notes
ADP ² Antibody-IRDye® QC-1	1.4 mg/mL solution in 100 mM KH ₂ PO ₄ (pH 8.5)*	1,000 assays (Part # 3013-1K) or 10,000 assays (Part # 3013-10K).
ADP ² Alexa Fluor® 594 Tracer	800 nM solution in 2 mM HEPES (pH 7.5) containing 0.01% Brij-35	1,000 assays (Part # 3013-1K) or 10,000 assays (Part # 3013-10K).
Stop & Detect Buffer B, 10X	200 mM HEPES (pH 7.5), 400 mM EDTA, and 0.2% Brij-35	The EDTA in the Stop & Detect Buffer B quenches the POLQ Enzyme Reaction by chelating Mg ²⁺ . The final concentrations of Mg ²⁺ and EDTA in the Complete Assay are 1 mM and 20 mM, respectively.
ATP	5 mM	The ATP supplied in this kit can be used for the POLQ Helicase ATPase Enzyme Reaction and the ATP/ADP standard curve.
ADP	5 mM	The ADP is used for the ATP/ADP standard curve.

Transcreener® ADP² TR-FRET Assay - SOLD SEPARATELY

Component	Composition	Notes
ADP ² Antibody - Terbium Conjugate	800 nM solution in HEPES-buffered saline	1,000 assays (Part # 3011-1K) or 10,000 assays (Part # 3011-10K).
ADP HiLyte647 Tracer	10 µM solution in 2 mM HEPES (pH 7.5) containing 0.01% Brij-35	1,000 assays (Part # 3011-1K) or 10,000 assays (Part # 3011-10K).
Stop & Detect Buffer C, 10X	500 mM HEPES (pH 7.5), 200 mM EDTA, and 0.2% Brij-35	The EDTA in the Stop & Detect Buffer B quenches the POLQ Enzyme Reaction by chelating Mg ²⁺ . The final concentrations of Mg ²⁺ and EDTA in the Complete Assay are 1 mM and 20 mM, respectively.
ATP	5 mM	The ATP supplied in this kit can be used for the POLQ Helicase ATPase

	Component	Composition	Notes
			Enzyme Reaction and the ATP/ADP standard curve.
	ADP	5 mM	The ADP is used for the ATP/ADP standard curve.

*The exact concentration may vary from batch to batch. Please refer to the Certificate of Analysis for an accurate concentration.

Before You Begin

- 3
 1. Read the entire protocol and note any reagents or equipment needed (see **Section 2.2**).
 2. Check the plate reader and verify that it is compatible with the Transcreener® ADP² Assay for the respective readout: Full list of compatible plate readers and settings.
 3. Please read and understand Transcreener® ADP² Assay Technical Manuals prior to use with this kit.

Protocol

- 4
 - The methods described below are for single-addition, endpoint detection: the POLQ Helicase ATPase Enzyme Reaction is quenched by the addition of EDTA along with the detection reagents (see **Figure 2**). The methods were designed for 384-well plates using a 10 µL POLQ Helicase ATPase Enzyme Reaction and 10 µL of detection/quench reagents (final volume 20 µL when the plates are read). The use of different plate densities or reaction volumes will require changes in reagent quantities (see **Section 5.1** for example reaction volumes).

The methods were optimized for initial velocity detection of ADP formation ($\leq 20\%$ conversion of ATP to ADP) by POLQ Helicase with ATP at the approximate K_m concentration (50 µM) and POLQ ssDNA at a saturating concentration (0.1 µM). These conditions will ensure sensitive detection of inhibitors that compete with ATP while minimizing the effects of compounds that inhibit POLQ via non-specific interactions with DNA. Significant changes in the ATP concentration may require optimization of the ADP² Antibody concentration (in the case of the FP and FI readout modes) or the ADP Tracer (in the case of the TR-FRET readout mode) to adjust the dynamic range as described in the Transcreener ADP² Assay Manual.

Note: Antibody (TR-FRET) and Tracer (FP and FI) concentrations remain constant in the 20 µL Complete Assay regardless of changes to other reaction conditions.

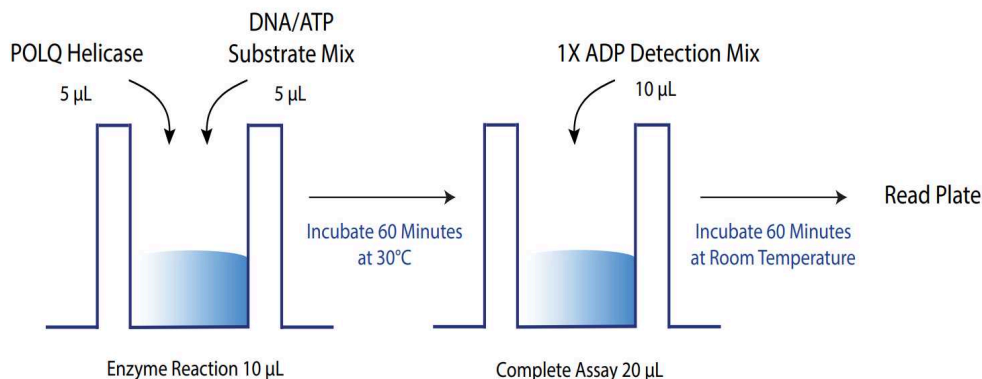


Figure 2. An Outline of the Procedure. The POLQ Helicase ATPase Enzyme Reaction is initiated by the addition of POLQ ssDNA and ATP. After the Enzyme Reaction incubation is completed, ADP detection reagents are added (Transcreener® ADP² Antibody and Tracer) along with EDTA to quench the POLQ Helicase ATPase reaction.

10 µL Enzyme Reaction Components

Component	Working Stock	Final Concentration in 10 µL
Enzyme Assay Buffer A, 10X	1X in Nuclease Free Water	1X (50 mM TRIS pH 7.5, 1 mM MgCl ₂ , and 0.01% Triton)
POLQ Helicase Enzyme, 0.1 mg/mL (1.1 µM)	2X in 1X Enzyme Assay Buffer A	250 pM – 500 pM*
ATP, 5 mM	100 µM in 1X Enzyme Assay Buffer A (with 0.2 µM POLQ ssDNA)	50 µM
POLQ ssDNA, 40 µM	0.2 µM in 1X Enzyme Assay Buffer A (with 100 µM ATP)	0.1 µM

*See Section 4.1 for Determining the Optimal Enzyme Concentration.

Table 1. POLQ Helicase ATPase Enzyme Reaction Components. Concentrations are provided for the standard protocol using 5 µL of POLQ Enzyme Mix and 5 µL of DNA/ATP Substrate Mix for the Enzyme Reaction.

Component	1X ADP Detection Mix - Add 10 µL Per Well					
	FP		FI		TR-FRET	
	Detection Mix Concentration	Complete Assay	Detection Mix Concentration	Complete Assay	Detection Mix Concentration	Complete Assay
Stop & Detect Buffer	1X	0.5X	1X	0.5X	1X	0.5X
ADP Tracer	4 nM	2 nM	8 nM	4 nM	500 nM	250 nM
ADP Antibody	55 µg/mL	27.5 µg/mL	50 µg/mL	25 µg/mL	8 nM	4 nM

Table 2. 1X ADP Detection Mix Components. The optimal concentrations for each of the detection reagents based on the preferred readout are shown. Changes to the concentrations may require re-optimization of the assay.

Determining the Optimal Enzyme Concentration

- 4.1 Using the enzyme concentration suggested in the POLQ Helicase Enzyme Certificate of Analysis should provide a robust signal that is within the linear range for ADP formation. However, for best results, we suggest performing an enzyme titration to identify the optimal enzyme concentration (EC_{50} to EC_{80}), especially when running the assay in a different buffer system or with a different ATP or POLQ ssDNA concentration. This example uses a 2X serial dilution; it should be performed at least in duplicate. If a compound screen is planned, you should include the solvent (e.g., DMSO) at its final assay concentration.

4.1.1 Enzyme Titration Steps

1. Prepare 1500 µL 1X Enzyme Assay Buffer A: dilute 150 µL of 10X Enzyme Assay Buffer A in 1350 µL Ultrapure Nuclease Free Water.
2. Prepare 344 µL of 16 nM POLQ Helicase Enzyme: dilute 5 µL of 1.1 µM POLQ Helicase Enzyme in 339 µL 1X Enzyme Assay Buffer A.
3. Add 10 µL of the POLQ Helicase Enzyme to well 1 (including replicates).
4. Add 5 µL of 1X Enzyme Assay Buffer A to wells 2-12, DO NOT add the Assay Buffer to well 1.
5. Transfer 5 µL from well 1 to well 2 and mix by pipetting, then transfer 5 µL from well 2 to well 3 and mix by pipetting; repeat this serial dilution process until well 11 has received POLQ Helicase Enzyme. Well 12 is to be used as a blank and should not include enzyme.
IMPORTANT: After mixing the last well (11) in the dilution series, remove 5 µL from that well only and discard, so that all the wells contain 5 µL final volume.
6. Prepare 500 µL of DNA/ATP Substrate Mix: dilute 2.5 µL 40 µM POLQ ssDNA and 10 µL 5 mM ATP in 487.5 µL 1X Enzyme Assay Buffer A.
7. Start the Enzyme Reaction by adding 5 µL of the DNA/ATP Substrate Mix to every well (1-12). Gently mix for 40 to 60 seconds on a plate shaker. Incubate at 30°C for 60 minutes.
8. Prepare 500 µL 1X ADP Detection Mix based on the concentrations provided in

Table 2: 50 μL 10X Stop & Detect Buffer B, 5 μL ADP² Alexa Fluor® 633 Tracer and 55 $\mu\text{g/mL}$ ADP² Antibody in Ultrapure Nuclease Free Water.

Note: The 1X ADP Detection Mix varies for the readout mode used. Table 2 lists the 1X ADP Detection Mix for each readout.

9. Add 10 μL of 1X ADP Detection mix to every well (1-12), in replicate.

10. Gently mix on a plate shaker for 40 to 60 seconds and then allow it to incubate at room temperature for 60 minutes before reading.

Note: The reagent volumes indicated above are sufficient for running the enzyme titration in duplicate plus excess for pipetting dead volume. Scaling of volumes can be performed if necessary.

For detection of inhibitors at single concentration or in dose response mode, we recommend selecting an enzyme concentration that produces a 50–80% change in signal (EC_{50} to EC_{80}) (see **Figure 3**). This will result in initial velocity conditions, which correspond to the linear phase of the reaction after conversion of values to ADP formation (see **Figure 6**). The EC_{50} is provided by common graphing programs; the EC_{80} enzyme concentration can be calculated from the EC_{50} , as follows:

$$\text{EC}_X = (X \div (100 - X))^{(1 \div |\text{hillslope}|)} \times \text{EC}_{50}$$

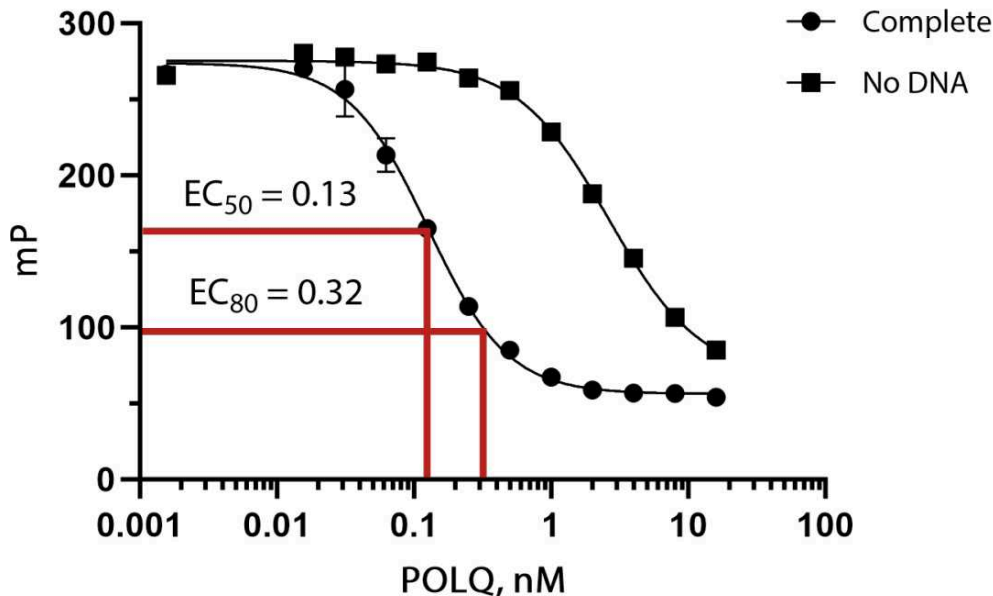


Figure 3. Enzyme Titration Curve. Example POLQ Helicase enzyme titration. The ideal range of enzyme concentrations is between EC_{50} and EC_{80} ; the specific concentration may vary depending on the enzyme lot.

Performing Single Compound Screening and Dose-Response Assays

4.2 4.2.1 Experimental Samples

1. Perform a serial dilution of test compounds with your method of choice. Add the enzyme to the test compounds at the desired concentration so that the total volume of this mixture is 5 μ L. Mix gently on a plate shaker for 40 to 60 seconds. Preincubate the Enzyme Inhibitor Mix for the desired time (typically at least 15 minutes) at room temperature to allow equilibration of the E-I complex

Note: Final concentration of test compounds should be based on the volume of the Enzyme Reaction.

2. Start the Enzyme Reaction by adding 5 μ L of the DNA/ATP Substrate Mix. It is recommended to incubate the Enzyme Reaction at 30°C for 60 minutes.

Note: The final volume of the Enzyme Reaction mixture should be 10 μ L for 384 well plates. See Section 5.1 for a list of other plate formats.

3. After the incubation, add 10 μ L of 1X ADP Detection Mix to the 10 μ L Enzyme Reaction and mix the 20 μ L Complete Assay using a plate shaker.

Note: The 1X ADP Detection Mix varies for the readout mode used. Table 2 lists the 1X ADP Detection Mix for each readout.

4. Incubate at room temperature (20–25°C) for 60 minutes and measure FP.

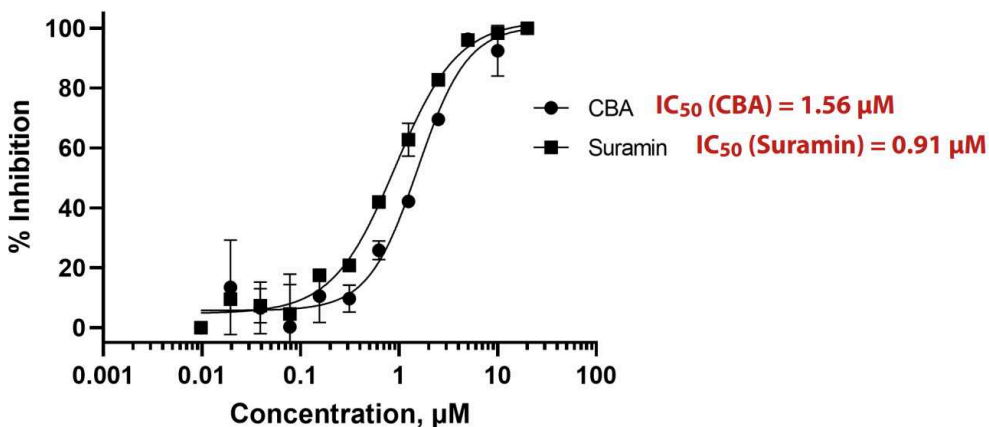
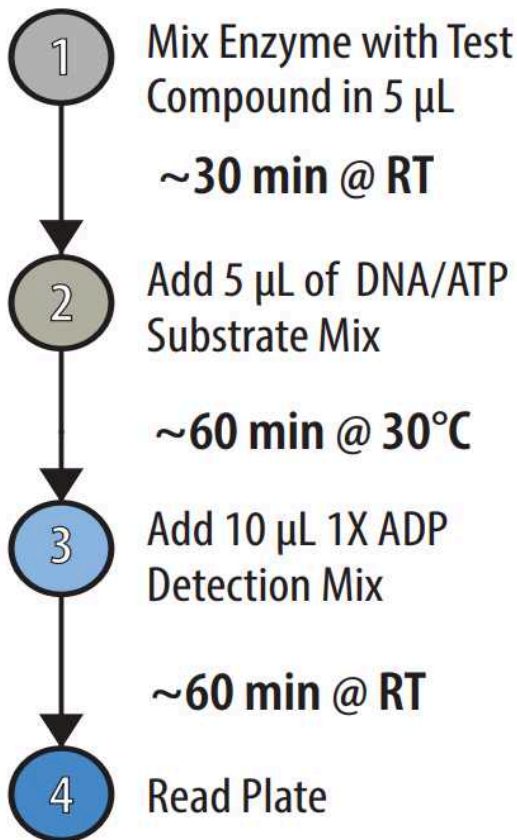


Figure 4. Dose-Response Curve. Example titration of probe inhibitors CBA and Suramin in the presence of POLQ Helicase Enzyme.

Setting Up a Standard Curve

4.3 Use of a standard curve for conversion of values to amount of ADP formed allows quantitative measurement of the enzyme activity and accurate IC₅₀ determinations; it is not typically done for screening at single concentrations. The standard curve mimics the Enzyme Reaction (as ATP concentration decreases, ADP concentration increases); the adenine nucleotide concentration remains constant. The ADP/ATP standard curve allows calculation of the concentration of ADP produced in the Enzyme Reaction and, therefore, the percent ADP conversion.

In this example, a 12-point standard curve was prepared using the concentrations of ADP and ATP shown in the following Table. Commonly, 8- to 12-point standard curves are used. Here we describe preparation of a standard curve from 50 μ M to 0.25 μ M ADP, which encompasses the appropriate range for this assay.

	% Conversion	ATP (μ M)	ADP (μ M)
	100	0	50
	50	25	25
	25	37.5	12.5
	15	42.5	7.5
	10	45	5
	7.5	46.25	3.75
	5	47.5	2.5
	3	48.5	1.5
	2	49	1
	1	49.5	0.5
	0.5	49.75	0.25
	0	50	0

Table 3. Concentration of ATP and ADP to prepare a 12-point standard curve

4.3.1 Standard Curve Steps

1. Prepare 1X Enzyme Assay Buffer A.
2. Dilute 5 mM ATP and ADP to 50 μ M in 1X Enzyme Assay Buffer A. The volume for both dilutions should be based on the total amount of volume to be utilized during preparation of the ATP/ADP percent conversion solutions.
3. Starting from 50 μ M of ADP, add proportional values of ATP and ADP with the

method of choice (i.e., Automatic Dispenser). For manual preparations, add reagents in separate microcentrifuge tubes to generate the desired percent conversions utilized in the standard curve. An example of the dilution scheme can be found in **Table 3**. Once all dilutions are completed, add 10 μL of each solution to the respective wells.

4. Afterwards, add 10 μL of 1X ADP Detection Mix. The 1X ADP Detection Mix varies for the readout mode used. **Table 2** lists the 1X ADP Detection Mix for each readout.

5. Finally, mix for 40 to 60 seconds, before measuring in the instrument of choice.

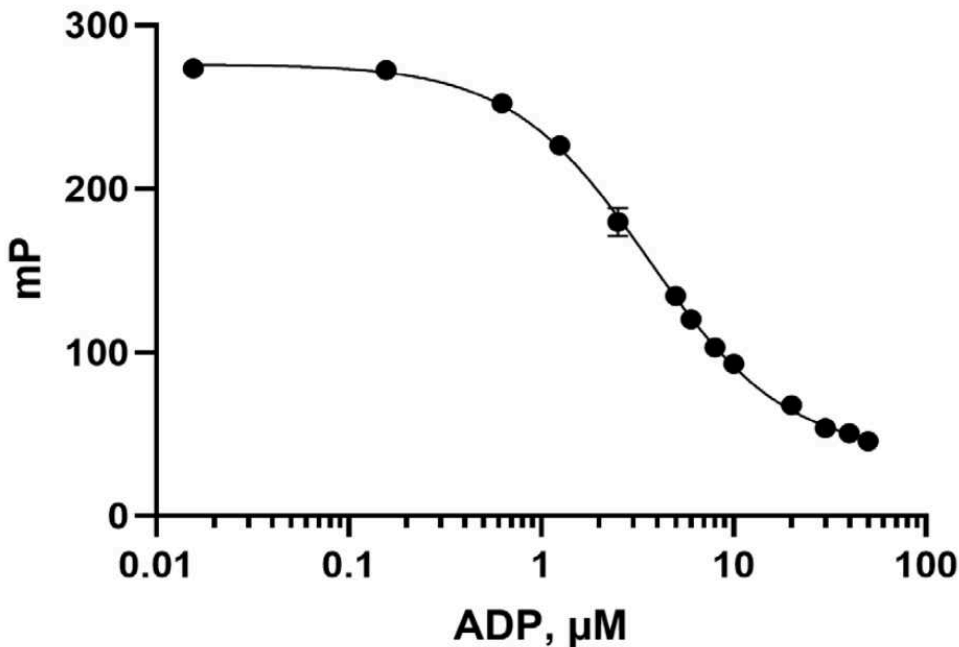


Figure 5. ADP Standard Curve. Standard curve using 1X ADP Detection Mix as shown in **Table 2**. (55 $\mu\text{g}/\text{mL}$ ADP² antibody, 4 nM ADP² Tracer)

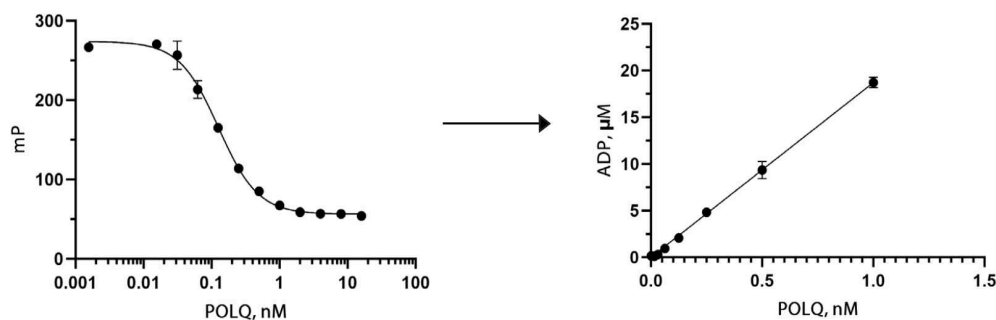


Figure 6. Enzyme Titration Curve Converted to ADP Formed. Raw polarization signal (mP) is converted to ADP formed using a standard curve as described in **Section 4.3**. Only the linear portion of the graph is shown; interpolation was performed using GraphPad Prism.

Measuring Assay Robustness with Z'

- 4.4 By taking into account both dynamic range and data variability at the high and low ranges of the assay, the Z' statistic provides a measure of what is of most interest when considering the suitability of an assay for HTS: the usable screening or "assay window." It is a dimensionless coefficient for the quality of the screening window that is relevant for any assay, regardless of detection method or readout, without the intervention of test compounds. As a guideline, a Z' value of 0.5 or greater is generally considered to be indicative of a very good screening window for a biochemical assay, thus the assay is an excellent assay. When running the POLQ Helicase ATPase Assay, run the controls with and without enzyme (no test compound) to achieve final results. Use the following formula to determine Z':

$$Z' = 1 - \frac{[(3 \times SD_{\text{No Enzyme}}) + (3 \times SD_{\text{Complete Reaction}})]}{|(\text{Mean}_{\text{No Enzyme}}) - (\text{Mean}_{\text{Complete Reaction}})|}$$

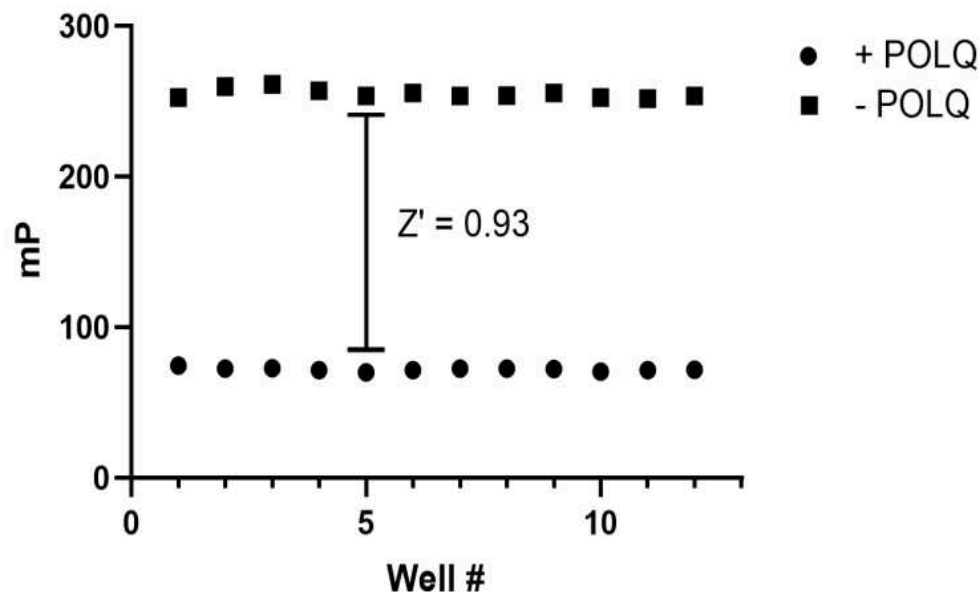


Figure 7. Z' Measurement. Complete Assay is performed with and without POLQ Helicase Enzyme (n=12). Z' is then calculated based on the formula shown in **Section 4.4**.

Appendix

5

Using the Assay with Different Volumes and Plate Formats

5.1

	Component	Total Volume	Enzyme Reaction Volume	1X ADP Detection Mixture Volume
	96 Well Low Volume Plate	50 µL	25 µL	25 µL
	384 Well Low Volume Plate	20 µL	10 µL	10 µL
	1536 Well Low Volume Plate	8 µL	4 µL	4 µL

Please check the working plate volumes from the manufacturer to ensure they are within the suggest volumes ranges of your plate.

Links to Applicable Application Notes

- 5.2
- [A Guide to Navigating Hit Prioritization After Screening Using Biochemical Assays](#)
 - [A Guide to Measuring Drug-Target Residence Times with Biochemical Assays](#)
 - [Using GraphPad Prism to Interpolate Data from a Standard Curve to Generate a Dose-Response](#)
 - [List of Commonly Used Plate Readers and Settings](#)

Contact Information

6

Email: TechSupport@bellbrooklabs.com

Phone: 608.443.2400

Toll-Free: 866.313.7881

FAX: 608.441.2967