

Nov 10, 2025

Enzolution™ TREX1 Assay System Technical Manual

DOI

<https://dx.doi.org/10.17504/protocols.io.bp2l6zz8dgqe/v1>

info¹

¹BellBrook Labs LLC

BellBrook Labs

Tech. support phone: +1 (608) 443-2400 email: techsupport@bellbrooklabs.com



info info

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bp2l6zz8dgqe/v1>

Protocol Citation: info 2025. Enzolution™ TREX1 Assay System Technical Manual. protocols.io <https://dx.doi.org/10.17504/protocols.io.bp2l6zz8dgqe/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 23, 2025

Last Modified: November 10, 2025

Protocol Integer ID: 230624

Keywords: enzymatic activity for trex1, purified human trex1, trex1 inhibitor, trex1 concentration, adjustment of the trex1 concentration, human trex1, degrading cytosolic dna, trex1, fp assay kit, cytosolic dna, dna substrate, stranded dna, stimulatory dna, preventing aberrant autoimmune response, ifn response in mammalian cell, assay, inhibitor dose response measurement, 45bp dsdna oligomer, interferon stimulatory dna, sting pathway, prime repair exonuclease, enzymatic activity, deoxynucleoside monophosphate, insensitive to dna, dna, releasing deoxynucleoside monophosphate, high throughput screening, 45bp dsdna oligomer from the listeria



Abstract

The Enzolution™ TREX1 Assay System is intended for use with the Transcreener® dAMP FP Assay Kit (Part #3028) to measure enzymatic activity for TREX1 (also known as Three Prime Repair Exonuclease 1). TREX1 is the major exonuclease responsible for degrading cytosolic DNA, and it plays an important role in preventing aberrant autoimmune responses in normal cells by acting as a checkpoint for the cGAS/STING pathway. It is a 3'-5' exonuclease that cleaves single bases from the 5' termini of double stranded or single stranded DNA, releasing deoxynucleoside monophosphates, including deoxy-AMP (dAMP). The Transcreener® dAMP FP assay enables sensitive detection of dAMP in a single addition, mix-and-read format that has been optimized and extensively validated for high throughput screening (HTS) and inhibitor dose response measurements using most multimode plate readers.

The Enzolution™ TREX1 Assay System provides all reagents required to screen and profile TREX1 inhibitors when used with the Transcreener® dAMP FP Assay Kit, including purified human TREX1 (amino acids 1-286, C-terminal 6xHis) and DNA Substrate. TREX1 is relatively non-specific for its DNA Substrate; the Enzolution™ TREX1 Assay System uses interferon stimulatory DNA (ISD), a 45bp dsDNA oligomer from the *Listeria monocytogenes* genome that induces a strong type I IFN response in mammalian cells. Note that we have optimized the assay using a saturating concentration of ISD, which makes the assay relatively insensitive to DNA-competitive inhibitors; use of lower ISD concentrations will require adjustment of the TREX1 concentration accordingly. Also, the protocol is configured for 384-well plates; use of different multiwell plate formats will require some additional optimization.


Introduction

- 1 The Enzolution™ TREX1 Assay System is intended for use with the Transcreener® dAMP FP Assay Kit (Part #3028) to measure enzymatic activity for TREX1 (also known as Three Prime Repair Exonuclease 1). TREX1 is the major exonuclease responsible for degrading cytosolic DNA, and it plays an important role in preventing aberrant autoimmune responses in normal cells by acting as a checkpoint for the cGAS/STING pathway. It is a 3'-5' exonuclease that cleaves single bases from the 5' termini of double stranded or single stranded DNA, releasing deoxynucleoside monophosphates, including deoxy-AMP (dAMP). The Transcreener® dAMP FP assay enables sensitive detection of dAMP in a single addition, mix-and-read format that has been optimized and extensively validated for high throughput screening (HTS) and inhibitor dose response measurements using most multimode plate readers.

The Enzolution™ TREX1 Assay System provides all reagents required to screen and profile TREX1 inhibitors when used with the Transcreener® dAMP FP Assay Kit, including purified human TREX1 (amino acids 1-286, C-terminal 6xHis) and DNA Substrate. TREX1 is relatively non-specific for its DNA Substrate; the Enzolution™ TREX1 Assay System uses interferon stimulatory DNA (ISD), a 45bp dsDNA oligomer from the *Listeria monocytogenes* genome that induces a strong type I IFN response in mammalian cells. Note that we have optimized the assay using a saturating concentration of ISD, which makes the assay relatively insensitive to DNA-competitive inhibitors; use of lower ISD concentrations will require adjustment of the TREX1 concentration accordingly. Also, the protocol is configured for 384-well plates; use of different multiwell plate formats will require some additional optimization.

Key Applications:

- Screening for TREX1 inhibitors
- Generating dose response curves and IC₅₀ values for TREX1 inhibitors
- Kinetic and mechanistic analyses

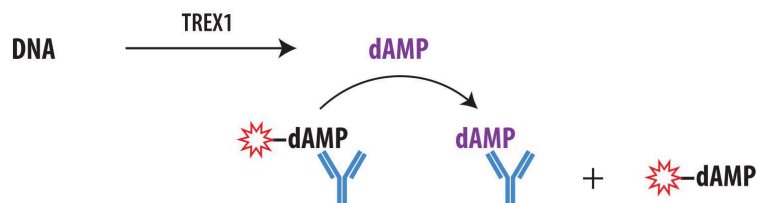

Product Specifications

- 2

Product	Quantity	Part #
Enzolution™ TREX1 Assay System	1,000 assays*	3029-1K
	10,000 assays*	3029-10K

*The exact number of assays depends on the enzyme reaction conditions. The kits are designed for use with 384-well plates, using 10 µL Enzyme Reaction and 20 µL Complete Assay volumes.

**Storage**

TREX1 should be stored at -80°C; other reagents can be stored at -20°C. Though we have confirmed that TREX1 is stable up to 2 freeze-thaw cycles, we recommend aliquoting the enzyme and snap-freezing for multiple uses to minimize loss of activity.

Use the reagents provided in this kit within 6 months from date of receipt.

Materials Provided

2.1

Component	Composition	Notes
TREX1 Enzyme	0.1 mg/mL (3.2 µM)* solution in 50 mM TRIS (pH 8.0), 500 mM NaCl with 20% glycerol	Amino acids 1-286, C-terminal 6xHis, 31.3 kDa. Sufficient enzyme is included in the kit to complete 1,000 assays (Part # 3029-1K) or 10,000 assays (Part # 3029-10K).
Interferon Stimulatory DNA	25 µM in H ₂ O	Supplied as a 45-bp annealed duplex interferon stimulatory DNA (ISD) oligomer.
TREX1 Assay Buffer, 10X	100 mM TRIS (pH 7.5), 75 mM MgCl ₂ , 0.05% BSA, and 0.1% Brij-35	Use the TREX1 Assay Buffer in the Enzyme Reaction and for inhibitor incubation. Changes to the assay buffer could affect TREX1 activity and/or detection of dAMP.
384-Well Low Volume Black Assay Plates	Corning #4514	Black polystyrene non binding assay plates in either a 3-pack (1,000+ Assays) or a 30-pack (10,000+ Assays). We strongly recommend the use of these plates as inconsistent results have been observed with other plate formats.

*The exact concentration may vary from batch to batch. Please refer to the Certificate of Analysis for an accurate concentration.

Materials Required But Not Provided

2.2

Component	Notes
Ultrapure Nuclease Free Water	Some deionized water systems are contaminated with nucleases that can degrade both nucleotide substrates and products, reducing assay performance. Use nuclease free water such as: Invitrogen Part # AM9930
Plate Reader	A multidetection microplate reader configured to measure FP of the dAMP Alexa Fluor® 633 tracer is required. Transcreener® FP Assays have been successfully used on the following instruments: BioTek Synergy™2 and Synergy™4; BMG Labtech PHERAstar® Plus and CLARIOstar® Plus; Molecular Devices SpectraMax™ Paradigm; Perkin Elmer EnVision® and ViewLux; and Tecan Infinite® F500, Safire2™, and M1000. Full list of compatible plate readers and settings.
Liquid Handling Devices	Use liquid handling devices that can accurately dispense a submicroliter volumes into 384-well plates.
Laboratory Incubator	An incubator model that is capable of maintaining temperature stability at 30°C is required.

Transcreener® dAMP Exonuclease Assay - SOLD SEPARATELY

Component	Composition	Notes
dAMP Antibody	3.5 mg/mL solution in PBS with 5% glycerol*	Mouse polyclonal antibody. Sufficient antibody is included in the kit to complete 1,000 assays (Part # 3028-1K) or 10,000 assays (Part # 3028-10K).
dAMP Alexa Fluor 633 Tracer	400 nM solution in 2 mM HEPES (pH 7.5) containing 0.01% Brij-35	Sufficient tracer is included in the kit to complete 1,000 assays (Part # 3028-1K) or 10,000 assays (Part # 3028-10K).
Stop & Detect Buffer B, 10X	200 mM HEPES (pH 7.5), 400 mM EDTA, and 0.2% Brij-35	The EDTA in the Stop & Detect Buffer B quenches the Enzyme Reaction by chelating Mg^{2+} . Therefore, the assay should work as an end-point assay as long as the EDTA is at least equimolar to the Mg^{2+} . The final concentrations of Mg^{2+} and EDTA in the Complete Assay are 7.5 mM and 20 mM, respectively.
dAMP	5 mM in H ₂ O (pH 7.0)	The dAMP in this kit can be used to generate a standard curve which can be used to convert mP values to dAMP product formed.

Before You Begin

1. Read the entire protocol and note any reagents or equipment needed (see **Section 2.2**).
2. Check the FP instrument and verify that it is compatible with the assay being performed (see **Transcreener® dAMP Assay Technical Manual Section 4.1**).
3. Please read and understand the Transcreener® dAMP Assay Technical Manual prior to using with this kit.

Protocol

4. The methods described below are for single-addition, endpoint detection in 384 well plates using 10 μ L TREX1 reactions and 10 μ L of detection/quench reagents (final volume 20 μ L when the plates are read). TREX1 reactions are initiated by the addition of the ISD Substrate and quenched by the addition of EDTA in the dAMP Detection Mix (see **Figure 2**) (TREX1 is a Mg-dependent enzyme). The use of different plate densities or reaction volumes will require changes in reagent quantities (see **Section 5.1** for example reaction volumes). Additionally, the methods were designed for initial velocity detection of dAMP formation by TREX1 over a range of 0.2 to 20 μ M dAMP; if desired, the dynamic range can be changed by adjusting the concentration of dAMP Ab, as described in the Transcreener® dAMP FP Assay Technical Manual.

10 μ L Enzyme Reaction Components

Component	As Provided	Working Stock	Final Concentration in 10 μ L
TREX1 Assay Buffer, 10X	10X	1X, Nuclease Free Water	1X (10 mM TRIS pH 7.5, 7.5 mM MgCl ₂ , 0.005% BSA, and 0.01% Brij-35)
TREX1 Enzyme	0.1 mg/mL (3.2 μ M)	2X in 1X TREX1 Assay Buffer	1X (0.1 nM - 1.2 nM)
Interferon Stimulatory DNA	25 μ M	0.70 μ M in 1X TREX1 Assay Buffer	0.35 μ M

Table 1. TREX1 Enzyme Reaction Components. Concentrations are provided for the standard protocol using 5 μ L of TREX1 Enzyme Mix and 5 μ L of ISD Substrate Mix for the Enzyme Reaction.

1X dAMP Detection Mix - Add 10 μ L Per Well

Component	As Provided	Detection Mix Concentration	Final Concentration in 20 μ L Complete Assay	Example Amounts for 10 mL Stock
dAMP Antibody*	3.5 mg/mL	20 μ g/mL	10 μ g/mL	57.1 μ L
dAMP Alexa Fluor® 633 Tracer	400 nM	2 nM	1 nM	50.0 μ L
Stop & Detect Buffer B, 10X	10X	1X	0.5X	1,000.0 μ L
Nuclease Free Water	-	-	-	8,892.9 μ L

*The exact concentration may vary from batch to batch. Please refer to the Certificate of Analysis for an accurate concentration.

Table 2. 1X dAMP Detection Mix Components. Volumes provided in the table are based on preparation of a 10 mL solution; adjust these appropriately for the desired volumes, including 10% extra for pipetting dead volumes.

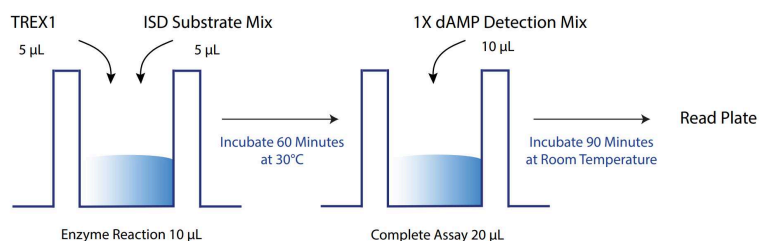


Figure 2. An Outline of the Procedure. The TREX1 Enzyme Reaction is initiated by the addition of ISD. After the Enzyme Reaction incubation is completed, dAMP detection reagents are added (Transcreener® dAMP Antibody and Tracer) along with EDTA to quench the TREX1 reaction.

Setting Up a Standard Curve

- 4.1 Use of a standard curve for conversion of mP values to amount of dAMP formed allows quantitative measurement of TREX1 activity and accurate IC₅₀ determinations; it is not typically done for screening at single concentrations. Here we describe preparation of a standard curve using 3X serial dilution from 200 μ M to 0.0034 μ M dAMP, which encompasses the appropriate range for this application.

Note: These concentrations represent the dAMP levels in the 10 μ L TREX1 Enzyme Reaction; they will be diluted 2-fold by addition of detection reagents. The reagent volumes indicated below are sufficient for running the standard curve in duplicate plus excess for pipetting dead volume.

4.1.1 Detailed Methods

1. Prepare 600 μ L 1X TREX1 Assay Buffer: dilute 60 μ L TREX1 Assay Buffer, 10X in 540

μL Ultrapure Nuclease Free Water.

2. Prepare 30 μL of 400 μM dAMP: dilute 2.4 μL 5 mM dAMP stock in 27.6 μL 1X TREX1 Assay Buffer.
3. Prepare 160 μL of ISD Substrate Mix: dilute 4.5 μL Interferon Stimulatory DNA in 155.5 μL 1X TREX1 Assay Buffer.
4. Prepare 300 μL 1X dAMP Detection Mix based on the concentrations provided in **Table 2**: 30 μL 10X Stop & Detect Buffer B, 1.5 μL dAMP Alexa Fluor® 633 Tracer and 20 μg/mL dAMP Antibody in Ultrapure Nuclease Free Water.

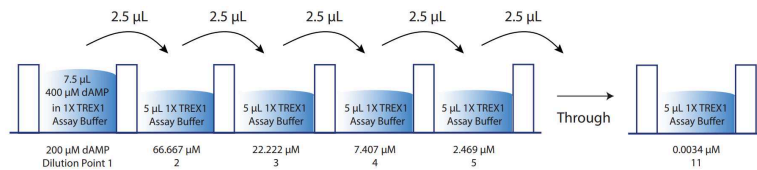


Figure 3. Performing a Serial Dilution. Example 3-fold serial dilution of dAMP to generate a standard curve.

5. Add 7.5 μL of the 400 μM dAMP to well 1 (including replicates).
6. Add 5 μL of 1X TREX1 Assay Buffer to wells 2-12, **DO NOT** add the 1X TREX1 Assay Buffer to well 1.
7. Transfer 2.5 μL from well 1 to well 2 and mix by pipetting, then transfer 2.5 μL from well 2 to well 3 and mix by pipetting; repeat this serial dilution process until well 11 has received dAMP. Well 12 is to be used as a blank and should correspond to 0 μM of dAMP on the standard curve.

IMPORTANT: After mixing the last well in the dilution series, remove 2.5 μL from that well only and discard, so that all the wells contain 5 μL final volume.

8. Add 5 μL of the ISD Substrate Mix to every well (1-12) then add 10 μL of 1X dAMP Detection mix to every well (1-12).
9. Gently mix on a plate shaker for 40 to 60 seconds and then allow it to incubate at room temperature for 90 minutes before reading.

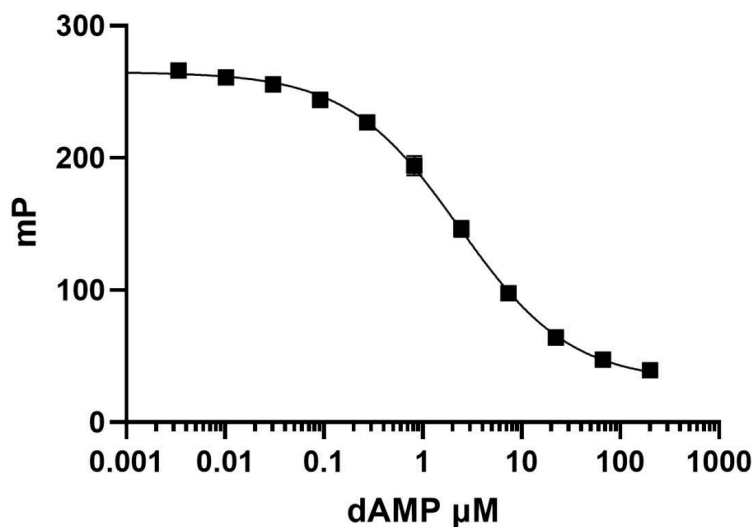


Figure 4. dAMP Standard Curve. Standard curve using 1X dAMP Detection Mix as described in **Step 4** and **Table 2**.

Determining the Optimal Enzyme Concentration

- 4.2 Using EC₈₀ enzyme concentration shown in the TREX1 Enzyme Certificate of Analysis should provide a robust signal that is within the linear range for dAMP formation. However, for best results, we suggest performing an enzyme titration to identify the optimal enzyme concentration, especially when running the assay in a different buffer system or with a different substrate concentration. The enzyme titration should be performed in duplicate and this example uses a 2X serial dilution. If a compound screen is planned, you should include the solvent (e.g., DMSO) at its final assay concentration. The Transcreener® dAMP FP Assay Kit has been shown to be able to tolerate up to 5% DMSO.

Note: TREX1 is a labile protein that is easily denatured; rapid or prolonged mixing should be avoided to preserve enzymatic activity.

4.2.1 Detailed Methods

1. Prepare 600 µL 1X TREX1 Assay Buffer: dilute 60 µL TREX1 Assay Buffer, 10X in 540 µL Ultrapure Nuclease Free Water.
 2. Prepare 160 µL of 20 nM TREX1 Enzyme: dilute 1 µL TREX1 Enzyme in 159 µL 1X TREX1 Assay Buffer with gentle mixing.
 3. Prepare 160 µL of ISD Substrate Mix: dilute 4.5 µL Interferon Stimulatory DNA in 155.5 µL 1X TREX1 Assay Buffer.
 4. Add 10 µL of the 20 nM TREX1 to well 1 (including replicates).
 5. Add 5 µL of 1X TREX1 Assay Buffer to wells 2-12, DO NOT add the 1X TREX1 Assay Buffer to well 1.
 6. Transfer 5 µL from well 1 to well 2 and mix by pipetting, then transfer 5 µL from well 2 to well 3 and mix by pipetting; repeat this serial dilution process until well 11 has received TREX1 Enzyme. Well 12 is to be used as a blank and should not include enzyme.
- IMPORTANT: After mixing the last well (11) in the dilution series, remove 5 µL from that well only and discard, so that all the wells contain 5 µL final volume.**
7. Start the enzyme reaction by adding 5 µL of the ISD Substrate Mix to every well (1-12). Gently mix for 40 to 60 seconds on a plate shaker. Incubate at 30°C for 60 minutes.
 8. Prepare 300 µL 1X dAMP Detection Mix based on the concentrations provided in **Table 2**: 30 µL 10X Stop & Detect Buffer B, 1.5 µL dAMP Alexa Fluor® 633 Tracer and 20 µg/mL dAMP Antibody in Ultrapure Nuclease Free Water.
 9. Add 10 µL of 1X dAMP Detection mix to every well (1-12), in replicate.
 10. Gently mix on a plate shaker for 40 to 60 seconds and then allow it to incubate at room temperature for 90 minutes before reading.

Note: The reagent volumes indicated above are sufficient for running the enzyme titration in duplicate plus excess for pipetting dead volume. Scaling of volumes can be performed if necessary.

For detection of inhibitors at single concentration or in dose response mode, we recommend selecting an enzyme concentration that produces an 50-80% change in FP signal (EC₅₀ to EC₈₀) (see **Figure 5**) and an assay window of at least 100 mP. This will result in initial velocity conditions, which correspond to the linear phase of the reaction after conversion of mP values to dAMP formation (see **Figure 6**). The EC₅₀ is provided by common graphing programs; the EC₈₀ enzyme concentration can be calculated from the EC₅₀, as follows:

$$EC_x = (X \div (100 - X))^{(1 \div |hillslope|)} \times EC_{50}$$

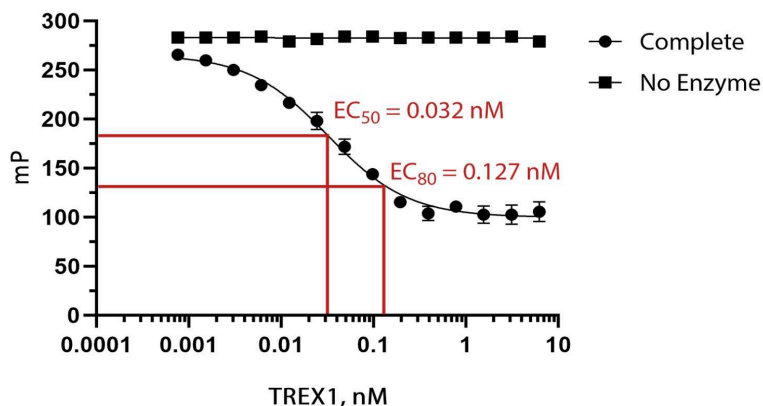


Figure 5. Enzyme Titration Curve. Example TREX1 Enzyme titration. The ideal range of enzyme concentrations is between EC_{50} and EC_{80} ; the specific concentration may vary depending on the enzyme lot.

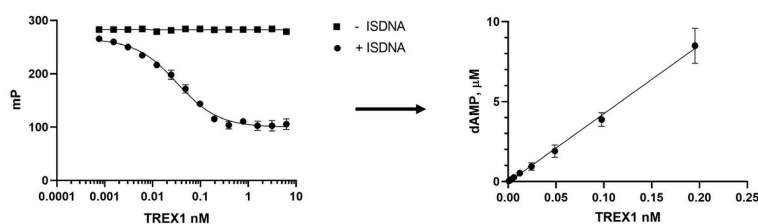


Figure 6. Enzyme Titration Curve converted to dAMP formed. Raw polarization signal (mP) is converted to dAMP formed using a standard curve as described in **Section 4.1**. Only the linear portion of the graph is shown; interpolation was performed using GraphPad Prism.

Performing Single Compound Screening and Dose-Response Assays

4.3 4.3.1 Detailed Methods

1. Perform a serial dilution of test compounds with your method of choice. Add the TREX1 Enzyme to the test compounds at the desired concentration so that the total volume of this mixture is 5 μ L. Mix gently on a plate shaker for 40 to 60 seconds. Preincubate the Enzyme Inhibitor Mix for the desired time (typically at least 30 minutes) at room temperature to allow equilibration of the E-I complex.

Note: Final concentration should be based on the volume of the Enzyme Reaction.

2. Start the enzyme reaction by adding 5 μ L of the ISD Substrate Mix. It is recommended to incubate the enzyme reaction at 30°C for 60 minutes.

Note: The final volume of the enzyme reaction mixture should be 10 μ L for 384 well plates. See **Section 5.1** for a list of other plate formats.

3. Following the 60 minute incubation, add 10 μ L of 1X dAMP Detection Mix to the 10 μ L Enzyme Reaction and mix the 20 μ L Complete Assay using a plate shaker for 40 to 60 seconds.

4. Incubate at room temperature (20–25°C) for 90 minutes and measure FP

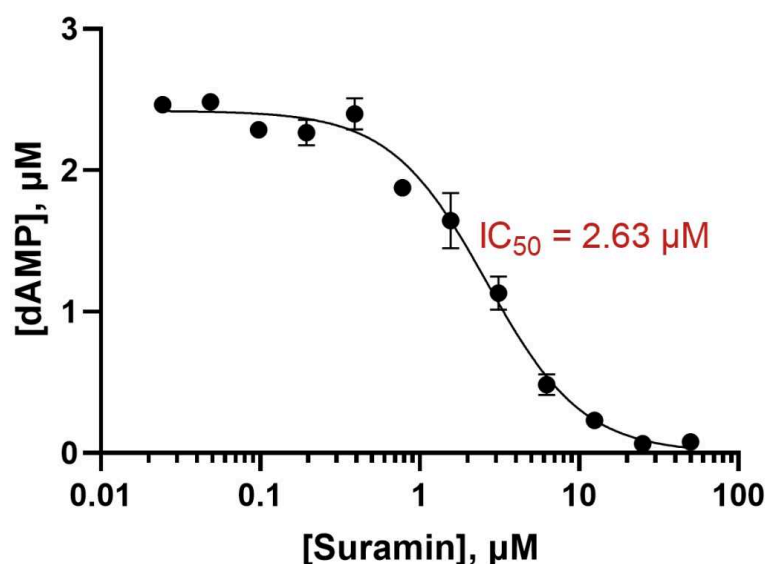
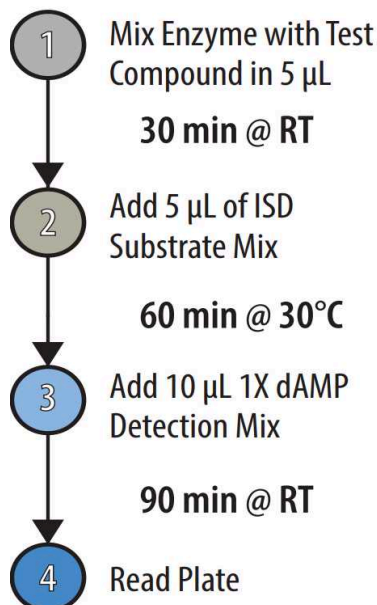


Figure 7. Dose-Response Curve. Example dose response curve with probe inhibitor Suramin.

Measuring Assay Robustness with Z'

- 4.4 By taking into account both dynamic range and data variability at the high and low ranges of the assay, the Z' statistic provides a measure of the overall quality of an assay for HTS based on the usable signal or "assay window." It is a dimensionless coefficient for the quality of the assay that is relevant for any assay, regardless of detection method or readout, without the intervention of test compounds. As a guideline a Z' value of 0.5 or greater is generally considered to be indicative of a very good screening window for a biochemical assay. Z' is determined using following formula, where complete reactions are untreated TREX1 enzymatic assays and No Enzyme are negative controls containing

all components except TREX1; assays should be performed at n =16 for statistical significance.

$$Z' = 1 - \frac{[(3 \times SD_{\text{No Enzyme}}) + (3 \times SD_{\text{Complete Reaction}})]}{|(\text{Mean}_{\text{No Enzyme}}) - (\text{Mean}_{\text{Complete Reaction}})|}$$

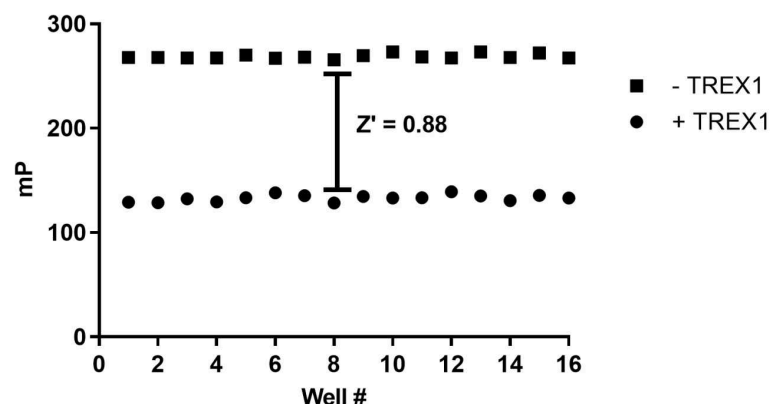


Figure 8. Z' Measurement. Complete assay is performed with and without TREX1 (n=16). Z' is then calculated based on the formula from 4.4.1.

Appendix

5

Using the Assay with Different Volumes and Plate Formats

5.1

Component	Total Volume	Enzyme Reaction Volume	dAMP Detection Mixture Volume
96 Well Low Volume Plate	50 µL	25 µL	25 µL
384 Well Low Volume Plate	20 µL	10 µL	10 µL
1536 Well Low Volume Plate	8 µL	4 µL	4 µL

Links to Applicable Application Notes

- 5.2
- [A Guide to Navigating Hit Prioritization After Screening Using Biochemical Assays](#)
 - [A Guide to Measuring Drug-Target Residence Times with Biochemical Assays](#)

- [List of Commonly Used Plate Readers and Settings](#)

Contact Information

6 **Email:** TechSupport@bellbrooklabs.com
 Phone: 608.443.2400
 Toll-Free: 866.313.7881
 FAX: 608.441.2967