


Oct 07, 2025

Enzyme-linked immunosorbent assay using the SARS-CoV-2 membrane glycoprotein analog M1s (sequence: ac-CTITVEELKKLLEQC-am) for detection of cognate COVID-19 antibodies as performed on high binding plates

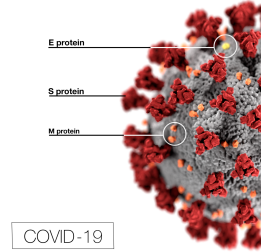

DOI

<https://dx.doi.org/10.17504/protocols.io.36wggpezyvk5/v1>

Brian Andrich Pollo¹, Danica S. Ching^{2,1}, Maria Isabel C. Idolor¹, Ruby Anne King³,
Fresthel Monica M. Climacosa^{4,1}, Salvador Eugenio C. Caoili¹

¹Biomedical Innovations Research for Translational Health Science (BIRTHS) Laboratory, Department of Biochemistry and Molecular Biology, University of the Philippines Manila;

²Institute of Clinical Epidemiology, National Institutes for Health, Manila;

³Department of Science and Technology - Philippine Council for Health Research and Development (DOST-PCHRD);

⁴Department of Medical Microbiology, College of Public Health, University of the Philippines Manila

Brian Andrich Pollo: corresponding author;

Coronavirus Method De...



Brian Andrich L. Pollo

University of the Philippines Manila

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.36wggpezyvk5/v1>

External link: <https://doi.org/10.6084/m9.figshare.30276484.v1>

Protocol Citation: Brian Andrich Pollo, Danica S. Ching, Maria Isabel C. Idolor, Ruby Anne King, Fresthel Monica M. Climacosa, Salvador Eugenio C. Caoili 2025. Enzyme-linked immunosorbent assay using the SARS-CoV-2 membrane glycoprotein analog M1s (sequence: ac-CTITVEELKKLLEQC-am) for detection of cognate COVID-19 antibodies as performed on high binding plates. [protocols.io https://dx.doi.org/10.17504/protocols.io.36wggpezyvk5/v1](https://dx.doi.org/10.17504/protocols.io.36wggpezyvk5/v1)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 06, 2025

Last Modified: October 07, 2025

Protocol Integer ID: 229142

Keywords: ELISA, SARS-CoV-2, COVID-19 antibodies, peptide antigen, membrane glycoprotein, serology, immunoassay, Protein A-peroxidase, TMB substrate, seroprevalence, immune monitoring, assay development, synthetic peptide, antibody detection, detecting antibody, using synthetic peptide antigen, linked immunosorbent assay, synthetic peptide antigen, peptide antigen in carbonate, immunosorbent assay, antibodies against the sar, antigen, antibody level, distinct absorbance profiles proportional to antibody level, immune response monitoring, synthetic peptide sequence, incubation with diluted serum, elisa, validation for sar, diluted serum

Funders Acknowledgements:

Dissertation grant

IDC211 grant

Abacov project

NIH faculty grant

CIP fellowship



Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to **protocols.io** is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with **protocols.io**, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

This protocol describes a peptide-based indirect enzyme-linked immunosorbent assay (ELISA) for detecting antibodies against the SARS-CoV-2 membrane glycoprotein analog M1s (synthetic peptide sequence: ac-CTITVEELKKLLEQC-am). High-binding polystyrene plates are coated overnight with the peptide antigen in carbonate-bicarbonate buffer, followed by blocking, incubation with diluted serum, detection using Protein A-peroxidase, and colorimetric readout with TMB substrate at 450 nm. The method offers a simple, reproducible, and cost-effective workflow for COVID-19 serological testing using synthetic peptide antigens. Expected results include distinct absorbance profiles proportional to antibody levels, supporting applications in seroprevalence analysis, immune response monitoring, and assay validation for SARS-CoV-2 antibody research.

Image Attribution

CDC/ Alissa Eckert, MS; Dan Higgins, MAM, Public domain, via Wikimedia Commons



Coat

- 1 overnight incubation at 4°C
100µL/well, Carbonate-bicarbonate buffer in water, pH 9.6

Wash

- 2 3x at 5 min interval
0.05% Tween 20 in 1X PBS (PBST)
freeze

Block

- 3 *leave plates at RT for 15min before blocking*
incubation at 37°C for 30 mins
200µL/well, 5% skimmed milk in PBST

Wash

- 4 3x at 5 min interval
0.05% Tween 20 in 1X PBS

Primary antibody

- 5 incubation at RT for 1hr
100µL/well, 1:100 sera in dilution buffer (0.05% skimmed milk in PBST)

Wash

- 6 3x at 5 min interval
0.05% Tween 20 in 1X PBS

Secondary antibody

- 7 incubation at RT for 1hr
50µL/well, 1:2000 Protein A-peroxidase (0.5ug/mL) in dilution buffer

Wash



- 8 3x at 5 min interval
0.05% Tween 20 in 1X PBS

Chromogenic substrate solution

- 9 incubation at RT for 30 min
50µL/well, 0.1mg/mL TMB in Phosphate Citrate Buffer in water, pH 6.0

Stop solution

- 10 50µL/well, 1M H₂SO₄
read at 450nm