

🌐 Evaluating endocytic rate in cells.

<https://dx.doi.org/10.17504/protocols.io.8epv5jjedl1b/v1>

¹KU Leuven

Marine Houdou

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.8epv5jjedl1b/v1>

External link: <https://www.thermofisher.com/order/catalog/product/T23366?>

gclid=CjwKCAjwzY2bBhB6EiwAPpUpZkuhTJcss8FLP963Hi9idgxNuMgwOubQVDObHI_Ru5cHPICunDbHdxoC24IQAvD_BwE
&ef_id=CjwKCAjwzY2bBhB6EiwAPpUpZkuhTJcss8FLP963Hi9idgxNuMgwOubQVDObHI_Ru5cHPICunDbHdxoC24IQAvD_Bw
E:G:s&s_kwcid=AL!3652!3!447292198772!!!g!!!10506731179!109642167771&cid=bid_pca_iva_r01_co_cp1359_pjt0000_bid0
0000_0se_gaw_dy_pur_con



Protocol Citation: Marine Houdou, Peter Vangheluwe, Nathalie Jacobs 2022. Evaluating endocytic rate in cells.. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.8epv5jjedl1b/v1>

Manuscript citation:

Houdou M, Jacobs N, Coene J, Azfar M, Vanhoutte R, Haute CVd, Eggermont J, Daniëls V, Verhelst SHL, Vangheluwe P, Novel Green Fluorescent Polyamines to Analyze ATP13A2 and ATP13A3 Activity in the Mammalian Polyamine Transport System. Biomolecules 13(2). doi: [10.3390/biom13020337](https://doi.org/10.3390/biom13020337)

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 03, 2022

Last Modified: May 31, 2024

Protocol Integer ID: 72253

Keywords: ASAPCRN, endocytic rate in cell, evaluating endocytic rate, endocytic rate, flow cytometry, cytometry readout, using tagged transferrin, tagged transferrin, cell

Abstract

Assess endocytic rate in cells using tagged transferrin (Alexa647) and flow cytometry readout.

Guidelines

- Make sure to **treat** the cells in **medium without FBS**.



Materials

0.25% Trypsin-EDTA: Gibco, 25200056

Albumin Fraction V: Carl Roth, 8076.4

Alexa647-Transferrin: T23366, invitrogen

Dulbecco's Phosphate Buffered Saline modified without calcium chloride and magnesium chloride (DPBS (-/-)): Gibco, D8537

Endocytosis inhibitors

- Dynasore: D7693, Sigma
- Genistein: ab120112, Abcam
- Pitstop 2: SML1169, Sigma

TrypLE Express Enzyme: Gibco, 12604021

Versene Solution: Gibco, 15040

Before start

- Prepare cell culture medium without FBS and keep it at 37°C.
- Prepare Flow Cytometry **(FC)** buffer made of 1% Albumin Fraction V in DPBS (-/-).
- Prepare Eppendorf tubes and FC tubes by labeling them and keeping them on ice.



- 1 Seed cells in **12 wellplate** that they reach **70-80% confluency** the day of the experiment.

Note

This protocol is made for a 12 wellplate format. Adaptations need to be done to scale up or down.

- 2 The day of the experiment, remove cell culture medium and **pre-treat** the cells in **medium without FBS** in a final volume of  500 μL /well , for

30m

 00:30:00 at 37°C, 5% CO₂ . The different conditions to consider are:

- 2.1 **No** pre-treatment for:

- blank samples
- Alexa647-transferrin-only samples.

In this case, change cell culture medium to cell culture medium without FBS.

- 2.2 **Pre-treatment** with endocytosis inhibitors cocktail made of:

-  100 μM Dynasore ,
-  50 μM Genistein and
-  50 μM Pitstop2 .

- 2.3 After pre-treatment, keep the plate **on ice** (4°C) for  00:15:00 .

15m

- 3 Add **Alexa647-Transferrin** at  50 μg /mL in each well, **except** for the **blank samples**.

Note

Alexa647-Transferrin is added to each well containing 500 μL of either culture medium without FBS or pre-treated medium including endocytosis inhibitors in culture medium without FBS.

- 4 Incubate  00:20:00 4°C (on ice)

20m



- 5 Incubate  00:20:00 37°C, 5% CO₂ . 20m
- 6 Harvest cells and prepare samples for Flow Cytometry (FC).
- 6.1 Discard medium.
- 6.2 Wash with  500 µL /well Versene or DPBS (-/-).
- 6.3 Discard Versene or DPBS (-/-).
- 6.4 Add  200 µL /well 0.25 % Trypsin or TrypLE. Incubate  00:05:00 at RT . 5m
- 6.5 Add  500 µL /well 1 % BSA-DPBS (-/-), collect the cells and transfer in Eppendorf tubes.
- 6.6 Centrifuge  00:05:00 at 2500 rpm and 4°C. 5m
- 6.7 Discard supernatants and resuspend cell pellet in  250-500 µL µL **1 % Albumin Fraction V in DPBS (-/-)**, depending on the size of the cell pellet.
- Note**
- Samples can't be too concentrated or too diluted to be ran on a flow cytometer.**
- If samples are too concetrated, cells will form clumps, they won't be single cells anymore and this will damage the flow cytometer.
 - If samples are too diluted, acquisition at the flow cytometer will take very long time.
- 6.8 Filter cell suspension through Nylon filter into FC tubes.
- 7 Keep on ice until acquisition at the flow cytometer. Record 10,000 live events per sample.

