




Jan 15, 2024

## Evaluating GPNMB ACD mutants by Western Blotting and immunofluorescence.

DOI

[dx.doi.org/10.17504/protocols.io.4r3l22yz4l1y/v1](https://dx.doi.org/10.17504/protocols.io.4r3l22yz4l1y/v1)

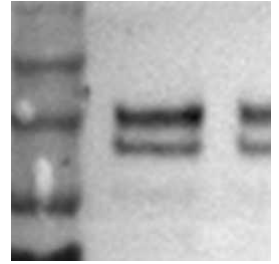

Erin Bogacki<sup>1</sup>, Patrick Lewis<sup>2,3</sup>, Susanne Herbst<sup>2,3</sup>

<sup>1</sup>NIH; <sup>2</sup>Royal Veterinary College;

<sup>3</sup>The Michael J. Fox Foundation for Parkinson's Research (MJFF) and the Aligning Science Across Parkinson's (ASAP) Initiative



Susanne Herbst

RVC

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l22yz4l1y/v1>

**Protocol Citation:** Erin Bogacki, Patrick Lewis, Susanne Herbst 2024. Evaluating GPNMB ACD mutants by Western Blotting and immunofluorescence.. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.4r3l22yz4l1y/v1>

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** November 13, 2023

**Last Modified:** May 31, 2024

**Protocol Integer ID:** 90843

**Keywords:** ASAPCRN, GPNMB, Western Blotting, Immunofluorescence, HEK293, gpnmb acd mutants by western blotting, gpnmb mutants by western blotting, evaluating gpnmb acd mutant, gpnmb mutant, immunofluorescence, immunofluorescent imaging, hek293 overexpression model, evaluation of cellular processing, cellular processing

**Funders Acknowledgements:**

The Michael J. Fox Foundation for Parkinson's Research (MJFF) and the Aligning Science Across Parkinson's (ASAP) Initiative  
Grant ID: .


## Abstract

This protocol describes the evaluation of cellular processing of GPNMB mutants by Western Blotting and Immunofluorescent imaging in a HEK293 overexpression model.



## Materials

### General

- HEK293T cells (ATCC CRL-3216)
- Eugene HD transfection reagent (E2311, Promega)
- pcDNA3.1-GPNMB-EGFP  pcDNA3.1-GPNMB-EGFP.png 104KB
- PBS, pH 7.4: #14190250, ThermoFisher Scientific

### Western Blotting

- Lysis buffer: 20 mM Tris-HCl (pH 7.5), 150 mM NaCl, 1 mM Na<sub>2</sub>EDTA, 1 mM EGTA, 1% (v/v) Triton X-100  
NOTE: add protease and phosphatase inhibitors fresh each time (eg Halt Protease and Phosphatase Inhibitor Cocktail (100X), #78440, ThermoFisher Scientific)
- Loading buffer: NuPAGE LDS sample buffer (#NP0007, ThermoFisher Scientific)
- Sample reducing agent: NuPAGE sample reducing agent (#NP0009, ThermoFisher Scientific)
- 4-12% Bis-Tris NuPAGE gels (eg #NP0321BOX, ThermoFisher Scientific)
- SDS-PAGE running buffer: MES running buffer (#NP0002, ThermoFisher Scientific)
- Trans-Blot® Turbo™ PVDF Transfer Packs: eg #1704157, BioRad
- TBS-T: 50 mM Tris-HCl, pH 7.5, 150 mM NaCl, 0.05% (v/v) Tween 20.
- Blocking and antibody dilution buffer: 5 % (w/v) non-fat milk powder in TBS-T
- Primary and secondary antibodies (see table 1 & 2 for antibody suggestions)


### Immunofluorescence

- 4 % (v/v) PFA/PBS: Dilute 16 % Paraformaldehyde Aqueous Solution (#15710, Electron Microscopy Sciences) to 4 % in PBS
- Blocking and antibody dilution buffer: 0.3 % Triton X-100, 5 % (v/v) FCS in PBS
- DAPI staining solution: 300 nM DAPI in PBS (#D1306, ThermoFisher Scientific or similar)
- Mounting medium: DAKO Fluorescence Mounting medium, # S3023, Agilent or similar
- Coverslips #1.5 (eg 631-0150, VWR)
- Slides (eg SuperFrost Plus, J1800AMNZ, Epredia)
- Primary and secondary antibodies (see table 3 & 4 for antibody suggestions)

## Troubleshooting



## Seed cells

- 1 Seed HEK293 cells.  
A) For Western Blotting, we recommend seeding  $2.5 \times 10^5$  cells per well of a 12-well culture plate.  
B) For Immunofluorescence, we recommend seeding  $1.2 \times 10^5$  cells per well of a 24-well culture plate. Seed cells on Poly-D-Lysine coated coverslips.  
Note: we routinely culture HEK293T cells in DMEM containing 10% FCS.  
Incubate in a tissue culture incubator  Overnight .






## HEK 293 cell transfection

10m

- 2 This protocol uses a DNA: Fugene ratio of 1:3. Prepare the transfection complexes as follows:

	A	B	C	D
		DNA	Fugene	Serum-free DMEM
12 well		1000 ng	3 ul	100 ul
24 well		500 ng	1.5 ul	50 ul

Preparation of transfection complexes (quantities are per well)

- 2.1 Add the required amount of plasmid DNA to serum-free DMEM. Mix briefly.
- 2.2 Add the required amount of Fugene HD Transfection Reagent.
- 2.3 Vortex and incubate for  00:10:00 at  Room temperature .
- 2.4 In the meantime, change the medium on the cells to fresh DMEM containing 10 % FCS
- 2.5 Add  50  $\mu$ L per 24-well or  100  $\mu$ L per 12-well drop-wise to the cells and incubate  Overnight

10m




## Western Blotting

25m

3 This section describes the sample preparation and analysis for Western Blotting.

3.1 Wash the cells gentle with PBS

3.2 Immediately add  100  $\mu$ L per well of ice-cold cell lysis buffer and place cells on ice.


3.3 Scrape cells with a cell scraper and harvest cell lysate into a 1.5 ml Eppendorf tube.



3.4 Incubate the cells  On ice for  00:10:00 , vortex occasionally

10m


3.5 Clear the cell lysate by spinning down at  16200 x g, 4°C for  00:15:00

15m

3.6 Transfer the post-nuclear supernatant to a fresh 1.5 ml Eppendorf tube and store at  -20 °C .

3.7 Prepare the cell lysates for Western Blotting by adding LDS sample buffer and denaturing agent and denature the samples at  80 °C for  00:08:00 .

8m


3.8 Run samples on 4-12 % Bis-Tris SDS-page. (approx.  00:35:00 at 160 V const.)

35m

3.9 Transfer proteins onto a PVDF membrane using the Turbo transfer system (BioRad) or similar.



3.10 Block membranes in 5% milk/TBS-T for  01:00:00 at  Room temperature .

1h



3.11 Incubate the membranes with primary antibodies at  4 °C  Overnight .

A	B	C	D	E
Target	Cat #	Supplier	Raised in	Dilution
GPNMB (N-terminal)	AF2550	R&D Systems	Goat	1:1000
GFP	MA5-15256	ThermoFisher Scientific	Mouse	1:1000
Actin	A1978	Sigma	Mouse	1:5000

Table 1: Primary antibodies for Western Blotting.

- 3.12 Wash the plots in TBS-T for  00:05:00 at  Room temperature . Repeat this step twice for a total of three washes.

5m

- 3.13 Dilute the secondary antibody in 5% milk/TBS-T and incubate the membranes with secondary antibodies at  Room temperature for  00:45:00 .

45m


A	B	C
Antibody suggestion		Dilution
anti-mouse-Peroxidase	eg, A3682, Sigma	1:10000
anti-goat-Peroxidase	eg, A5420, Sigma	1:10000

Table 2: Secondary antibodies for Western Blotting.



- 3.14 Develop the blots using an appropriate developer. Full-length GPNMB-EGFP is detected as a double band at ~125 kDa. A cleaved C-terminal GPNMB fragment can be detected with the anti-GFP antibody at ~35 kDa.

## Immunofluorescence

2h 20m

- 4 Gently wash coverslips with PBS and fix in 4 % PFA/PBS for  00:15:00 min.

15m

- 4.1 Gently wash the cells with PBS. Replace the PBS and add the permeabilisation/blocking. Incubate coverslips for a minimum of  00:20:00 min at  Room temperature .


20m

4.2 In the meantime, place a piece of Parafilm onto your bench and label if required. This will act as a flat surface to stain the coverslips on.

4.3 Prepare the antibody staining solution in Blocking and staining buffer. Find a suggestion of antibodies for counterstaining below:

	A	B	C	D	E
	<b>Target</b>	<b>Cat #</b>	<b>Supplier</b>	<b>Raised in</b>	<b>Dilution</b>
	LAMP-1	H4A3	DSHB	Mouse	1:100
	TGN46	13573-1-AP	Proteintech	Rabbit	1:100

Table 3: Primary antibodies for immunofluorescence.

4.4 Pipette a  45 µL drop of the antibody staining solution onto the Parafilm and invert the coverslip onto the staining solution so that the cells face downwards.

4.5 Incubate for  01:00:00 hr in the dark.


1h

4.6 Wash the coverslips three times with PBS.

4.7 Prepare a staining solution containing the secondary antibody:

A	B
<b>Antibody suggestion</b>	<b>Dilution</b>
anti-mouse-AF647	1:1000
anti-rabbit-AF586	1:1000

Table 4: Secondary antibodies for Immunofluorescence.

4.8 Pipette a  45 µL drop of the antibody staining solution onto the Parafilm and invert the coverslip onto the staining solution so that the cells face downwards.

4.9 Incubate for  00:45:00 min in the dark.

45m



- 4.10 Wash the coverslips once in PBS, stain with DAPI (or other nuclear stain), and mount onto glass slides.
- 4.11 GPNMB can be observed predominantly at the trans-Golgi network but can also be seen at lysosomal compartments.