



Oct 04, 2023

## Ex vivo mouse brain patch clamp recordings and Fura-2 imaging

DOI

[dx.doi.org/10.17504/protocols.io.eq2lyj8zrlx9/v1](https://dx.doi.org/10.17504/protocols.io.eq2lyj8zrlx9/v1)

taylor.panczyk<sup>1</sup>

<sup>1</sup>Northwestern University



taylor.panczyk

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.eq2lyj8zrlx9/v1>

**Protocol Citation:** taylor.panczyk 2023. Ex vivo mouse brain patch clamp recordings and Fura-2 imaging. **protocols.io**  
<https://dx.doi.org/10.17504/protocols.io.eq2lyj8zrlx9/v1>

**License:** This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** October 03, 2023



**Last Modified:** May 31, 2024

**Protocol Integer ID:** 88723

**Keywords:** ASAPCRN, mouse brain patch, whole cell patch clamp, mouse brain, electrophysiology, brain slice, attached configuration recording, recording, mouse, quick pulses of negative pressure, imaging in this protocol, quick pulse, recording chamber, types of recording

## Abstract

In this protocol we detail the steps to perform ex-vivo brain slices electrophysiology and Fura-2 recordings. For both types of recordings, the brain slices with the desired nucleus must be placed into the recording chamber and patch clamp high resistance seal must be performed. For cell attached configuration recording, the seal must be kept unbroken. For whole cell configuration, quick pulses of negative pressure to break into the cell. For Fura-2 recordings, a whole cell patch clamp must be performed to allow the Fura-2 to enter the cell.

## Materials

### Solutions to prepare:

#### Internal solutions:

*These are prepared prior to the experiment day and aliquoted in 1.5 ml tubes, and stored at -20°C until the day of experiment.*

#### K-based internal solution for whole cell

Compound	M.W.	Conc (mM)
KMeSO <sub>4</sub>	150.20	126.00
EGTA	380.40	1.00
KCl	74.55	14.00
HEPES	238.30	10.00
Mg-ATP	551.10	2.00
Na <sub>3</sub> -GTP	523.20	0.50
CaCl <sub>2</sub>	110.98	0.50
MgCl <sub>2</sub>	203.32	3

Adjust pH to 7.3 with KOH

#### Internal solution for Fura 2

Compound	M.W.	Conc (mM)
KMeSO <sub>4</sub>	150.20	135.00
EGTA	380.40	5.00
KCl	74.55	5.00
Na-PCr	255.10	10.00
HEPES	238.30	5.00
Mg-ATP	551.10	2.00
Na <sub>3</sub> -GTP	523.20	0.50

	CaCl <sub>2</sub>	110.98	0.50

Adjust pH to 7.3 with KOH

### Electrophysiology solutions:

*These are prepared as 10' stock solutions (below) and diluted for use on the day of the experiment.*

#### Stock aCSF

Solution 10x

1L

Compound	MW	g	conc (mM)
NaCl	58.44	79.33	1357.5
KCl	74.55	1.86	25
sodium bicarbonate	84.01	21	250
sodium phosphate	120	1.5	12.5

### Physiological glucose

aCSF 1x

1L

A	B	C	D
Compound	MW	g	conc (mM)
NaCl	diluted from 10x		125
KCl	diluted from 10x		2.5



	A	B	C	D
	<b>sodium bicarbonate</b>	<b>diluted from 10x</b>		<b>25</b>
	<b>sodium phosphate</b>	<b>diluted from 10x</b>		<b>1.25</b>
	<b>CaCl<sub>2</sub></b>	<b>1M solution</b>	2ml	<b>2</b>
	<b>MgCl<sub>2</sub></b>	<b>1M solution</b>	1ml	<b>1</b>
	<b>Glucose*</b>		0.63	<b>3.5</b>

\* add glucose to working solution (do not include in 10x stock)

### High glucose aCSF

Solution 1x

1L

	A	B	C	D
	<b>Compound</b>	<b>MW</b>	<b>g</b>	<b>conc (mM)</b>
	<b>NaCl</b>	<b>diluted from 10x</b>		<b>135.75</b>
	<b>KCl</b>	<b>diluted from 10x</b>		<b>2.5</b>
	<b>sodium bicarbonate</b>	<b>diluted from 10x</b>		<b>25</b>
	<b>sodium phosphate</b>	<b>diluted from 10x</b>		<b>1.25</b>
	<b>CaCl<sub>2</sub></b>	<b>1M solution</b>	2ml	<b>2</b>
	<b>MgCl<sub>2</sub></b>	<b>1M solution</b>	1ml	<b>1</b>
	<b>Glucose*</b>		4.5	<b>25</b>

\* add glucose to working solution (do not include in 10x stock)


## Troubleshooting



## Prepare patch pipettes

- 1 Turn on the Sutter P-1000 puller and enter the desired pull protocol.
- 2 Insert a thick-walled borosilicate glass capillary and press pull.
- 3 Pipette resistance must be of 2.5 to 5 megaohms.

## Setting up patch rig and environment

- 4 Turn on the MultiClamp 700B Amplifier, Axon Digidata 1550B digitizer, micromanipulator, computer tower and the associated software. Note: amplifier and digitizer must be turned on prior to opening software.
- 5 Turn on O<sub>2</sub>/CO<sub>2</sub> tank and bubble aCSF solution.
- 6 Take an aliquot of internal solution from the -20° fridge.
- 7 Fill syringe with internal solution (or aCSF for cell attached recordings), place a filter on the end of the syringe, and place a MicroFil Pipette Filler on the end of the filter.
- 8 Turn on the pump and circulate recording aCSF solution through chamber.
- 9 Adjust and set the rate of the pump to 3-4 mL/min.
- 10 Turn on water heater and set to desired temperature (~  34 °C )

## Examine slices and patching cells

- 11 Transfer brain slice from incubation beaker to the recording chamber.
- 12 Secure down slice with a harp (slice anchor).
- 13 Locate and focus the desired brain region under the 4x objective.
- 14 Change the microscope lens to the 60x objective.
- 15 Slowly focus on healthy neurons in slices for patching.
- 16 For whole cell configuration, fill a glass micropipette one-third full of internal solution. For cell attached configuration fill it with aCSF. Ensure there is no residual internal solution on exterior of glass micropipette, as this may introduce salts into the micromanipulator and add additional noise to recordings. Remove any air bubbles by gently flicking the glass micropipette.
- 17 Gently place the glass micropipette onto the wire electrode and tighten.
- 18 Position the electrode using a micromanipulator.
- 19 Under the 60x objective, bring the tip of the glass pipette above the slice.
- 20 Apply a positive pressure and maintain it.
- 21 Approach the cell diagonally. The positive pressure should create a small dimple on the cell.
- 22 Once a dimple is formed, zero the pipette voltage, release the positive pressure, and apply a small amount of negative pressure. The resistance should begin to increase rapidly.





- 23 As the resistance increases, clamp the cell at your resting potential of interest (typically -60 mV).
- 24 After a giga-ohm seal is formed, leave it this way to record in cell attached configuration or apply a few quick pulses of negative pressure to break into the cell to record in whole cell configuration.
- 25 Start recording.

## Fura 2 imaging

- 26 Transfer slices to a 2PLSM and turn on the 780-nm wavelength laser.
- 27 Add Fura-2 (100  $\mu$ M) and Alexa Fluor 594 hydrazide (25  $\mu$ M) to the internal solution.
- 28 Fill the patch pipettes with the internal solution.
- 29 Perform a whole cell patch clamp recording as previously explained.
- 30 Wait 10 minutes until the cell is completely filled with the internal solution containing Fura-2.
- 31 Choose a dendrite from the patched neuron 20-100  $\mu$ m far from soma.
- 32 Draw a straight line ROI over the chosen dendrite.
- 33 Start line scan acquisitions with 0.195- $\mu$ m pixels and 12-s dwell time.