



Mar 13, 2025

Expansion-assisted iterative fluorescence *in situ* hybridization (EASI-FISH) for morphoFISH

 Forked from [Expansion-assisted iterative fluorescence in situ hybridization \(EASI-FISH\) in *Drosophila* CNS](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.5jyl824n6l2w/v1>

Monique Copeland¹, Tiago Ferreira², Mark Eddison³


¹Scientific Operations, Janelia Research Campus; ²Janelia Research Campus;

³Project Technical Resources, Janelia Research Campus



Mark Eddison

Janelia Research Campus

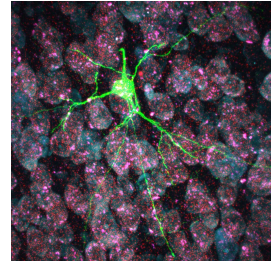

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.5jyl824n6l2w/v1>

Protocol Citation: Monique Copeland, Tiago Ferreira, Mark Eddison 2025. Expansion-assisted iterative fluorescence in situ hybridization (EASI-FISH) for morphoFISH. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.5jyl824n6l2w/v1>

Manuscript citation:

Discovery of neuronal cell types by pairing whole cell reconstructions with RNA expression profiles

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 09, 2024

Last Modified: March 13, 2025

Protocol Integer ID: 109476

Keywords: In situ hybridization, HCR, Expansion Microscopy, EASI-FISH, MorphoFISH, Mice, iterative fluorescence in situ hybridization, expansion microscopy, thick slices of the mouse brain, assisted iterative fluorescence, multiplexed in situ hybridization, situ hybridization, mouse slices with gfp, mouse brain, cleared mouse slice, labeled neuron, microscopy, hybridization chain reaction, modified gel recipe, gel recipe, neuron

Funders Acknowledgements:

HHMI

Grant ID: N/A

Abstract

A straightforward, robust, and reliable protocol (EASI-FISH) that utilizes expansion microscopy and the hybridization chain reaction for multiplexed *in situ* hybridization in thick slices of the mouse brain has recently been described (Wang et al., 2021). Below details a modified version of this protocol for cleared mouse slices with GFP-labeled neurons fixed under morphoFISH perfusion (Ferreira et al., 2024). This uses a modified gel recipe (TReX-1000) for greater gel stability. We also include an updated design for the Z1 sample holder.

Attachments



[Table of materials_M...](#)

10KB



[LIGHTSHEETHOLDER_V1](#)

8...

14.2MB

Image Attribution

Mark Eddison

Materials

Solutions for Day 1: Labelling RNA

-  MOPS (Fine White Crystals/Molecular Biology) **Fisher Scientific Catalog #bp308100** . Store  Room temperature .
- Melphalan-X (Stock  2 mg/mL ; working  1 mg/mL). Store  -20 °C .
-  Acryloyl-X, SE **Thermo Fisher Scientific Catalog #A20770** , Stock  10 mg/mL ; working  0.1 mg/mL . Store  -20 °C .
-  Press-to-Seal[®]; Silicone Isolator with Adhesive, eight wells, 9 mm diameter, 0.5 mm deep **Thermo Fisher Catalog #P24743**
-  Poly-L-Lysine solution 10ml **TEDPELLA Catalog #18026** ,  1.6 mL +  3.2 µL
 PHOTO-FLO 200 SOLUTION **Electron Microscopy Sciences Catalog #74257** . Store  4 °C .
- PBS-Triton (0.1%)
- 0.2 ml PCR tubes (USA Scientific; 1402-4700).
- RNase Away (Thermo Scientific; 7003).

Solutions for Day 2: Gelation and Proteinase K digestion

- Trex-1000. Store  -20 °C .
- 10%  Ammonium Persulfate **Merck MilliporeSigma (Sigma-Aldrich) Catalog #A3678** . Store  -20 °C .
- 10%  NNN'N'-Tetramethylethylenediamine **Merck MilliporeSigma (Sigma-Aldrich) Catalog #T22500** . Store  -20 °C .
- 0.5%  4-Hydroxy-TEMPO **Merck MilliporeSigma (Sigma-Aldrich) Catalog #176141**). Store  -20 °C .
-  Proteinase K, Molecular Biology Grade - 2 ml **New England Biolabs Catalog #P8107S** ,  800 U/ml). Store  -20 °C .
-  50 millimolar (mM) Proteinase K Buffer: Store  Room temperature .

- Small paint brush.

Solutions for Day 3: Hybridization

- Hybridization Buffer (Molecular Instruments, Store $-20\text{ }^{\circ}\text{C}$).
- Probe Wash Buffer (Molecular Instruments, Store $-20\text{ }^{\circ}\text{C}$).
- DNA oligo probes (Designed and made by Molecular Instruments, Store $-20\text{ }^{\circ}\text{C}$).
- DAPI (Sigma D9534)/  PBS Fisher Scientific Catalog #BP24384 at [M] 500 ng/ml .

Solutions for Day 5: Hybridization Chain Reaction (HCR)

- Amplification Buffer (Molecular Instruments, Store $4\text{ }^{\circ}\text{C}$).
- Fluorescent Hairpins (448 or 546 from Molecular Instruments or 669 conjugated in lab, Store $-20\text{ }^{\circ}\text{C}$).
- 5x SSCT (5x SSC, 0.1% Tween in Nuclease Free Water). Store Room temperature .
- 0.5x SSCT (0.5x SSC, 0.1% Tween in Nuclease Free Water). Store Room temperature .
-  Invitrogen GFP Polyclonal Antibody Alexa Fluor™ 488 Thermo Fisher Scientific Catalog #A-21311 .
-  UltraPure™ BSA (50 mg/mL) Thermo Fisher Scientific Catalog # AM2616 .

Solutions for Stripping Probes and Hairpins for Multiplexing

 DNase I, RNase-free Qiagen Catalog #79254 .

Note

Do **not** use RDD buffer supplied.

DNase1 buffer

	A	B
	Tris-HCL pH 8.0	10 mM
	MgCl2	2.5 mM

	A	B
	CaCl ₂	0.5 mM

Recipes and Reagents

200mM MOPS Buffer (10X Stock):

🧪 1046.5 mg in 🧪 25 mL in NFW, pH to 7.7 with 1M 10 Normality (N) NaOH. Store 🌡️ -20 °C .

Melphalan stock (2.5mg/ml)

🔗 Melphalan **Cayman Chemical Company Catalog #16665**

🔗 DMSO, Anhydrous **Thermo Fisher Catalog #D12345**

- Dissolve 🧪 2.5 mg per ml in anhydrous DMSO (Invitrogen, D12345).
- Dissolve, heat to 🌡️ 37 °C and vortex vigorously and place on a shaker.
- May take an hour to dissolve.
- Aliquot in 🧪 800 µL batches.
- Store in a desiccated environment at 🌡️ -20 °C .

Acryloyl-X (AcX) stock (10 mg/ml)

🔗 Acryloyl-X, SE **Thermo Fisher Scientific Catalog #A20770**

- Dissolve in anhydrous DMSO.
- Aliquot in 🧪 200 µL batches.
- Any extra aliquot in 🧪 5 µL batches (for extra AcX).
- Store in a desiccated environment at 🌡️ -20 °C .
- Don't re-use AcX after thawing.

Melphalan-X (2mg/ml):



- Combine an equal concentration of Acryloyl-X ($[M] 10 \text{ mg/mL}$) and Melphalan ($[M] 2.5 \text{ mg/mL}$) (1-part AcX to 4-parts Melphalan (ie. $\text{200 } \mu\text{L} : \text{800 } \mu\text{L}$).
- Incubate Overnight at Room temperature with shaking.
- Store in $\text{50 } \mu\text{L}$ aliquots in a desiccated environment at -20°C .
- Use at $[M] 1 \text{ mg/mL}$ by 1:1 dilution with $[M] 20 \text{ millimolar (mM)}$ MOPS.

Trex-1000:

$[M] 4.04 \text{ Mass Percent}$ Sodium Acrylate* – made from Acrylic Acid as it is made at variable purity

40% Acrylamide Solution **Bio-Rad Laboratories Catalog #1610140**

2% Bis Solution **Bio-Rad Laboratories Catalog #1610142**

Nuclease-free water **Ambion Catalog #AM9932**

10X PBS (ThermoFisher; AM9625)

4.04M Sodium Acrylate stock solution

- In a fume hood, place 5.5 mL of acrylic acid into a 50mL tube. Place tube in a Room temperature water bath (e.g. a beaker).
- Add 4.5 mL water.
- Add 7.2 mL $[M] 10 \text{ Mass Percent}$ NaOH gradually to prevent excessive heating and evaporation/boiling.
- Remove tube from hood (at this point most of the acrylic acid has been converted to non-volatile sodium acrylate).
- Add $[M] 1 \text{ Mass Percent}$ NaOH (nominally 1 mL) gradually until the pH is between 7.5 and 8 using a pH meter, at Room temperature . Do not use pH test strips.
- Add water up to a final volume of 20 mL .

Note

Acrylic acid has a pKa of 4.76 – at $\text{pH } 7.75$ this solution has about $[M] 4 \text{ millimolar (mM)}$ remaining buffering capacity.

For 9.4 mls Trex-1000 (not including APS +TEMED + 4HT):



	A	B
	4.04M Sodium Acrylate	2500 ul
	40% Acrylamide	3571 ul
	2% N,N MethylBisacrylamide	500 ul
	10x PBS	1000 ul
	Nuclease Free Water	1829 ul

10% APS:

	A	B
	APS	100 mg
	H2O	900 ul

10% Temed:

	A	B
	TEMED	100 ul
	H2O	900 ul

0.5% 4HT:

	A	B
	4HT	5 mg
	H2O	995 ul

Note

Store APS, TEMED and 4HT in  150 µL aliquots in PCR tubes @  -20 °C .

50mM ProK/SDS Buffer (50ml)

⊗ EDTA (0.5 M, pH 8.0, nuclease-free) Thermo Fisher Scientific Catalog #AM9260G

⊗ 10% SDS solution Thermo Fisher Scientific Catalog #15553027

	A	B
	Tris-HCL- pH8	50 mM
	EDTA	1 mM
	TritonX	0.5 %
	NaCl	50 mM
	SDS	0.3 %

RNase-Free DNase1

- Add  550 µL DNAase1 Buffer to DNase1 powder. Mix.
- Aliquot  50 µL per PCR tube, store @  -20 °C .

DNase1 buffer (50ml)

	A	B
	1 M Tris-HCL > 10 mM	500 ul
	1 M MgCl2 > 2.5 mM	125 ul
	1 M CaCl2 > 0.5 mM	25 ul
	Nuclease Free Water	49.350 ml

HCR Wash Buffers (50 ml)

5x SSCT:



⊗ SSC, RNase-free, 20× Ambion Catalog #AM9763

	A	B
	20x RNase free-SSC (ThermoFisher, AM 9763)	12.5 ml
	10% Tween	500 ul
	Nuclease Free Water	37 ml

0.5x SSCT:

	A	B
	20x RNase free-SSC	1.25 ml
	10% Tween	500 ul
	Nuclease Free Water	48.25 ml



Preparation of slices for EASI-FISH

- 1 To visualize the GFP neurons, rehydrate the slice in PBS for 10 mins.
- 2 Under a dissection scope and using a sharp knife, cut ROI with GFP cells from brain slice. The area should be about 2mm x 4mm and have a nick in the right hand corner to keep track of orientation. Take a photo of before and after to keep track of slice and cells.
- 3 Store ROI slice in 70% EtOH until ready to use

Day 1: Labelling RNA

- 4 Transfer each ROI slice to a 0.2ml PCR tube
- 5 Rehydrate with 2 × 5 min wash in  150 µL PBT (0.1%). 
- 5.1 Rehydrate with 2 x  00:05:00 wash in  150 µL PBT (0.1%) (1/2). 
- 5.2 Rehydrate with 2 x  00:05:00 wash in  150 µL PBT (0.1%) (2/2). 
- 6 Incubate slice 1 x  00:30:00 in  150 µL  20 millimolar (mM) MOPS Buffer. 

- 7 Thaw Melphalan-X (MelphX) and Acryloyl-X (Ac-X) solutions.
- 8 Using a P20 pipette, take off as much MOPS buffer from the brains as possible. 
- 9 Dilute Melphalan-X stock 1:1 with MOPS Buffer (to 1mg/ml).



10 Add 1/100 Ac-X (**10 mg/mL**) to Melphalan-X working solution. Vortex, mix, and spin.



11 Add **50 µL** of Melphalan-X/AcX solution to each PCR tube and gently mix.



12 Incubate **Overnight** @ **37 °C** .

2m



13 Prepare gel chambers for gelation the next day. Wipe a non-charged slide with RNase away. Adhere a gasket (4-6 wells max) and coat the glass surface of the chamber with **1 µL** poly-Lysine using a P20 pipette tip. Air dry and repeat.



Chambers.jpg

Day 2: Gelation and Proteinase K digestion

14 Wash slice 2 × 2 min **150 µL** PBT and 1 × 2 mins **150 µL** PBS.



14.1 Wash slice 2 x **00:02:00** **150 µL** PBT and 1 x **00:02:00** **150 µL** PBS.

4m

14.2 Wash slice 2 x **00:02:00** **150 µL** PBT and 1 x **00:02:00** **150 µL** PBS.

4m

15 Thaw TREx-1000 and 4HT, Temed and APS. Vortex well and keep **On ice** .



16 Gently stick down the slice in the center of the chamber, once stuck down, carefully add a drop of PBS to prevent dehydration.



17 Mix together Trex-1000 and 4HT, Temed and APS at a ratio of 94:2:2:2. Vortex.



**Note**

Each chamber needs ~  120 μL of Trex-1000, make excess. (ie. for two chambers make  300 μL gel solution).

18 Remove PBS from chamber with pipette tip and carefully wick away remaining PBS with a tissue.

19 Pipette  40 μL of gel solution, on top of the slice, to each chamber. Incubate slide in the fridge ( 4 $^{\circ}\text{C}$) for  00:10:00 .

10m



20 Take off gel solution and repeat step 22.

Note

Place waste gel solution in an Eppendorf. Before discarding, polymerize the gel waste @  37 $^{\circ}\text{C}$

21 Take off the gel solution and gasket surface adhesive. Add a final  40 μL of gel solution and gently place a cover slip over the chamber. Gently press to seal the coverslip and incubate @  4 $^{\circ}\text{C}$ for  00:10:00 .

10m

**Note**

Adding  5 μL of gel solution to the center of the underside of the coverslip can help prevent air bubbles when sealing.

22 Polymerize the gel @  37 $^{\circ}\text{C}$ for  01:30:00 to  02:00:00 .

3h 30m

23 Cool gels on the bench for a few minutes.



24 Take off the chamber lid and gasket with a razor blade. Trim the gels into a rectangle and nick the top right-hand corner to track the orientation of the sample.

25 Take off the gel from the slide with a paintbrush that has been wetted with a small amount of ProK Buffer and transfer it to a  2 mL Eppendorf.

26 Incubate each gel with  1 mL ProK Buffer and  10 μ L (1/100) ProK Enzyme @  37 °C  Overnight .

2h



Day 3: Hybridization

27 Wash gels 4 $\times$ 15 min with  1 mL PBS at  Room temperature .



27.1 Wash gels 4 $\times$  00:15:00 in  1 mL PBS at  Room temperature (1/4).

15m

27.2 Wash gels 4 $\times$  00:15:00 in  1 mL PBS at  Room temperature (2/4).

15m

27.3 Wash gels 4 $\times$  00:15:00 in  1 mL PBS at  Room temperature (3/4).

15m

27.4 Wash gels 4 $\times$  00:15:00 in  1 mL PBS at  Room temperature (4/4).

15m

28 DNase treat sample before hybridization (see Step 48) for Cyto-DAPI staining/segmentation

29 Thaw and mix hybridization (hyb) and probe wash buffer.



30 Incubate gel in  500 μ L hyb buffer for  00:30:00 @  37 °C .

30m



31 Dilute probes 1/100 ([IM] 10 Mass Percent), in 300 µL hyb buffer per gel. Vortex.
Incubate @ 37 °C .



32 Incubate gels with probes overnight @ 37 °C , no shaking necessary.



33 Put probe wash buffer and PBS @ 37 °C .

Day 4: Probe Washing

34 Wash 3 × 30 min 750 µL Probe Wash Buffer @ 37 °C .



34.1 Wash 3 x 00:30:00 750 µL Probe Wash Buffer @ 37 °C (1/3).

30m

34.2 Wash 3 x 00:30:00 750 µL Probe Wash Buffer @ 37 °C (2/3).

30m

34.3 Wash 3 x 00:30:00 750 µL Probe Wash Buffer @ 37 °C (3/3).

30m

35 Wash 3 × 30 min 1 mL PBS @ 37 °C .



35.1 Wash 3 x 00:30:00 1 mL PBS @ 37 °C (1/3).

30m

35.2 Wash 3 x 00:30:00 1 mL PBS @ 37 °C (2/3).

30m

35.3 Wash 3 x 00:30:00 1 mL PBS @ 37 °C (3/3).

30m

36 Wash 3 × 1hr 1 mL PBS @ 37 °C .



36.1 Wash 3 x 01:00:00 1 mL PBS @ 37 °C (1/3).

1h



36.2 Wash 3 x 01:00:00 1 mL PBS @ 37 °C (2/3).

1h

36.3 Wash 3 x 01:00:00 1 mL PBS @ 37 °C (3/3).

1h

37 Keep gels in PBS at Room temperature Overnight .

1h



Day 5: Hybridization Chain Reaction (HCR)

38 Incubate gel in 500 μ L Amplification buffer for at least 00:30:00 @ Room temperature

30m



39 Snap cool hairpins with PCR machine @ 95 °C for 00:01:30 and cool @ Room temperature for 00:30:00 .

31m 30s



Note

Choice of far-red dyes: We found that JF669 tends to bleach faster in morphoFISH tissue. For this reason, we prefer to use AF647-conjugated hairpins.

40 Mix hairpins h1 and h2 @ 1/100 in 300 μ L Amp Buffer per gel. Vortex.



41 Incubate gel with hairpins for 03:00:00 @ Room temperature in the dark.

3h



42 Wash gels 2 $\times$ 20 min in 750 μ L 5X SSCT @ Room temperature .



42.1 Wash gels 2 x 00:20:00 in 750 μ L 5X SSCT @ Room temperature (1/2).

20m

- 42.2 Wash gels 2 x  00:20:00 in  750 μ L 5X SSCT @  Room temperature (2/2). 20m
- 43 Wash gels 2 $\times$ 40 min in  1 mL 0.5X SSCT @  Room temperature . 
- 43.1 Wash gels 2 x  00:40:00 in  1 mL 0.5X SSCT @  Room temperature (1/2). 40m
- 43.2 Wash gels 2 x  00:40:00 in  1 mL 0.5X SSCT @  Room temperature (2/2). 40m

Day 6: Mount and Image

- 44 Wash 2 $\times$ 30 min with  1 mL PBS-Triton (0.1%). 
- 44.1 Wash 2 x  00:30:00 with  1 mL PBS-Triton (0.1%) (1/2). 30m
- 44.2 Wash 2 x  00:30:00 with  1 mL PBS-Triton (0.1%) (2/2). 30m
- 45 Wash 2 $\times$ 30 min and 1 $\times$ 1hr with  1 mL PBS. 
- 45.1 Wash 2 x  00:30:00 with  1 mL PBS (1/2). 30m
- 45.2 Wash 1 x  01:00:00 with  1 mL PBS (2/2). 1h
- 46 DAPI stain gels for  01:00:00 with  1 mL PBS/DAPI ( 500 ng/ml). 1h
- 47 Mount gels (sample up) on an 8mm poly-lysine coated coverslip superglued to a Z.1 light-sheet sample holder and image. 



48 For long-term storage, keep gels in  750 μ L 5X SSC @  4 °C

Stripping Probes and Hairpins for Multiplexing

2h 45m

49 Incubate gel for  00:30:00 in  1 mL of DNase1 Buffer @  37 °C .



50 Add  450 μ L of DNase Buffer to  50 μ L DNase1. Mix.



51 Incubate gel in DNase1 for  02:00:00 @  37 °C .

2h



52 Wash 4 $\times$ 15 min with  1 mL PBS.



52.1 Wash 4 $\times$  00:15:00 with  1 mL PBS (1/4).

15m

52.2 Wash 4 $\times$  00:15:00 with  1 mL PBS (2/4).

15m

52.3 Wash 4 $\times$  00:15:00 with  1 mL PBS (3/4).

15m

52.4 Wash 4 $\times$  00:15:00 with  1 mL PBS (4/4).

15m

53 Hybridize with next round of probes (Day 3, step 34).



Protocol references

Wang et al. (2021) EASI-FISH for thick tissue defines lateral hypothalamus spatio-molecular organization. Cell 184 6361-6377.e24, <https://pubmed.ncbi.nlm.nih.gov/34875226>

Hugo GJ et al. (2022) Visualizing cellular and tissue ultrastructure using Ten-fold Robust Expansion Microscopy (TREx). eLife 11:e73775, <https://doi.org/10.7554/eLife.73775>

Ferreira AT et al. (2024) Discovery of neuronal cell types by pairing whole cell reconstructions with RNA expression profiles. bioRxiv 2024.12.30.630829, <https://doi.org/10.1101/2024.12.30.630829>