

Mar 01, 2019

Version 1

Experiment protocol: a syringe-filter based DNA extraction V.1

DOI

dx.doi.org/10.17504/protocols.io.yqzfvx6

Sungwoo Bae¹

¹National University of Singapore



Sungwoo Bae

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.yqzfvx6>

Protocol Citation: Sungwoo Bae 2019. Experiment protocol: a syringe-filter based DNA extraction. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.yqzfvx6>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: March 01, 2019



Last Modified: March 01, 2019

Protocol Integer ID: 20985

Keywords: dna extraction, based dna extraction, dna, syringe, experiment protocol, filter

Troubleshooting

- 1 Prepare the following solutions
 - a. TE buffer (10 mM Tris-HCl and 1 mM EDTA)
 - b. TE/lysozyme (TL) buffer (10 mM Tris-HCl, 1 mM EDTA, and 7.5 mg/mL lysozyme)
 - c. TE/proteinase K/SDS (TPS) buffer (10 mM Tris-HCl, 1 mM EDTA, 300 µg/mL proteinase K, and 1% SDS[w/v])
- 2 Filtration

The 10 mL sample or 100µl was passed through a 0.2 µm pore-sized syringe filter in the direction of the filtration.
- 3 TE buffer addition

A 1 mL of TE buffer was injected into the same filter, still in the direction of the filtration; the flow through was pooled into a sterile 15 mL tube.
- 4 Backwashing step

Backwashing was done in the opposite direction of filtration with the flow through pooled into the same 15 mL tube; the tubes were mixed by inversions for 5 mins in between each back-washing step:

 - a. A 1 mL of TL buffer was flushed through the filter then 300 µL of TL was added into the pooled flow-through.
 - b. Next, 1 mL of TPS buffer was flushed through the filter.
 - c. Lastly, 1 mL of TPS buffer and 3 times of 1 mL air was flushed and mixed to the flow through.
 - d. The filter is then discarded
- 5 Magnetic beads addition

A 50 µL of silica-coated magnetic beads solution (Dynabeads MyOne SILANE, Invitrogen) was added to the flow-through and was mixed by inversions for 5mins.
- 6 Washing

The beads were magnetized to the tube walls then the supernatant was discarded.
- 7 DNA elution and collection
 - a. The DNA was then eluted from the beads with 50 µL of TE buffer.
 - b. The magnetic beads were collected, and the DNA samples were recovered.
 - c. The recovered DNA was then diluted 3x with TE buffer to lessen possible inhibitions on downstream applications. The diluted DNA can also be stored in -20°C.