

Jun 30, 2023

Version 4

# Explant Surgery: Chronic recoverable Neuropixels in mice V.4

 In 1 collection

DOI

[dx.doi.org/10.17504/protocols.io.bp2l6113dvqe/v4](https://dx.doi.org/10.17504/protocols.io.bp2l6113dvqe/v4)


Emily A Aery Jones<sup>1</sup>

<sup>1</sup>Stanford University



Emily A Aery Jones

Stanford University

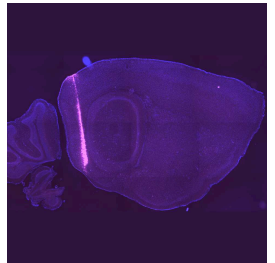

## Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bp2l6113dvqe/v4>

**Protocol Citation:** Emily A Aery Jones 2023. Explant Surgery: Chronic recoverable Neuropixels in mice. **protocols.io**  
<https://dx.doi.org/10.17504/protocols.io.bp2l6113dvqe/v4> Version created by **Emily A Aery Jones**

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



**Protocol status:** Working

**We use this protocol and it's working**

**Created:** June 30, 2023

**Last Modified:** June 30, 2023

**Protocol Integer ID:** 84255

**Keywords:** electrophysiology, silicon probe, electrode, Neuropixels, chronic recoverable neuropixels in mice, chronic recoverable neuropixel, implanted neuropixel probe, neuropixel probe, neuropixel, mice, explant surgery, probes for future use

**Funders Acknowledgements:**

Simons Collaboration on the Global Brain

Grant ID: 542987SPI

Stanford School of Medicine Dean's Postdoctoral Fellowship

A.P. Giannini Foundation Postdoctoral Fellowship

## Abstract

This protocol collection explains how to build a low-cost, lightweight system to implant 1 Neuropixels 1.0 probe or 2 Neuropixels 2.0 probes into mice, record during freely moving behavior, then recover the probes for future use. This protocol explains how to recover a previously implanted Neuropixel probe and clean for future use. See full collection for more details.

## Materials

*Tools:*

- Dovetail holder & stereotax adapter (Sensapex uMp-NH & uMp-NPR-200)
- Tiny flathead screwdriver (Wiha, 0.8 × 40)
- Scissors
- Headbar stereotactic adapters (see *Implant surgery* protocol)
- optional: ultrasonic cleaner (e.g. Amazon B075FTFD3R)

*Consumables:*

- Isoflurane
- Tergazyme (1g per 100mL, Alconox 1325, made within 8h of surgery)
- dH2O
- Dow-Sil DS-2025 Silicone Cleaning Solvent
- 100% isopropyl alcohol (IPA)

## Troubleshooting



## Prepare mouse

5m

- 1 Set O<sub>2</sub> flow rate to 1-1.5L/min and isoflurane to 3%. Place mouse in anesthetic chamber.

### Safety information

Use vacuum systems connected to isoflurane lines and work on a downdraft table to move uninhaled isoflurane away from the surgeon.

- 2 When breathing has slowed to 1Hz, move animal to the toothbar. Switch isoflurane from chamber to nosecone. Wait until unresponsive to pedal reflex test, then lower to 1.5% isoflurane.

### Note

Monitor and record anesthetic depth throughout procedure. Ears and toes should be pink, respiration should be once per 1-2 seconds, and pedal reflex should be lost. If respiration slows or extremities start to lose color, lower isoflurane levels. If respiration speeds up or reflex is present, increase isoflurane levels and halt the procedure until animal is fully anesthetized again.

- 3 Set the nosecone to the position you recorded from the implant surgery. Attach mouse to headbar stereotactic adapters.

## Remove probe

25m

- 4 Remove all the tape to expose the screws and probe.
- 5 Cut the ground wire above the pin.
- 6 Attach the Sensapex holder and set the stereotax to the angle used in the implant surgery. Tax to the dovetail cap. A fully open Sensapex holder can be taxed in from anterior or dorsal to the dovetail holder.
- 7 Adjust the screws so the holder grips the dovetail cap.

- 8 Unscrew the wings from the body. Screws can be saved for future surgeries.
- 9 Slowly tax up to bring the probe & body piece away from the animal. Set back to 0 degrees and adjust the arm so the probe is away from the surgical field.
- 10 Lower the probe into tergazyme and let sit for 30 minutes, then rinse with dH2O.

## Perfuse

20m

- 11 Remove the mouse from the stereotax and prepare for perfusion.



Remaining implant after single 1.0 probe is explanted. After perfusion, the assembly can be removed from the skull in 1 piece by grasping the headbar and rotating it away from the skull.

## Re-use probe

- 12 Probes may have residual Dow-Sil on them or dried gunk. Examine probe shanks under the stereoscope and look for changes in reflection of light along the shank. To remove Dow-Sil, soak shank in Dow-Sil DS-2025 Solvent for 10 minutes, then dip several times in 100% IPA to remove solvent.
- 13 If pieces still remain or 2.0 probe shanks are sticking together, dip into ultrasonic cleaner for 10 second intervals in 100% IPA, examining between each dip and stopping once shank is clean or at least improved. This will quickly clear small residual pieces and will dry out larger pieces so they can be removed manually.
- 14 If the previous step still wasn't sufficient, move to manual cleaning and shank separation using the 30g needle. Mount the probe on the stereotax, shanks facing down and PCB facing you as if you were about to insert the probe during a surgery. Visualize through the stereoscope, focused so you can see the whole shank. Brace your hand against a



solid surface and your second hand to steady it. Hold the syringe parallel to the table and bring the needle perpendicular to the shanks. Insert the bent tip between the probe shanks high on the probe and gently move downwards until they separate. The shanks are designed to safely flex towards and away from you along this axis, but can break if flexed left or right, so move the needle up and down only, following the shanks. It's okay if the needle catches on the shanks and pulls them toward or away from you; keep moving slowly down or up, never sideways, until it releases.

- 15 Assemble following the *Assembly* protocol, skipping the steps that modify the probe (sharpening, soldering wire to the PCB, attaching the body piece). Extend the ground wire back to full length and attach a new pin.