



Dec 02, 2022

Expression and purification protocol of the Human (Homo sapiens) LC3B Ubiquitin-like modifier

 Forked from [Expression and purification protocol of the Human \(Homo sapiens\) mCherry-LC3B Ubiquitin-like modifier](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.j8nlkw82dl5r/v1>

Dorotea Fracchiolla¹, Liv Jensen²

¹Team Hurley; ²University of California, Berkeley



Liv Jensen



Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.j8nlkw82dl5r/v1>

Protocol Citation: Dorotea Fracchiolla, Liv Jensen 2022. Expression and purification protocol of the Human (Homo sapiens) LC3B Ubiquitin-like modifier. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.j8nlkw82dl5r/v1>

Manuscript citation:

<https://doi.org/10.1083/jcb.201912098>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: December 02, 2022

Last Modified: May 31, 2024

Protocol Integer ID: 73497

Keywords: expression, purification, recombinant, Microtubule-associated proteins 1A/1B light chain 3B, LC3B, mCherry-tag, ASAPCRN, tagged human recombinant ubiquitin, human recombinant ubiquitin, lc3b ubiquitin, purification protocol of the human, protein conjugation, associated protein, protein, like modifier lc3b, purification protocol, purification, purification procedure, amino acid, phosphatidyl ethanolamine, microtubule, like modifier this protocol, map1lc3b

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to [protocols.io](#) is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with [protocols.io](#), can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract


This protocol describes expression and purification procedures for obtaining mCherry-tagged human recombinant Ubiquitin-like modifier LC3B (MAP1LC3B, Microtubule-associated proteins 1A/1B light chain 3B) lacking five C-terminal amino acids to allow in vitro protein conjugation to target PE, Phosphatidyl Ethanolamine.

Image Attribution


© Dorotea Fracchiolla

Guidelines

Affinity purification

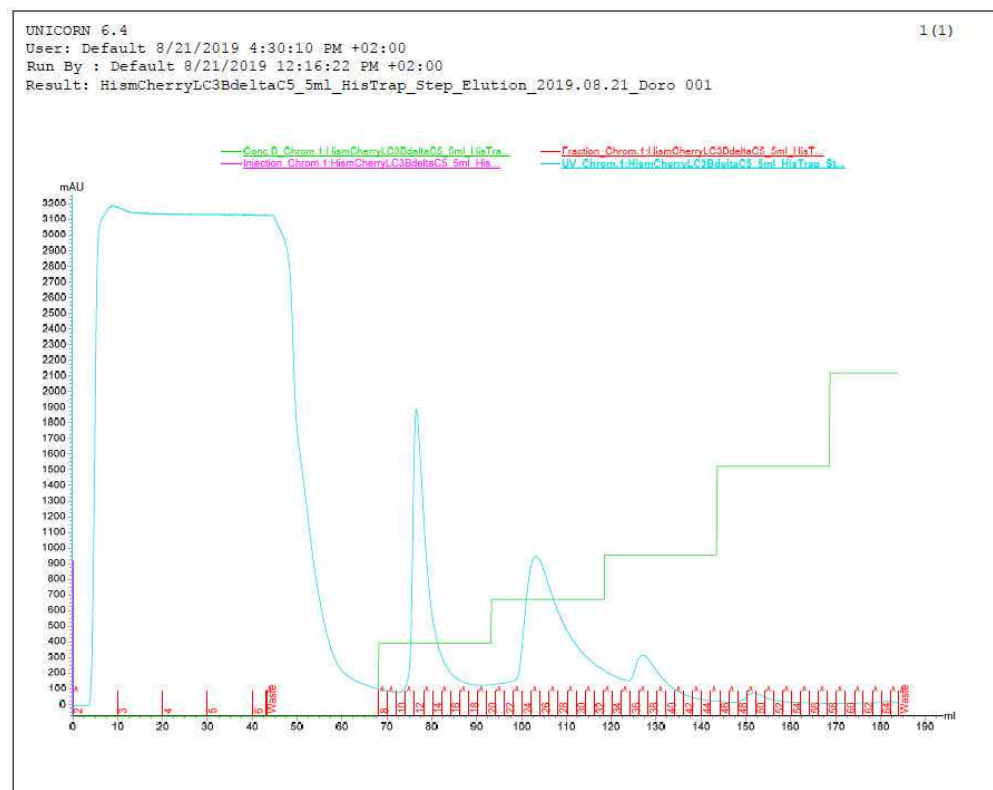


Chromatograph of His-tag affinity purification for mCherry-hLC3B.

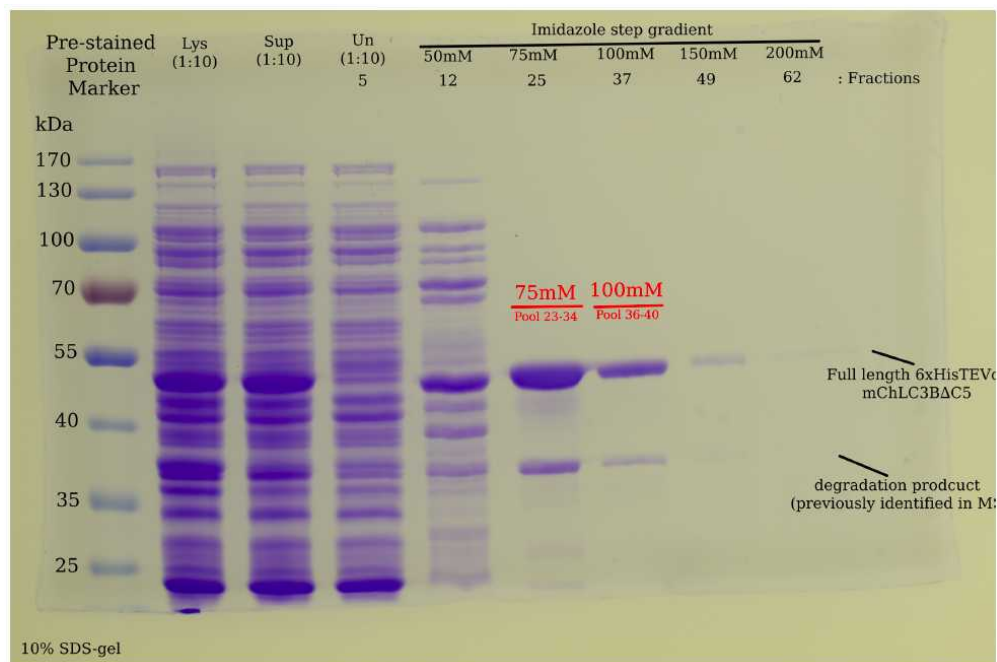


Coomassie BB stained gel of His-tag affinity purification for mCherry-hLC3B.

Size Exclusion Chromatography

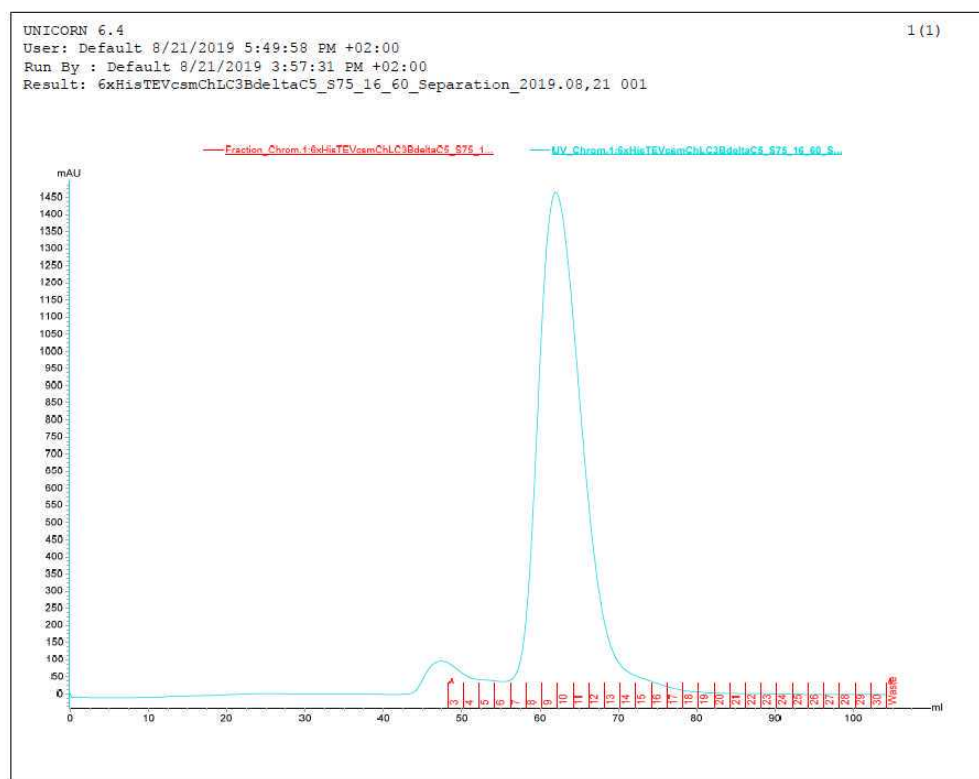

Chromatograph of Size Exclusion purification for mCherry-hLC3B.

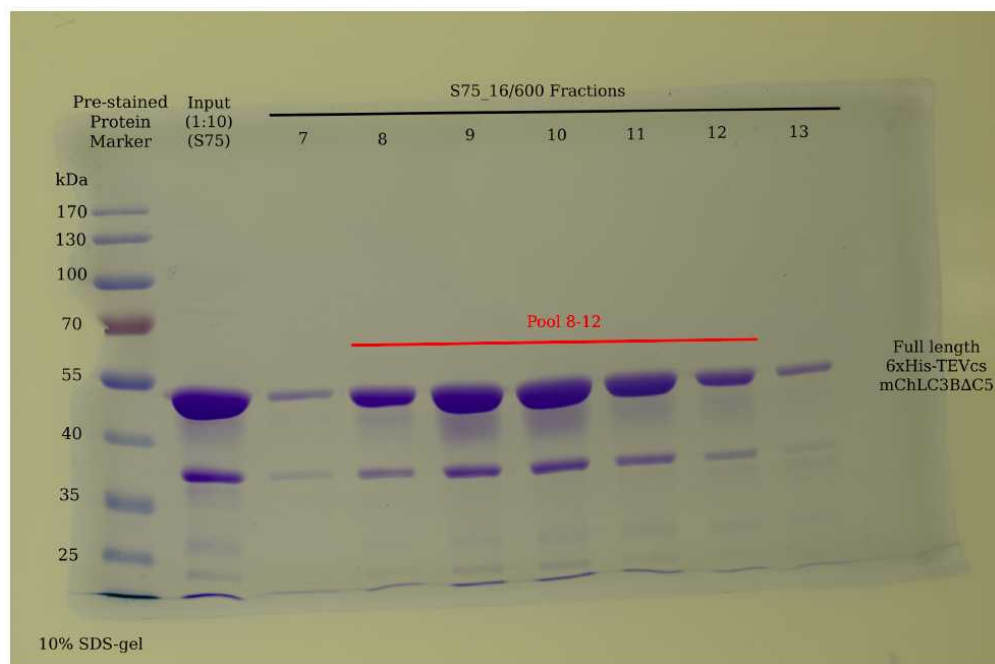

Coomassie BB stained gel of Size Exclusion Chromatography purification for mCherry-hLC3B.



Materials

Materials and Reagents

- Escherichia coli Rosetta pLyss cells
- Luria Bertani (LB) medium with antibiotics (final conc. 50µg/ml Ampicillin, 34µg/ml Chloramphenicol)
- IPTG (isopropyl-β-D-thiogalactopyranoside)
- 37°C shaker incubator
- sterile flasks/sterile pipettes
- tip sonicator
- **Lysis Buffer:** 50mM Hepes pH=7.5, 300mM NaCl, 10mM Imidazole, 2mM MgCl₂, 2mM β-mercaptoethanol, 1mM Pefablock, Complete Protease Inhibitors (EDTA-free CIP tablet, Roche), DNase (Sigma).
- **Buffer A:** 50mM Hepes pH=7.5, 300mM NaCl, 10mM Imidazole (filtered and degassed) + 1mM β-mercaptoethanol;
- **Buffer B:** 50mM Hepes pH=7.5, 300mM NaCl, 300mM Imidazole (filtered and degassed) + 1mM β-mercaptoethanol;
- **Size Exclusion Chromatography (SEC) Buffer:** 25mM Hepes pH=7.5, 150mM NaCl, 1mM DTT (Dithiothreitol).

Note: all purification buffers are filtered and degassed. Reducing agents (β-mercaptoethanol and Dithiothreitol) are added after degassing step.

Columns:- HT 5ml column (GE Healthcare)
- S75_16/60 (GE Healthcare)

Gels:10% SDS-PAGE

Safety warnings

! Please refer to the Safety Data Sheets (SDS) for health and environmental hazards.

Before start

General information

Insert: Homo sapiens LC3B, [NP_073729.1](#); Expression system: E.Coli Rosetta pLyss; plasmid origin: Sascha Martens Lab, Addgene 169168, lab internal construct database number SMC948; backbone: pET-Duet1; plasmid resistance: Ampicillin; tags & cleavage sites: N-term 6xHis, followed by Tobacco Etch Virus (TEV) cleavage site, mCherry tag, LC3B ORF. Ext coeff: 41830 M⁻¹ cm⁻¹, MW 47,89 kDa.



Protein Expression

16h 15m

- 1 Transform plasmid DNA (Addgene_190237) into E.Coli Rosetta pLyss cells and plate on Ampicillin LB agar plate for Overnight at 37 °C .
- 2 The following day, inoculate a 5 mL LB + Amp/Cam pre-culture with 1-2 colonies and grow Overnight at 37 °C shaking.
- 3 The following day, use 5 mL pre-culture to inoculate 1 L LB medium + Amp/Cam at 37 °C until an OD₆₀₀ (Optical Density) of 0.6 is reached.
- 4 Induce protein expression with 400 micromolar (μM) IPTG and grow for a further 16:00:00 at 18 °C shaking. 16h
- 5 Pellet cells at 4000 rpm, 4°C, 00:15:00 in a Sorvall RC6+ centrifuge (Thermo Scientific), discard supernatant and resuspend pellets in ice cold lysis buffer (25 ml/1 lt culture). 15m
- 6 Flash freeze resuspended pellets in liquid nitrogen and store at -80 °C until purification.

Protein Purification

- 7 Perform His-Trap affinity purification followed by Size Exclusion Chromatography.

Purification of LC3B Ubiquitin-like modifier

30m 30s

- 8 Cells are lysed via freeze/thaw cycles and sonication: thaw pellet corresponding to 1 L culture by freeze/thawing in Room temperature water bath. All following steps are to be executed at 4 °C or On ice .



- 9 Lyse cells by sonicating them with an immersion Tip Sonicator (2x  00:00:30). Note: adjust times and intensity according to the available instrument. 
- 10 Clear lysate by spinning in a Beckman centrifuge, at  40000 x g, 4°C, 00:30:00 , Ti45 Rotor . 
- 11 Inject soluble fraction onto a 5ml HT column operating at  4 °C pre-equilibrated in Buffer A at 1ml/min flow rate.
- 12 Wash column for  5 µL with Buffer A at 2 ml/min flow rate to remove unspecific bound proteins. 
- 13 Elute protein at 300mM Imidazole concentration. Perform elution at 1ml/min flow rate.
- 14 Perform TEV cleavage with 1mg/ml TEV protease overnight at 4°C while dialyzing against elution buffer containing no imidazole.
- 15 Apply cleavage reaction to equilibrated HT column, and collect the flow-through.
- 16 Measure protein absorbance at A₂₈₀ with a Spectrophotometer against dialysis buffer. 
- 17 Aliquot protein, flash freeze in liquid Nitrogen and store at  -80 °C .