

Jul 05, 2017

Extraction method C (FMS)

 [GigaScience](#)

 In 1 collection

DOI

[dx.doi.org/10.17504/protocols.io.inbcdan](https://doi.org/10.17504/protocols.io.inbcdan)

Faezah Mohd Salleh¹, Jazmin Ramos-Madriral¹, Fernando Penaloza¹, Shanlin Liu¹, Mikkel-Holger S Sinding¹, Riddhi P Patel¹, Renata Martins¹, Dorina Lenz¹, Jorns Fickel¹, Christian Roos¹, Mohd Shahir Shamsir¹, Mohammad Shahfiz Azman¹, Burton K Lim¹, Stephen J Rossiter¹, Andreas Wilting¹, M Thomas P Gilbert¹

¹Natural History Museum of Denmark, Copenhagen, Denmark

GigaScience Press



Faezah Mohd Salleh

OPEN  ACCESS



DOI: [dx.doi.org/10.17504/protocols.io.inbcdan](https://doi.org/10.17504/protocols.io.inbcdan)

External link: <https://doi.org/10.1093/gigascience/gix053>

Protocol Citation: Faezah Mohd Salleh, Jazmin Ramos-Madriral, Fernando Penaloza, Shanlin Liu, Mikkel-Holger S Sinding, Riddhi P Patel, Renata Martins, Dorina Lenz, Jorns Fickel, Christian Roos, Mohd Shahir Shamsir, Mohammad Shahfiz Azman, Burton K Lim, Stephen J Rossiter, Andreas Wilting, M Thomas P Gilbert 2017. Extraction method C (FMS). **protocols.io** <https://dx.doi.org/10.17504/protocols.io.inbcdan>

Manuscript citation:

Salleh FM, Ramos-Madriral J, Peñaloza F, Liu S, Mikkel-Holger SS, Riddhi PP, Martins R, Lenz D, Fickel J, Roos C, Shamsir MS, Azman MS, Burton KL, Stephen JR, Wilting A, Gilbert MTP, An expanded mammal mitogenome dataset from Southeast Asia. GigaScience 6(8). doi: [10.1093/gigascience/gix053](https://doi.org/10.1093/gigascience/gix053)

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

Created: June 26, 2017

Last Modified: November 09, 2017


Protocol Integer ID: 6563


Abstract

This protocol provides an efficient DNA extraction and purification of historical museum hides, which potentially have been chemically tanned.


Materials


MATERIALS

 MinElute PCR Purification Kit **Qiagen Catalog #28004**

 Chloroform

 Sodium Hypochlorite Solution

 70% Ethanol

 Digestion Buffer consisting of 10 mM tris(hydroxymethyl)aminomethane hydrochloride (Tris-HCl) (pH 8.0), 10 mM NaCl, 2% w/v sodium dodecyl sulfate (SDS)

Before start

Separate PCR-free facility


Extraction

- 1 Prior to extraction, individually vortex skin samples in a 10% commercial sodium hypochlorite solution (bleach) solution to decontaminate surface.
- 2 To remove the bleach subsequently vortexed the samples in 70% ethanol.
- 3 Finally vortex the samples in H₂O.
- 4 Add the skin samples to 1 mL digestion buffer consisting of 10 mM tris(hydroxymethyl)aminomethane hydrochloride (Tris-HCl) (pH 8.0), 10 mM NaCl, 2% w/v sodium dodecyl sulfate (SDS), 5 mM CaCl₂, 2.5 mM ethylenediaminetetraacetic acid (EDTA) (pH 8.0), 40 mM dithiothreitol (DTT), and 10% Proteinase K [Gilbert et al. 2007].

Note

Gilbert, M. T. P., Tomsho, L. P., Rendulic, S., Packard, M., Drautz, D. I., Sher, A., ... & Campos, P. F. (2007). Whole-genome shotgun sequencing of mitochondria from ancient hair shafts. *science*, 317(5846), 1927-1930


- 5 Incubate the samples for 12 hours at 56°C.

 12:00:00 Incubation at 56°C

- 6 In order to purify the DNA from contaminants, first, mix 1mL supernatant with 1mL phenol.

 1 mL Phenol


- 7 Vortex the sample for 20 sec.

 00:00:20 Vortexing


- 8 Gently rotate the sample for 5 min.

 00:05:00 Gentle rotation


- 9 Centrifuge the sample at 3000 g for 3 min.

 00:03:00 Centrifugation


- 10 Remove approximately 1mL aqueous liquid and mix with 1mL chloroform.

 1 mL Chloroform

11 Vortex the mixture for 30 sec.

 00:00:30 Vortex

12 Rotate the mixture for 5 min.

 00:05:00 Rotation

13 Centrifuge the mixture at 3000 g for 3 min.

 00:03:00 Centrifugation

14 Remove approximately 1mL aqueous liquid and purify using the MinElute PCR Purification kit (Qiagen, Valencia, CA) according to manufacturer's instruction with a slight modification: Firstly, modify the PB buffer according to [Allentoft et al. 2015].

Note

Allentoft, M. E., Sikora, M., Sjögren, K. G., Rasmussen, S., Rasmussen, M., Stenderup, J., ... & Malaspinas, A. S. (2015). Population genomics of Bronze Age Eurasia. *Nature*, 522(7555), 167-172

15 Secondly, increase the volume of PB binding buffer to 10x.

16 Apply the buffer to the spin columns following the method developed by [Dabney et al. 2013], use a Zymo-Spin V reservoir (Zymo Research, Irvine, CA) to pass the large buffer volume through the MinElute column.


Note

Dabney J, Knapp M, Glocke I, et al (2013) Complete mitochondrial genome sequence of a Middle Pleistocene cave bear reconstructed from ultrashort DNA fragments. *Proc Natl Acad Sci U S A* 110:15758–15763


17 Prior to the final centrifugation, add 15 µL of EB buffer to the column.

 15 µL EB Buffer

18 Incubate for 15 minutes at 37°C.

 00:15:00 Incubation at 37°C

19 Centrifugate at 6000 g for 1 min.

 00:01:00 Centrifugation



- 20 Quantify the extracted DNA using a Qubit fluorometer with a dsDNA high sensitivity (HS) assay (Life Technologies, Carlsbad, CA).