



Jun 09, 2018

Extraction of genomic DNA from diatoms by the modified method described in (Jacobs et al., 1992)

 [PLOS One](#)

DOI

dx.doi.org/10.17504/protocols.io.qh6dt9e

Artyom M. Marchenkov¹, Darya P. Petrova¹, Alexey A. Morozov¹, Yulia R. Zakharova¹, Michael A. Grachev¹, Alexander A. Bondar²

¹Limnological Institute of Siberian Branch of Russian Academy of Sciences;

²Institute of Chemical Biology and Fundamental Medicine of Siberian Branch of Russian Academy of Sciences



Artyom Marchenkov

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

[Create free account](#)

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.qh6dt9e>

External link: <https://doi.org/10.1371/journal.pone.0203161>

Protocol Citation: Artyom M. Marchenkov, Darya P. Petrova, Alexey A. Morozov, Yulia R. Zakharova, Michael A. Grachev, Alexander A. Bondar 2018. Extraction of genomic DNA from diatoms by the modified method described in (Jacobs et al., 1992). **protocols.io** <https://dx.doi.org/10.17504/protocols.io.qh6dt9e>

Manuscript citation:

Marchenkov AM, Petrova DP, Morozov AA, Zakharova YR, Grachev MA, Bondar AA (2018) A family of silicon transporter structural genes in a pennate diatom *Synedra ulna* subsp. *danica* (Kütz.) Skabitsch. PLoS ONE 13(8): e0203161. doi: [10.1371/journal.pone.0203161](https://doi.org/10.1371/journal.pone.0203161)

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 28, 2018

Last Modified: June 09, 2018

Protocol Integer ID: 12574

Keywords: diatom, DNA extraction , protocol for diatom dna extraction, diatom dna extraction, freshwater diatom, sb ras for freshwater diatom, extraction of genomic dna, dna extraction, extraction, genomic dna, dna

Abstract

A protocol for diatom DNA extraction used in the Limnological Institute SB RAS for freshwater diatoms.

Troubleshooting

Cells collecting

- 1 The cells were sedimented on polycarbonate filters (5- μ m pores) (Whatman, USA), briefly rinsed with cold Diatom Medium (DM), harvested by centrifugation for 2 min at 16,100 g at 4 °C and then stored at -70 °C.

Cells lysis

- 2 A 4.5 mL volume of lysis buffer (20 mM Tris-HCl, pH 8.0; 100 mM NaCl; 50 mM Na₂EDTA), 20% SDS (250 μ L) and proteinase K (250 μ L, 10 mg/mL) were added to 1 g of frozen cells, which were then incubated for 2 h at 60 °C with intermittent mixing. Then 250 μ L of proteinase K (10 mg/mL) were added again, and the cells were incubated for 40 min under the same conditions.

Extraction

- 3 The lysate was mixed for 10 min with 8 mL of phenol saturated with TE buffer, pH 8.0 using the tube rotator. The extract was centrifuged (10 min, 16,100 g, 20 °C), then the aqueous phase was taken and again mixed with an equal volume of phenol: chloroform: isoamyl alcohol mixture (25:24:1). It was then mixed and centrifuged under the same conditions. The aqueous phase was mixed with a chloroform: isoamyl alcohol mixture (24:1), mixed and centrifuged under the same conditions. Aqueous phase was used for the subsequent stages.

DNA precipitation

- 4 DNA was precipitated by adding 1/10 volume of 3 M sodium acetate, pH 5.2, and 1 volume of isopropanol to the aqueous phase. The solution was matured at 0 °C for 1 h and then centrifuged (30 min, 16,100 g, 4 °C). The supernatant was removed, and the precipitate was twice rinsed with 70% ethanol, dried in air and dissolved in sterile TE buffer (10 mM Tris-HCl; 1 mM Na₂EDTA, pH 8.0).