

Oct 17, 2023

## F/2 medium at 27 PSU of salinity from Sea Red salts

DOI

[dx.doi.org/10.17504/protocols.io.n92ldzyrxv5b/v1](https://dx.doi.org/10.17504/protocols.io.n92ldzyrxv5b/v1)



Estelle Bigeard<sup>1</sup>

<sup>1</sup>Station Biologique de Roscoff, France

Ecology of Marine Plank...

Roscoff Culture Collection

1 more workspace



Estelle Bigeard

Station Biologique de Roscoff, UMR7144 CNRS Sorbonne Univers...

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n92ldzyrxv5b/v1>

**Protocol Citation:** Estelle Bigeard 2023. F/2 medium at 27 PSU of salinity from Sea Red salts. **protocols.io**  
<https://dx.doi.org/10.17504/protocols.io.n92ldzyrxv5b/v1>

**License:** This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** January 07, 2022

**Last Modified:** October 17, 2023

**Protocol Integer ID:** 56659

**Keywords:** roscoff parasites culture collection, marine parasite, sea red salt, red sea salt, salinity, culture microalgae, fungi dinomyce, marine water enrichment solution, dinoflagellate, amoebophrya culture collection, psu of salinity, parasite

**Funders Acknowledgements:**

ANR AQUARADOX, PARALEX, EPHEMER

Grant ID: SU & CNRS

## Abstract

F/2 medium (Guillard's Marine Water Enrichment Solution) is used to maintain in culture microalgae (for example, dinoflagellates) and their marine parasites (Alveolates, Syndiniales & perkinsids, Fungi *Dinomyces*).

Most of these strains have been sampled from estuaries, at a 27 PSU of salinity. Adding soil extract is sometimes necessary to increase their growth. Since 2007, Roscoff Parasites Culture Collection was maintained in this medium.

In order to maintain the *Amoebophrya* Culture Collection with a medium which would be stable over time, we tested then validated since October 2022 the use of the Red Sea salt to prepare F/2 medium at a salinity of 27 PSU.

## Materials

### Equipments

Autoclave  
Laminar flow cabinet (or biosafety cabinet)  
Stainless Steel Filter Holder  
Peristaltic pump

### Materials

RED SEA Salt 22 kg - Sel pour aquarium – Available on the website « Aqua Store » - Ref R11065  
<https://www.aqua-store.fr/sel-aquarium-eau-de-mer/1238-red-sea-salt-22-kilos-livraison-incluse-7290100771990.html>

Guillard's (F/2) Marine Water Enrichment Solution 50 ×, liquid, plant cell culture tested - Merck Millipore (Sigma) - G9903

Stericup® Filter Unit - Merck Millipore - SCGVU10RE  
or Filter Unit PES 0,2µm 1000mL – Nalgene – 567-0020  
or 500 ml rapid-flow sterile bottle top filter with 0.2µ PES Membrane (Nalgene) - Thermo Fisher Scientific - 595-4520  
linked to a DURAN 1000ml glass bottle (use the same clean bottle)

10L Polycarbonate carboys (Nalgene) - ThermoScientific - 2251-0020

1L Polycarbonate bottle (Nalgene) - ThermoScientific - 2015-1000

Pipette

### Flow Cytometry

*SYBR*® Green I Nucleic Acid Gel Stain - Euromedex - 17590-1mL

## Troubleshooting

## Before start

### Red Sea Salt

Red Sea Salt is naturally harvested directly from the waters of the Red Sea. Natural seawater contains over 70 chemical elements. Most of them influence water parameters, but only a few play a significant role in its overall chemical stability.

The Red Sea contains moderately high yet balanced levels of Fundamental Elements (Calcium, Magnesium, Bicarbonates). These three "Fundamental Elements" have a major impact on the stability of pH, alkalinity, and the ionic strength of seawater. There are no nitrates or phosphates (algae nutrients), no toxic levels of heavy metals, nor chemical binders.

More informations to the link : <https://g1.redseafish.com/fr/red-sea-salts/red-sea-salt/>

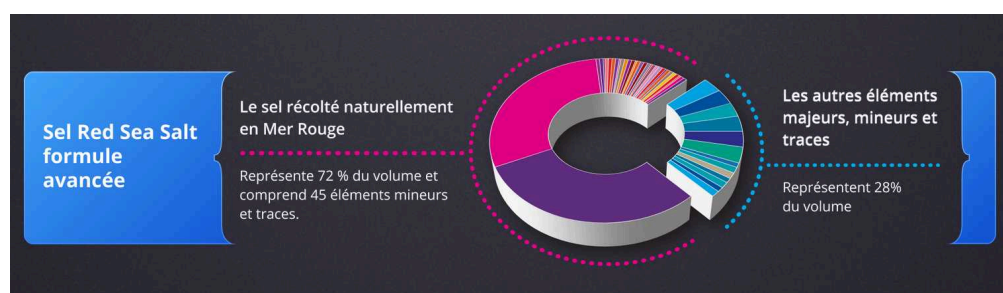


Figure 1 : Red Sea Salt Composition

## Preparation of the (Natural, Controlled) Seawater Solution

- 1 **Be careful** : Always use clean and non-toxic utensils for mixing.

**Note:** Due to possible salt stratification during transport, the entire salt should be mixed before sampling and dissolve it into water.

Use Milli-Q water as distilled water may contain some trace elements such as Ca, Mg, HCO<sub>3</sub>.

Check the salinity of the mixed seawater using a specific seawater refractometer or a conductivity meter. Ensure that the measuring equipment is properly calibrated before its use.

Transparent containers are used as opaque containers cannot be autoclaved at 121°C.

Salinity	pH	Alkalinity (°dKH)	Ca (mg/L)	Mg (mg/L)	K (mg/L)	Dose (g/L)
31.0 ppt	8,2-8,4	6,8-7,2	365-385	1090-1150	330-350	33,4
33,5 ppt	8,2-8,4	7,3-7,7	390-410	1170-1230	360-380	36,0
35,5 ppt	8,2-8,4	7,8-8,2	420-440	1250-1310	380-400	38,3

Figure 2 : Composition and salinity of seawater based on concentration.

### Seawater Salinity 27 PSU

To obtain seawater with a salinity of 27 PSU using Red Sea Salt, weigh 294g of salt and transfer it to a 10L Nalgene container.

Add 10 liters of Milli-Q water at room temperature.

Shake the carboy.

Autoclave at 121°C for 20 minutes.

**Comment :** There is no need to mix before autoclaving until complete dissolution. Heating will dissolve salts, even if salts will precipitate again during cooling.

## Preparation of Culture Medium under Sterile Conditions

- 2 Think to thaw F/2 medium (Stored at -20°C) at 4°C.

Add 20mL of F/2 medium per liter of seawater at least the day before filtration, up to 3 days in advance.

Perform a sterile filtration using a 0.22µm filter unit, up to a maximum of 5L.

Check the salinity using a salinity probe or another suitable device.

Store at room temperature.



## Sterility control using Flow Cytometry

- 3 Collect a sub-sample of 1 mL from 2 or 3 1L Nalgene culture bottles in a cytometry tube. Add 10 $\mu$ l of glutaraldehyde (25% grade) followed by 1% final concentration of SyBr Green for the DNA staining, wait 5 minutes at dark. Bacteria population can be estimated using a flow cytometer, based on their FSC and Green fluorescence under a 488 laser excitation.

If bacteria are detected, the medium should be filtered and tested again.