

Nov 13, 2025

## 🌐 Fast-ATACseq protocol – Stanford Bone Marrow TMC

DOI

<https://dx.doi.org/10.17504/protocols.io.3byl464b8go5/v1>

Patricia Favaro<sup>1</sup>, Warren Reynolds<sup>1</sup>, Reema Baskar<sup>1</sup>, Sean Bendall<sup>1</sup>

<sup>1</sup>Stanford University

Human BioMolecular Atlas Program (HuBMAP) Method Development Community  
Tech. support email: [Jeff.spraggins@vanderbilt.edu](mailto:Jeff.spraggins@vanderbilt.edu)



Patricia Favaro

Stanford University

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



**DOI:** <https://dx.doi.org/10.17504/protocols.io.3byl464b8go5/v1>

**Protocol Citation:** Patricia Favaro, Warren Reynolds, Reema Baskar, Sean Bendall 2025. Fast-ATACseq protocol – Stanford Bone Marrow TMC. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.3byl464b8go5/v1>

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** November 13, 2025

**Last Modified:** November 17, 2025

**Protocol Integer ID:** 232241

**Keywords:** Fast-ATAC-seq, HuBMAP, Tissue Map Center (TMC), Bone Marrow, bulk atac, atacseq protocol, enriched cell, purified cell, cell suspension, atac library, dna purification, atac, mobilized peripheral blood, cell, cocktail of surface antibody, surface antibody, staining media, seq analysis, peripheral blood

**Funders Acknowledgements:**

HuBMAP NIH

Grant ID: U54HL165445

## Abstract

Frozen CD34<sup>+</sup> enriched cells from mobilized peripheral blood were thawed and suspended in cell-staining media with TruStain FC blocker for 10 min at room temperature prior to staining. Next, the cocktail of surface antibodies was added (**Table S1**), and cells were stained in the dark for 30 min on ice and then washed in cell-staining media. Prior to data acquisition, cell suspensions were spiked with 7-AAD (BioLegend) to label non-viable cells. The sorted populations (**Figures S3B and S5D**) were used for bulk ATAC-seq analysis. For bulk Fast-ATAC-seq, cells were processed as described previously (**M.R. Corces et al., Nat. Genet., 48 (2016)**). Briefly, approximately 50,000 FACS-purified cells were pelleted and used for the transposition reaction for each technical replicate, followed by DNA purification. Amplification and purification of the transposed fragments were performed as described previously (**J.D. Buenrostro et al., Curr. Protoc. Mol. Biol., 109 (2015)**) with modified indexing primers (**J.D. Buenrostro et al., Nature, 523 (2015)**). Libraries were quantified using qPCR. All Fast-ATAC libraries were sequenced on an Illumina NovaSeq through Novogene. Each population analyzed included two technical replicates and two biological replicates.

## Troubleshooting

## Details

---

- 1 Table S1 (novel EMP and MP FACS Panel with the cocktail of surface antibodies, Figures S3B and S5D): <https://doi.org/10.1016/j.celrep.2025.115913>
  
- 2 ATAC-seq protocol:  
M.R. Corces et al., Nat. Genet., 48 (2016) - [10.1038/ng.3646](https://doi.org/10.1038/ng.3646)  
J.D. Buenrostro et al., Curr. Protoc. Mol. Biol., 109 (2015) - [.1002/0471142727.mb2129s109](https://doi.org/10.1002/0471142727.mb2129s109)  
J.D. Buenrostro et al., Nature, 523 (2015) - [10.1038/nature14590](https://doi.org/10.1038/nature14590)

## Acknowledgements

This work was supported by HuBMAP NIH grant U54HL165445. The study was also supported by a research grant from bluebird bio, which also provided the mPB used in this experiment.