

Oct 16, 2019

## Fatty acid extraction and derivatisation

DOI

[dx.doi.org/10.17504/protocols.io.79jhr4n](https://dx.doi.org/10.17504/protocols.io.79jhr4n)

Igem Dusseldorf<sup>1</sup>

<sup>1</sup>Heinrich-Heine Universität Düsseldorf



Igem Dusseldorf

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.79jhr4n>

**Protocol Citation:** Igem Dusseldorf 2019. Fatty acid extraction and derivatisation. **protocols.io**  
<https://dx.doi.org/10.17504/protocols.io.79jhr4n>

**Manuscript citation:**

Browse, John, Peter J. McCourt, and Christopher R. Somerville. "Fatty acid composition of leaf lipids determined after combined digestion and fatty acid methyl ester formation from fresh tissue." *Analytical biochemistry* 152.1 (1986): 141-145.

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



**Protocol status:** Working

**We use this protocol and it's working**

**Created:** October 15, 2019

**Last Modified:** October 16, 2019

**Protocol Integer ID:** 28683

**Keywords:** fatty acid extraction, pot fatty acid extraction, acid extraction, fatty acid, bacterial samples for preparation, derivatisation of plant, bacterial sample, extraction, weber lab for plant seedling, cyanobacteria, derivatisation this protocol, ms analysis, derivatisation, plant seedling, fatty

## Abstract

This protocol enables direct, one-pot fatty acid extraction and derivatisation of plant and bacterial samples for preparation for GC-MS analysis.

The protocol was established by the Weber lab for plant seedlings and seeds, but was successfully used for cyanobacteria and *E.coli* as well.

## Guidelines

All steps can be done at RT

## Materials

- C17 internal standard (stock = 1 mg/ml)
- MeOH//3N HCl (CAS: 7647-01-0)
- hexane (CAS: 110-54-3)
- 1% NaCl
- clean glass tubes

## Troubleshooting

## Safety warnings

! acrid, always wear safety goggles and gloves, don't use plastic

## Before start

Always use clean glass tubes and avoid using washing detergent




## Start

1 samples with 4 oD-units ( e.g.: oD = 1, you will need 4 ml culture )

you should make 4 or more replicates

2 centrifuge in a clean glass tube @ 4500 x g

 00:10:00

3 discard supernatant

4 Freeze @ -80 °C until further use

## Extraction solution

5 200 µl C17 internal standard (1 mg/ml hexane) + 9,8 ml MeOH/3N HCl


caution! fill in a beaker and take the needed amount out of this beaker, don't take out directly from original container to avoid contaminating stock solution.

**always work with gloves**


## Extraction

6 add 1 ml extraction solution to samples and blank  
For each new batch of samples, include a blank

7 heat @ 90°C  
after 5 minutes re-tighten lids of the glass tubes!

 01:00:00

8 let cool down @ RT

 00:15:00

9 add 1 ml hexane


5m

**Note**


hexane is very volatile, that's why working quickly is necessary

10 add 1 ml 1% NaCl

11 vortex

 00:00:30

12 spin down @ 2000 rpm

 00:05:00

13 transfer hexane phase in to GC vial with screw cap

14 dilute samples ( 10  $\mu$ l sample + 90  $\mu$ l hexane )  
wash hamilton between samples with hexane

all of these samples can be frozen @ -20 until further analysis