

Oct 15, 2020

Version 2

FCMPASS - Acquisition and gating of light scatter reference materials V.2



DOI

<https://dx.doi.org/10.17504/protocols.io.bnhimb4e>

Joshua Welsh¹, Jennifer Jones¹

¹Translational Nanobiology Section, Laboratory of Pathology, Center for Cancer Research, National Cancer Institute, National Institutes of Health

Translational Nanobiolo...



Jennifer Jones

NIH

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bnhimb4e>

Protocol Citation: Joshua Welsh, Jennifer Jones 2020. FCMPASS - Acquisition and gating of light scatter reference materials. protocols.io <https://dx.doi.org/10.17504/protocols.io.bnhimb4e> Version created by **Joshua A Welsh**

Manuscript citation:

Welsh J A, Jones J C, Small Particle Fluorescence and Light Scatter Calibration Using FCMPASS Software, Current Protocols in Cytometry, 94, e79. doi: [10.1002/cpcy.79](https://doi.org/10.1002/cpcy.79)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](https://creativecommons.org/licenses/by/4.0/), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 15, 2020

Last Modified: June 29, 2023

Protocol Integer ID: 43274

Keywords: flow cytometry, calibration, extracellular vesicles, data from light scatter reference material, light scatter reference materials this protocol, gating of light scatter reference material, fcmpass software, fcmpass software package, light scatter reference material, use with the fcmpass software, fcmpass, performing small particle calibration, small particle calibration,

Disclaimer

This protocol summarizes key steps for a specific type of assay, which is one of a collection of assays used for EV analysis in the NCI Translational Nanobiology Section at the time of submission of this protocol. Appropriate use of this protocol requires careful, cohesive integration with other methods for EV production, isolation, and characterization. By using the FCMPASS software you agree to the following terms and conditions.

Terms & Conditions of use for FCMPASS software.

Definitions: The term "SOFTWARE" throughout this agreement means the machine readable, binary, object code form, and the related documentation for FCMPASS, a software package that is designed to allow flow cytometer calibration for small particles. The term "RECIPIENT" means the party that downloads the software. The term "PROVIDER" means the National Cancer Institute (NCI), a participating institute of the National Institutes of Health (NIH), and an agency of the United States Government. By downloading or otherwise receiving the SOFTWARE, RECIPIENT may use the SOFTWARE subject to RECIPIENT's agreement to the following terms:

1. THE SOFTWARE SHALL NOT BE USED IN THE TREATMENT OR DIAGNOSIS OF HUMAN SUBJECTS. RECIPIENT is responsible for compliance with all laws and regulations applicable to the use of the SOFTWARE.
2. RECIPIENT shall not distribute the SOFTWARE, in whole or in part without express advance written approval of PROVIDER.
3. The SOFTWARE may be used for research, academic, and educational purposes only. The SOFTWARE may not be used for commercial purposes. RECIPIENT will not license or sell or use the SOFTWARE for commercial purposes or applications.
4. The SOFTWARE that is distributed pursuant to this Agreement has been created by United States Government employees. In accordance with Title 17 of the United States Code, section 105, the SOFTWARE is not subject to copyright protection in the United States. Other than copyright, all rights, title and interest in the SOFTWARE shall remain with the PROVIDER.
5. RECIPIENT shall not modify, extend, decompile, make derivatives of, merge, publish, reverse engineer or distribute the SOFTWARE without written permission from PROVIDER.
6. RECIPIENT may publish or otherwise publicly disclose the results of using the SOFTWARE. RECIPIENT agrees to acknowledge PROVIDER's contribution of the SOFTWARE in all written publications containing any data or information regarding or resulting from use of the SOFTWARE.
7. THE SOFTWARE IS PROVIDED "AS IS" AND ANY EXPRESS OR IMPLIED WARRANTIES, INCLUDING, BUT NOT LIMITED TO, THE IMPLIED WARRANTIES OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE AND NONINFRINGEMENT ARE DISCLAIMED. IN NO EVENT SHALL THE PROVIDER OR THE INDIVIDUAL DEVELOPERS BE LIABLE FOR ANY DIRECT, INDIRECT, INCIDENTAL, SPECIAL, EXEMPLARY, OR CONSEQUENTIAL DAMAGES (INCLUDING, BUT NOT LIMITED TO, PROCUREMENT OF SUBSTITUTE GOODS OR SERVICES; LOSS OF USE, DATA, OR PROFITS; OR BUSINESS INTERRUPTION) HOWEVER



CAUSED AND ON ANY THEORY OF LIABILITY, WHETHER IN CONTRACT, STRICT LIABILITY, OR TORT (INCLUDING NEGLIGENCE OR OTHERWISE) ARISING IN ANY WAY OUT OF THE USE OF THIS SOFTWARE, EVEN IF ADVISED OF THE POSSIBILITY OF SUCH DAMAGE. PROVIDER makes no representations that the use of SOFTWARE will not infringe any patent or proprietary rights of third parties.

8. RECIPIENT may, on an as-needed basis, send to PROVIDER reports regarding the application of the SOFTWARE and the effectiveness and problems encountered in using the SOFTWARE, without disclosing RECIPIENT's confidential information. Information from general reports may be used by the PROVIDER to enhance the capabilities of the SOFTWARE. Reports can be forwarded to the PROVIDER at one of the following addresses: joshua.welsh@nih.gov or jennifer.jones2@nih.gov
9. No indemnification for any loss, claim, damage, or liability is intended or provided by either Party under this Agreement. Each Party shall be liable for any loss, claim, damage, or liability that said Party incurs as a result of said Party's activities under this Agreement, except that Provider, as an agency of the United States, assumes liability only to the extent as provided under the Federal Tort Claims Act (28 U.S.C. Chapter 171 Sections 2671-2680).
10. RECIPIENT agrees not to claim, infer, or imply endorsement by the United States Government, or any of its organizational units, contractors or employees. RECIPIENT agrees not to use any trademarks, service marks, trade names, logos or product names of NCI or NIH to endorse or promote products derived from the SOFTWARE without specific, prior and written permission.
11. Title in the SOFTWARE shall remain with the PROVIDER. It is understood that nothing herein will be deemed to constitute, by implication or otherwise, the grant to either party by the other of any license or other rights under any patent, patent application or other intellectual property right or interest. PROVIDER reserves the right to distribute the SOFTWARE to others and to use it for PROVIDER's own purposes. The United States Government explicitly retains all rights to use the SOFTWARE for any purpose, to have it used on the Government's behalf or to allow others to use it.


Abstract

This protocol outlines the steps required to collect data from light scatter reference materials for use with the FCMPASS software. This is one of a number of protocols in the pipeline for performing small particle calibration using the fcompass software package.

Materials

MATERIALS


 Flow Cytometer

 Vortex

 Low Protein Binding Collection Tubes (1.5 mL) **Thermo Fisher Scientific Catalog #90410**

 DPBS **Invitrogen - Thermo Fisher Catalog #14190**

 NIST-Traceable Polystyrene Beads (3000 Series) **Thermo Fisher Scientific**

 Falcon® 5 mL Round Bottom Polystyrene Test Tube with Snap Cap Sterile Individually Wrapped
Corning Catalog #352003

- 1 Calculate the stock traceable size calibration reference bead particle concentration using percent solids value and particle density provided by the manufacturer and the following formula, where N_P is the concentration (particles mL⁻¹), $W_V\%$ is the percent solids, ρ is the particle density (g mL⁻¹), and D is the average diameter (μm).

$$N_P = \frac{(W_V\% \times 6 \times 10^{12})}{(\pi \rho D^3)}$$

Note

For example, 100 nm polystyrene beads at 1% with 1.05 g mL⁻¹ would be calculated using:

$$1.82 \times 10^{13} = \frac{(0.01 \times 6 \times 10^{12})}{(\pi \times 1.05 \times 0.1^3)}$$

A online calculator can also be found at this [link](#).

- 2 Thoroughly vortex the traceable size calibration reference bead stock bottles to homogenize the mixtures before dispensing 1 drop (~50 μL) into separate 500 μL low-protein binding Eppendorf.
- 3 Using the working stock from step 2, make up 500 μL solution at 1×10^7 particles mL⁻¹.

Note

It is recommended that serial dilutions are used and volumes of no less than 10 μL to avoid pipetting errors. The optimal particle concentration at which to run the reference materials will vary depending on several factors, including the flow rate, beam height, and electronic sampling rate. If running for the first time, it is recommended that serial dilutions are performed to determine the optimal concentration for preparation of the beads.


- 4 On the flow cytometer, set the triggering threshold to the most sensitive light scatter detector and ensure the parameter is using log-scaling (not linear or biexponential).
- 5 Running DPBS, lower the triggering threshold until the noise floor of the instrument becomes visible. This is most clearly when using a histogram.
- 5.1 Plotting the trigger-channel height parameter against time and monitoring while running DPBS is a good indication for determining whether an instrument is clean. If the spread

of noise (and event rate) decreases over time, it is indicative that the instrument was dirty and is becoming cleaner.

Note


There are a number of cleaning solutions. Some shared resource lab managers have a preference. These include bleach, contrad 70, micro 90, surfanol

- 5.2 The extent to which the opto-electronic noise of an instrument can be sampled will vary between instruments. Legacy flow cytometers will tolerate a couple of 1000-2000 events/second whilst allowing room to sample desired events, while high-speed jet-in-air sorters are capable of sample 10,000+ events per second.
- 5.3 Triggering using a light scatter parameter on the opto-electronic noise of the instrument has benefits in determining and tracking the lower limit of detection, as well as being informative for buffer + reagent controls where background fluorescence will show clear shifts due to many events being triggered from sampling the noise. The use of this method comes at the cost of having high event rates and therefore larger files. Before utilizing this method the instrument should be validated to determine: 1) its ability to detect and accurately process particles, 2) the event rate at which single small particles are detected, and 3) the degree to which the opto-electronic noise can be sampled without creating artefacts or reducing the ability to detect genuine events.
- 5.4 On some instruments that utilize peristaltic pumps there can appear to be an increase and decrease of the baseline corresponding to the turnover of the pump. This is a result of the threshold being set close to (but above) the electronic noise, resulting in the increase and decrease in trigger events in light scatter. This can be overcome by lowering the threshold so that the noise is being sampled regardless of the peristaltic pump turnover or increasing the threshold and therefore decreasing the instrument's limit of sensitivity.
- 6 Analyze each bead sample at the same acquisition settings until >5000 bead events are recorded.
- 6.1 It is preferable to analyze and store bead populations individually. This will minimize population overlap, aggregates, background noise, and artifacts.

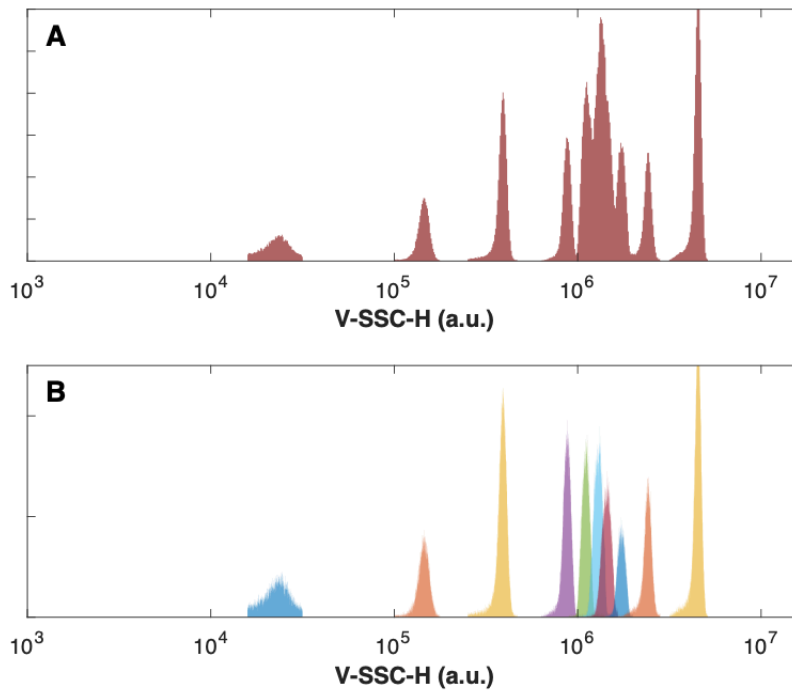

Panel A demonstrates the cumulative distribution of the gated populations when mixed together. While some populations are clearly distinguished some are not. The areas where bunching of populations occurs is dependent upon the cytometer and is useful in determining the collection angle. Panel B illustrates overlaid and colored gated bead population from Oanel A.

- 7 Gate each bead population using the parameter Height vs. Area in a dot-plot to remove doublets/aggregates and then use a histogram on the light scatter parameter (Height) to obtain statistics for each population. The light scatter parameter should use log scaling.

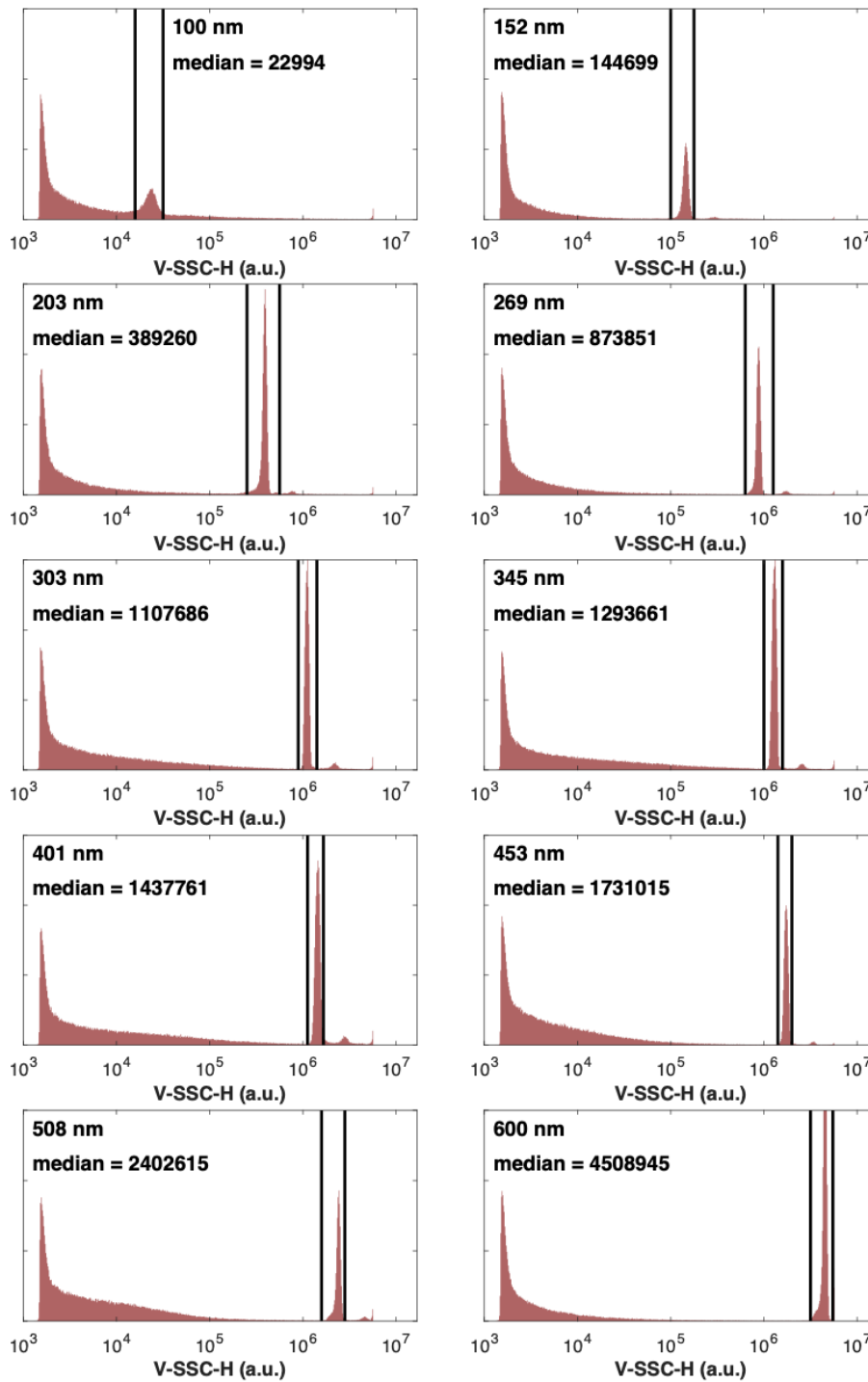

Gating light scatter reference beads. Each panel shows the gating of polystyrene NIST-traceable reference beads ranging in mean diameter from 100 to 600 nm. The median light scatter statistic of the gated population is given in each panel

8 Obtain the median statistic for each of the bead populations.



- 8.1 By default, flow cytometers trigger the acquisition of an event using the pulse height parameter. In cases where a trigger threshold is being defined (e.g. SSC), it is recommended that the pulse-height is used so that the limit of detection can be defined in calibrated units. There is no consensus within the small particle community over the use of pulse height vs. area. We recommend that, in general, if the parameter being calibrated was not used as a trigger channel the pulse area statistic should be used due to the tendency for low signal intensities to be linear and therefore a more reliable method for extrapolation.