

Aug 28, 2024

Fear Conditioning

DOI

<https://dx.doi.org/10.17504/protocols.io.261ge56bdg47/v1>

Sabina Marciano^{1,2}, Roberta Marongiu^{1,2}

¹Department of Neurological Surgery, Weill Cornell Medical College, New York, NY 10065;

²Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD 20815, USA



Jacquelyn Haytayan

Weill Cornell Medicine

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.261ge56bdg47/v1>

Protocol Citation: Sabina Marciano, Roberta Marongiu 2024. Fear Conditioning. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.261ge56bdg47/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: April 04, 2024

Last Modified: September 23, 2024

Protocol Integer ID: 97806

Keywords: ASAPCRN, motor impairment, mouse, behavior, fear learning, memory in mice, conditioning behavioral, mice

Funders Acknowledgements:

Aligning Science Across Parkinson's

Grant ID: 020608

Abstract

Behavioral assay to measure fear learning and memory in mice.

Materials

- MedAssociates fear conditioning chambers
- white floor inserts
- black ceiling inserts
- peppermint and limonene spray



Fear Acquisition

- 1 Habituate mice to the room. Set up fear conditioning chambers as well as protocols (listed below) in VideoFreeze software.
- 2 Place individual mice in a lit fear conditioning apparatus scented with peppermint spray.
- 3 Protocol: 3m 20s
 - 00:02:00 baseline followed by 5 tone-shock pairings where the shock follows the tone
 - shock = 0.5mA, 00:00:01 is delivered after the end of the tone (75dB, 2kHz, 00:00:19)
 - empty trace intervals between the shocks last 00:01:00
- 4 After the protocol runs for each mouse, place the mouse back in the home cage and clean the fear conditioning chamber thoroughly.
- 5 Contextual test is performed 24:00:00 after the fear acquisition test. Cued Test is performed at least 02:00:00 after the contextual test. 1d 2h

Contextual Test

- 6 Goal: Measure the subjects' fear response to the environment where fear conditioning was received
- 7 24 hours after fear acquisition: Mice are placed in a lightened FC apparatus scented with peppermint spray.
- 8 Protocol: Neither cue nor shock are delivered during this test session (Duration 00:07:30) 7m 30s

Cued Test



9 Goal: Measure the subjects fear response to the CS, and ideally no other features of the training environment. The key is to alter the chamber so the subjects are exposed to a new set of environmental cues.

10 Mice are placed in a lightened FC apparatus which is altered by placing a smooth floor insert on shock grid and black triangles on top.

Limonene spray is used, instead of peppermint scent.

11 Protocol:

- Baseline: 2 min followed by five tone-cues
- No shock is delivered after the end of the tone (75dB, 2 kHz,  00:00:19)

19s

Primary Analyses

12 Fear conditioning acquisition

- % Component freezing for baseline, each tone (if used), shock, and interstimulus interval
- Graphs for each component should be graphed separately
- Average motion

Context

- % Component freezing per minute and total
- Per minute and total will be on separate graphs
- Average motion per minute and total

Cued

- %Component freezing for baseline, each tone, and interstimulus interval
- Graphs for each component should be graphed separately
- Average motion for baseline, each tone, and interstimulus interval