

Sep 08, 2021

FindingNemo Library 2: Modified RAD004

DOI

<https://dx.doi.org/10.17504/protocols.io.bxv4pn8w>



Inswasti Cahyani¹, John Tyson², Nadine Holmes¹, Josh Quick³, Nicholas Loman³, Matthew Loose¹

¹University of Nottingham; ²University of British Columbia; ³University of Birmingham



Inswasti Cahyani

University of Nottingham

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bxv4pn8w>

Protocol Citation: Inswasti Cahyani, John Tyson, Nadine Holmes, Josh Quick, Nicholas Loman, Matthew Loose 2021. FindingNemo Library 2: Modified RAD004. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.bxv4pn8w>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: September 01, 2021

Last Modified: September 08, 2021

Protocol Integer ID: 52892


Keywords: ultra-long sequencing, cohex, glass bead, nanopore, MinION, UHMW DNA, Monarch, Circulomics, phenol, SDS, CTAB, GM12878, Whatman, PromethION, Nanobind, modified rad004 protocol, uhmw dna, nanopore sequencer, principle of the rad004 kit, rad004 kit, ul library protocol, findingnemo library, rad004, based adapter ligation kit, dna, adapter ligation kit, findingnemo

Abstract

This sub-protocol is designed to prepare library from extracted ultra-high molecular weight (UHMW) DNA to obtain ultra-long (UL) reads on Nanopore sequencers. The UL library protocol we tested here is based on ONT's rapid kit, *i.e.*, **SQK-RAD004**, a transposase based adapter ligation kit.

This **Modified RAD004** protocol consistently produced N50 > 100 kb from a good input quality of UHMW DNA, for when ULK001 is not accessible/available. Transposase-based reaction is done in a large volume of up to 1 ml.

The working principle of the RAD004 kit is shown in the diagram below:

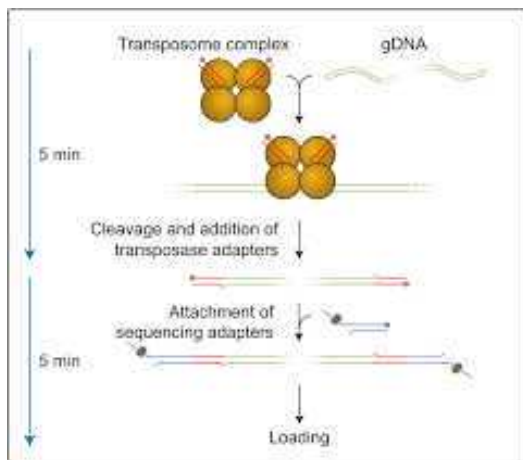

source: <https://store.nanoporetech.com/uk/rapid-sequencing-kit.html>



Guidelines

Acknowledgements

The FindingNemo protocol series was developed by Inswasti Cahyani for the Long Read Club with significant contributions from John Tyson and Nadine Holmes, also Josh Quick, and continuous discussion and support of Matt Loose and Nick Loman. We would also like to thank Giron Koetsier (NEB) and Kelvin Liu (Circulomics) for lending their expertise and advance product trials.

Please follow on Twitter for latest updates and results:

@NininUoN

@mattloose



Materials

Chemicals/Compounds

-  Tris-HCl pH 8.0 **Thermo Scientific Catalog #J22638-AE**
-  Nuclease-free Water **Thermofisher Catalog #AM9920**
-  NaCl (5 M) RNase-free **Thermo Fisher Scientific Catalog #AM9759**
-  Magnesium Chloride **Fisher Scientific Catalog #AC223210010**
-  Triton X-100 **Merck MilliporeSigma (Sigma-Aldrich) Catalog #T8787-50ML**
-  Glycerol **Bio Basic Inc. Catalog #GB0232.SIZE.500ml**

Made-up Buffer

4x MuA Buffer

- 100 mM Tris-HCl pH 8.0
- 40 mM MgCl₂
- 440 mM NaCl
- 0.2% TritonX-100
- 40% Glycerol

Kits

-  Rapid Sequencing Kit **Oxford Nanopore Technologies Catalog #SQK-RAD004**

Disposables

-  DNA LoBind Tubes, 1.5 mL **Eppendorf Catalog #0030108051**
-  DNA LoBind 2.0ml PCR Clean Eppendorf Tubes **Eppendorf Catalog #0030 108.078**
- Wide-bore (or cut off) P1000 and P200 tips



Safety warnings

- ⚠ When handling phenol always wear PPE, keep a solution of 50% (w/v) PEG-400 nearby to treat the burn in the case of accidental splashes.

Before start

Things to observe at all times:

- Excessive and vigorous pipetting and vortexing should be avoided as these may shear the DNA.
- Make up buffers with nuclease-free water to avoid introducing nucleases to solutions.
- Avoid unnecessary heating and freezing; isolated DNA should be stable for storage in the fridge for months

Library Prep Notes

- 1 Extracted UHMW DNA is often difficult to quantify due to its viscosity. However, accurate measurement of DNA concentration is crucial for calculating optimum ratio of the transposase enzyme to the DNA molecules.

We provide a protocol section for quantifying UHMW DNA in our 'FindingNemo' protocol master file.

Properly quantified DNA can then be processed for this library prep.

Both cell number and DNA concentration/amount are used to calculate the amount of transposase (FRA) and adapter (RAP).

We base this protocol on SQK-RAD004 but follow the SQK-ULK001 protocol for the optimum ratio of transposase amount to human genomic DNA:

6 µl FRA to 6 million human cells (or around 40 µg DNA)

For other species, the genome size has to be taken into account and the FRA to DNA ratio optimised, *e.g.*, we had optimised a non-human cell line of **6.2 Gb genome** at:

2.5 µl FRA to 1 million non-human cell (around 12-15 µg DNA)

- 2 The following table lists the replaced reagents between ULK001 and RAD004.

	A	B	C
	Reagent	ULK001	RAD004 (modified)
	Dilution buffer	FDB	MuA buffer (see Materials)
	Transposase	FRA*	FRA^
	Sequencing adapter	RAP-F	RAP

(*) FRA enzyme in ULK001 kit

(^) FRA enzyme in RAD004 kit

Transposase Reaction

- 3 In a 2 ml tube, dilute UHMW DNA to a concentration of around 50 ng/µl in a total volume of 750 µl (with water or elution buffer if required).
Mix well with a P1000 wide-bore tip.



**Note**

- DNA concentration can still range from 20-50 ng/μl to have optimum transposase reaction.
- If input DNA amount is less than 20 μg (1-3 million cells used), halve all the reaction volumes, *i.e.*, 375 μl total DNA volume instead of 750 μl as in the table below.
- It is important to have as homogeneous DNA as possible at this step so the transposase can access and cut the DNA molecules with an even distribution. It is OK to pipette thoroughly but gently.

	A	B	C	D	E
	Cell No. (million)	Approx. DNA amount (μg)	Total DNA volume (μl)	DNA concentrati on (ng/μl)	Total reaction volume (μl)
	6	>20-40	750	20-50	1000
	5				
	4				
	3	5-20	375		500
	2				
	1				

- 4 In a 1.5 ml tube, dilute the corresponding amount of transposase (FRA) with MuA buffer to a total volume of 250 μl (or 125 μl if doing half-reaction). More details in the table below.

	A	B	C	D	E
	Cell No. (million)	Approx. DNA amount (μg)	FRA (μl)	4X MuA Buffer (μl)	Total reaction volume (μl)
	6	>20-40	6	244	1000
	5		5	245	
	4		4	246	
	3	5-20	3	122	500
	2		2	123	

	A	B	C	D	E
	1		1	124	

- 5 Mix the diluted FRA by vortexing for 2-3 seconds.
- 6 Using a P1000 wide-bore tip, add the diluted FRA to the DNA sample.
Stir the reaction with the pipette tip whilst expelling the diluted FRA to ensure an even distribution.
Mix thoroughly by gentle pipetting.

 On ice

- 7 Incubate the reaction as follows:

 23 °C  00:10:00

 70 °C  00:05:00

 Room temperature  00:10:00 at least

Note

It is important that the room temperature at the fragmentation step (first incubation step) does not fall below 20°C to ensure optimum reaction condition. The use of a water bath or heating block is recommended.

Adapter Ligation

- 8 Add the corresponding volume of sequencing adapter (RAP) as in the table below.

	A	B	C	D	E
	Cell No. (million)	Approx. DNA amount (µg)	FRA (µl)	RAP (µl)	Total reaction volume (µl)
	6	>20-40	6	5	1000
	5		5	4.2	
	4		4	3.3	



	A	B	C	D	E
	3	5-20	3	2.5	500
	2		2	1.7	
	1		1	0.8	

Note

- Use a P1000 wide-bore tip to pipette mix. Visually check to ensure the reaction is thoroughly mixed.
- Tube inversion can be used to aid mixing.

9 Incubate for 30 minutes at 23°C.



23 °C



00:30:00