



Jun 20, 2025

Version 4

🌐 Flaviviruses (West Nile, Zika, Dengue serotypes 1-4) NS2B-NS3 protease fluorescence dose response and single point assay V.4

DOI

<https://dx.doi.org/10.17504/protocols.io.q26g7yebkgwz/v4>

Haim Barr^{1,2}, Noa Lahav^{1,2}

¹Weizmann Institute of Science; ²ASAP Discovery Consortium

ASAP Discovery



Noa Lahav

Weizmann Institute

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.q26g7yebkgwz/v4>

Protocol Citation: Haim Barr, Noa Lahav 2025. Flaviviruses (West Nile, Zika, Dengue serotypes 1-4) NS2B-NS3 protease fluorescence dose response and single point assay. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.q26g7yebkgwz/v4> Version created by **Mary-Ann Xavier**

Manuscript citation:

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 04, 2025

Last Modified: June 20, 2025

Protocol Integer ID: 219547

Keywords: Flaviviridae, Fluorescence assay, NS2B-NS3, ZIKV, WNV, DENV, ASAP, flavivirus infection, flavivirus, ns3 protease fluorescence dose response, compounds top assay concentration, biochemical assay, compounds assay concentration, dengue, top assay concentration, ns3 protein, treatments for viral infectious disease, assay buffer, direct enzyme activity measurement method, viral infectious disease, enzyme, zika, assay reliability, point assay, assay, total assay volume, stock to assay, ns3 protease, fluorescence intensity, fluorescence, protein

Funders Acknowledgements:

National Institutes of Health/National Institute Of Allergy and Infectious Diseases (NIH/NIAID)

Grant ID: U19AI171399

Disclaimer

The content is solely the responsibility of the authors and does not necessarily represent the official views of the National Institutes of Health.

Abstract

This is a **functional, biochemical assay** used to identify treatments for viral infectious diseases related to **Flavivirus infection**, (specifically **West Nile, Zika, and Dengue**) and targets the conserved **NS2B-NS3 protein**. In this current version, we added the Addgene id.

Utilizing a direct enzyme activity measurement method, the experiment was performed in a 384-well plate reading the fluorescence intensity. This assay tested the mode of action of inhibition. Initial screening was conducted with 2 compound concentrations to determine %inhibition, followed by dose response for IC₅₀ measurement (as specified below). In each run a reference compound was added to verify the assay reliability and robustness.

Assay was developed at the Weizmann Institute of Science, as a part of the ASAP Discovery Consortium.

Experiment Concentrations (From Stock to Assay)

	A	B	C	D	E
	Reagent	Stock	Concentration Loaded into CERTUS	Final Concentration in Assay Plate	Units
	Substrate	10000	10	5	μM
	DENV NS2B-NS3	217000	200	100	nM
	ZIKV NS2B-NS3	225000	200	100	nM
	WNV NS2B-NS3	222000	200	100	nM

Assay Buffer

	A	B	C	E
	Reagent	Stock	Final concentration	Units
	Tris pH=8.5	1000	20	mM
	Glycerol	50	10	%
	Triton	10	0.01	%



Compound Plate Design for Dose Response:

Total assay volume: 20 μ L

Compounds top assay concentration: 100 μ M

Dilution factor: 2

Dose response points: 12

Number of replicates: 2

Backfill with DMSO: yes

Compounds Plate Design for 2-Point Assay:

Total Assay Volume: 20 μ L

Compounds Assay Concentration: 100 μ M and 50 μ M

Dilution Factor: 2

Dose Response Points: 2

Number of Replicates: 2

Backfill with DMSO: Yes



Materials

Assay Buffer Reagents (Concentration listed are Stock Solution Concentrations)

1. [M] 50 Mass / % volume

⊗ Glycerol - for molecular biology, ≥99% **Merck MilliporeSigma (Sigma-Aldrich) Catalog #G5516**

2. [M] 1000 millimolar (mM)

⊗ Tris(2-carboxyethyl)phosphine hydrochloride **Merck MilliporeSigma (Sigma-Aldrich) Catalog #75259**

3. [M] 10 % (v/v) Triton ⊗ Triton X-100 **Merck MilliporeSigma (Sigma-Aldrich) Catalog #T8787**

*Note: all components are added fresh to the assay buffer before each experiment

Additional Reagents:

West Nile Virus (WNV) Reagents:

[M] 222000 nanomolar (nM) WNV NS2B/NS3 Enzyme

- WNV NS2B/NS3 was originally [M] 222000 nanomolar (nM) and was diluted to [M] 200 nanomolar (nM) with freshly made **Assay Buffer** before each experiment

[M] 10000 nanomolar (nM) WNV Enzyme Substrate

- Enzyme Substrate** was ⊗ Boc-Gly-Arg-Arg-AMC acetate salt **Biosynth Catalog #FB110553**
- Substrate stock** was created by dissolving the substrate powder in **DMSO** to create [M] 10 millimolar (mM) Substrate Stock aliquots and stored at 🌡️ -80 °C . Before each experiment, the Substrate Stock was diluted to [M] 10 micromolar (μM) Substrate with freshly made **Assay Buffer**.

Zika (ZIKV) Virus Reagents:

[M] 225000 nanomolar (nM) ZIKV NS2B/NS3 Enzyme

- ZIKV NS2B-NS3 was originally [M] 225000 nanomolar (nM) and was diluted to [M] 200 nanomolar (nM) with freshly made **Assay Buffer** before each experiment

[M] 10000 nanomolar (nM) ZIKV Enzyme Substrate

- Enzyme Substrate** was ⊗ Boc-Gly-Arg-Arg-AMC acetate salt **Biosynth Catalog #FB110553**
- Substrate stock** was created by dissolving the substrate powder in **DMSO** to create [M] 10 millimolar (mM) Substrate Stock aliquots and stored at 🌡️ -80 °C . Before each experiment, the **Substrate Stock** was diluted to [M] 10 micromolar (μM) Substrate with freshly made **Assay Buffer**.

Dengue (DENV) Reagents:

[M] 217000 nanomolar (nM) DENV NS2B/NS3 Enzyme

- DENV NS2B/NS3 was originally [M] 217000 nanomolar (nM) and was diluted to [M] 200 nanomolar (nM) with freshly made **Assay Buffer** before each experiment

**[M] 10000 nanomolar (nM) DENV Enzyme Substrate**

- **Enzyme Substrate** was  Bz-Nle-KRR-AMC (hydrochloride) **Cayman Chemical Company Catalog #27710**
- **Substrate Stock** was created by dissolving the substrate powder in **DMSO** to create **[M] 10 millimolar (mM) Substrate Stock aliquots** and stored at  **-80 °C** . Before each experiment, the **Substrate Stock** was diluted to **[M] 10 micromolar (μM) Substrate** with freshly made **Assay Buffer**.

Expected result

A literature reference inhibitor (**compound 83**, Journal of Medicinal Chemistry **2015** 58 (23), 9354-9370 <https://doi.org/10.1021/acs.jmedchem.5b01441>) was added to each run of NS2B-NS3 to give IC50 values of ~2 μM for DENV-2 and WNV and ~3 μM for ZIKA .

Protocol materials

-  ZIKV VE_Ganxian2016 NS2B-NS3 protease fusion (non-cleavable linker, N-term Twin-Strep-tag) **addgene Catalog #228661**
-  DENV-1 strain US/Hawaii/1944 NS2B-NS3 protease fusion with non cleavable linker **addgene Catalog #228648**
-  DENV-2 Thailand/16681/84 NS2B-NS3 protease fusion (non-cleavable linker) **addgene Catalog #228650**
-  WNV NY99 NS2B-NS3 protease fusion (non-cleavable linker) K→A mutation to reduce autoproteolysis **addgene Catalog #228660**
-  DENV-3 Philippines/H87/1956 NS2B-NS3 protease fusion (non-cleavable linker) **addgene Catalog #228651**
-  DENV-4 strain H241 NS2B-NS3 protease fusion (non-cleavable linker) **addgene Catalog #228652**

Safety warnings

-  Please be sure to wear proper Personal Protective Equipment (PPE) while performing this experiment.

Before start

Dummy plate: In each step of the protocol, dispense the reagents into an **empty** plate before dispensing them into the assay compounds plate (this will detect and prevent dispensing issues).

Note: Inhibitor compounds stock concentration is 20 mM. Compounds are pre-dispensed into 384 plates and stored at -20°C until use.



Flavivirus NS2B-NS3 protease expression and purification

- 1 Zika virus NS2B-NS3 fused protease expression and purification of a construct suitable for biochemical assay. Using the plasmid



ZIKV VE_Ganxian2016 NS2B-NS3 protease fusion (non-cleavable linker, N-term Twin-Strep-tag) **addgene Catalog #228661**

Protocol

NAME

Zika NS2B-NS3 fusion construct for assay small scale expression and purification protocol

CREATED BY

korvus.wang

Preview

- 2 West Nile virus NS2B-NS3 fused protease expression and purification of the construct suitable for biochemical assay.



WNV NY99 NS2B-NS3 protease fusion (non-cleavable linker) K→A mutation to reduce autoproteolysis **addgene Catalog #228660**

Protocol

NAME

West Nile virus NS2B-NS3 protease fusion construct small scale expression and purification protocol

CREATED BY

korvus.wang

Preview



- 3 Dengue virus serotype 1 virus NS2B-NS3 fused protease expression and purification of the construct suitable for biochemical assay.



DENV-1 strain US/Hawaii/1944 NS2B-NS3 protease fusion with non cleavable linker **addgene Catalog #228648**

Protocol

NAME

**Dengue virus serotype 1 NS2B-NS3 protease fusion construct
small scale expression and purification protocol**

CREATED BY

korvus.wang

Preview

- 4 Dengue virus serotype 2 virus NS2B-NS3 fused protease expression and purification of the construct suitable for biochemical assay.



DENV-2 Thailand/16681/84 NS2B-NS3 protease fusion (non-cleavable linker) **addgene Catalog #228650**

Protocol

NAME

**Dengue virus serotype 2 NS2B-NS3 protease fusion construct
small scale expression and purification protocol**

CREATED BY

korvus.wang

Preview



- 5 Dengue virus serotype 3 virus NS2B-NS3 fused protease expression and purification of a construct suitable for biochemical assay.



DENV-3 Philippines/H87/1956 NS2B-NS3 protease fusion (non-cleavable linker) **addgene Catalog #228651**

Protocol



NAME

Dengue virus serotype 3 NS2B-NS3 protease fusion construct using parallel rapid protein expression and purification: 100 mL cultures

CREATED BY

korvus.wang

Preview

- 6 Dengue virus serotype 4 virus NS2B-NS3 fused protease expression and purification of a construct suitable for biochemical assay.



DENV-4 strain H241 NS2B-NS3 protease fusion (non-cleavable linker) **addgene Catalog #228652**

Protocol

NAME

Dengue virus serotype 4 NS2B-NS3 protease fusion construct: 100 mL cultures

CREATED BY

korvus.wang

Preview

Determine which Flavivirus is needed and prepare solutions

7 Prepare solutions based on the materials section.

	A	B	C	D	E
	Reagent	Stock	Loaded into CERTUS	Final in assay plate	units
Choice of NS2B-NS3 Enzyme Protein					
	DENV NS2B-NS3	217000	200	100	nM
	ZIKV NS2B-NS3	225000	200	100	nM
	WNV NS2B-NS3	222000	200	100	nM
Choice of Viral Substrate					
	WNV/ZIKV Substrate	10000	10	5	μM
	DENV Substrate	10000	10	5	μM
Assay buffer					
	Tris pH=8.5	1000	20	20	mM
	Glycerol	50	10	10	%
	Triton	10	0.01	0.01	%



Prepare 384-well Plate for experiment

2h 31m

- 8 **PRIME** the dispenser with **Assay Buffer** by selecting the **PRIME** button until the tubes are filled completely.
- 8.1 **DISPENSE**  10 μL Assay Buffer to Columns **1** and **23** of assay plate
Note: These will represent the **inhibitor control columns** (Contain: Substrate + Assay Buffer + DMSO, no enzyme, no compounds)
- 8.2 **EMPTY** the dispenser tubes.
▪ Discard Assay Buffer discharged from the dispenser tubes.
- 9 **PRIME** the dispenser with  200 nanomolar (nM) Flavivirus NS2B/NS3 by selecting the **PRIME** button until the tubes are filled completely.
- 9.1 **DISPENSE**  10 μL  200 nanomolar (nM) Flavivirus NS2B/NS3 to columns **2** through **22** and column **24**.
Note:
▪ Be sure to cycle dispensing on an empty plate first, to detect and avoid dispensing issues.
▪  200 nanomolar (nM) Flavivirus NS2B/NS3 is two times the assay concentration. The final concentration of the Flavivirus NS2B/NS3 in the assay plate is  100 nanomolar (nM) .
▪ Columns 2 and 24 are **neutral control columns** (Contain: Enzyme + substrate + DMSO, no compounds)
- 9.2 **EMPTY** the dispenser tubes .
▪ Discard the  200 nanomolar (nM) Flavivirus NS2B/NS3 discharged from the tubes
- 10 **CENTRIFUGE** plate  1500 rpm, 00:01:00 to remove bubbles
- 11 **INCUBATE** plate  02:00:00 at  Room temperature
▲ Make sure the plate is protected from light!
- During Incubation:** wash and prepare the dispenser to the next step
- 12 **PRIME** the dispenser with  200 μL  10 micromolar (μM) Flavivirus Substrate



- **Note:** Be sure to cycle dispensing on an empty plate first, to detect and avoid dispensing issues.

13 **DISPENSE**  10 μL of **[M] 10 micromolar (μM) Flavivirus Substrate** to **Columns 1-24** (the full plate).

- **Note:** **[M] 10 micromolar (μM) Flavivirus Substrate** is two times the assay concentration. The final concentration of the Flavivirus Substrate in the assay plate is **[M] 5 micromolar (μM)**.

14 **CENTRIFUGE** plate  1500 rpm, Room temperature, 00:01:00 to remove bubbles

1m

15 **INCUBATE** plate for  00:30:00 at  Room temperature

30m

▲ Make sure the plate is protected from light!

Recommended: Clean the dispenser during incubation

Read Plate Fluorescence

- 16 **READ** and **RECORD** the plate Relative fluorescence units (RFU) via the "**Flavivirus protocol**" on the **PERAstar FS Control Software**.
- Software is a standard Fluorescence Assay set for Optimal excitation wavelength 360 nm, emission wavelength 470 nm, and a Gain of 200.

Equipment

PERAstar FS

Microplate reader

BMG LABTECH

0471B0001A

https://www.bmglabtech.com/en/pherastar-fsx/?utm_term=pherastar%20plate%20reader&utm_campaign=usa.roi.products&utm_source=FwOoFR_5EUHUaAlkREALw_wcB



Expected result

gain 300 should yield ~20,000 RFU in full reaction; ~7000 RFU in Buffer control