

Dec 01, 2023

Flow cytometry

DOI

dx.doi.org/10.17504/protocols.io.ewov1qm47gr2/v1

Pranay Srivastava¹

¹Massachusetts General Hospital, ASAP



Fang Zhou

Massachusetts General Hospital, Aligning Science Across Park...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.ewov1qm47gr2/v1>

Protocol Citation: Pranay Srivastava 2023. Flow cytometry. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.ewov1qm47gr2/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 30, 2023

Last Modified: May 31, 2024

Protocol Integer ID: 91669

Keywords: ASAPCRN, flow cytometry this protocol, immune cell profile inn spleen, flow cytometry, cell

Abstract

This protocol is to assess immune cell profile inn spleen (Chauhan et al. 2018).

Troubleshooting

- 1 Spleen was removed in a 35 mm petri plate with 5 ml RPMI 1640 and digested mechanically and passed through 70 µm filter screen.
- 2 The cell suspension was centrifuged, and the pellet was incubated in RBC lysis buffer. The resulting cell suspension was washed in 1xPBS and blocked with Fc Block (Biolegend, Cat# 101302, 1 µl/50 µl).
- 3 The cells were incubated with MC1R (Invitrogen, Cat# PIPA521911, 1.39 µg) antibody followed by fluorophore conjugated primary antibodies for extracellular markers (Biolegend Cat# 101235, CD11b-BV421(0.25 µg); Cat# 127641, Ly6G-BV-650 (0.25 µg); Cat# 128041, Ly6C-BV785 (0.125 µg); Cat# 100516, CD4-APC (0.25 µg); Cat# 100751, CD8a-BV510 (0.5 µg); Cat# 152405, CD19-PerCP-Cy5.5 (0.25 µg); Cat# 102036, CD25-BV605 (0.3 µg)) and AF488 (Invitrogen, Cat# A11034, 1:200).
- 4 Zombie dye (Biolegend, Cat# 423101, 1µl/sample) was used to differentiate between live and dead cells. Helper T cells and cytotoxic T cells were identified by CD4+ and CD8+, respectively. CD4+CD25+ cells were used to mark Tregs. CD19+ cells were used as marker for B cells. Monocytes were identified as CD11b+Ly6G-Ly6C^{high} cells and neutrophils were marked by CD11b+Ly6C-Ly6G+. For monocytes and neutrophils, CD11b-positive cells were first extracted from the live cell subset by expansion with SSC-A, followed by Ly6C. After expansion with Ly6C and Ly6G, we excluded Ly6G-positive cells. Monocytes were identified as CD11b+Ly6G-Ly6C high cells, and neutrophils were marked by CD11b+Ly6C-Ly6G+.
- 5 SORP 5 Laser BD Fortessa X-20 (BD Bioscience) and FlowJo v10.7.1 (Becton Dickson & Company) software was used for data acquisition.