


Sep 30, 2024

🌐 Fluorescein diacetate assay - for plastic degrading enzymes in algae

DOI

<https://dx.doi.org/10.17504/protocols.io.n2bvj3j9blk5/v1>

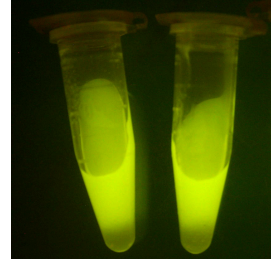

Joao Vitor Molino¹

¹University of California, San Diego

Lab



Joao Vitor Molino

Ronin Institute, University of California, San Diego

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n2bvj3j9blk5/v1>

Protocol Citation: Joao Vitor Molino 2024. Fluorescein diacetate assay - for plastic degrading enzymes in algae. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.n2bvj3j9blk5/v1>



License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 21, 2023

Last Modified: September 30, 2024

Protocol Integer ID: 91215

Keywords: Fluorescein diacetate FDA assay, plastic degrading enzymes in algae fluorescein, enzymes in algal culture, algae fluorescein, plastic degrading enzyme, enzymes in microalgae, degrading enzyme, enzymatic activity of plastic, sampling algal culture, hydrolysis assay, using fluorescein diacetate, algal culture, enzyme, fluorescein diacetate, enzyme activity of cell, enzyme activity, applications in environmental biotechnology, environmental biotechnology, enzymatic activity, fluorescence, assay with phosphate buffer, activity assay, based activity assay, cell contamination in the sample, microalgae, cell contamination, assay, hydrolysis, plastic waste management, clear supernatant, green glow

Funders Acknowledgements:

APEX

Grant ID: EE0009671 UCSD

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to **[protocols.io](#)** is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with **[protocols.io](#)**, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

Fluorescein diacetate (FDA) hydrolysis assays can be used to measure the enzyme activity of cells in a sample. A bright yellow-green glow is produced and is strongest when enzymatic activity is greatest. This can be quantified using a spectrofluorometer or a spectrophotometer (**REF**).

In this protocol, we describe a method for assessing the enzymatic activity of plastic-degrading enzymes in algal cultures using Fluorescein Diacetate (FDA) hydrolysis. The assay involves sampling algal cultures, centrifuging to obtain clear supernatants, and performing a fluorescence-based activity assay. Key steps include ensuring no cell contamination in the samples, preparing the assay with phosphate buffer at optimal pH, and conducting kinetic readings using a plate reader. This protocol is crucial for identifying efficient plastic-degrading enzymes in microalgae, with applications in environmental biotechnology and plastic waste management

Guidelines

Objective: This assay measures the enzymatic activity of plastic-degrading enzymes in algal cultures using fluorescein diacetate (FDA) hydrolysis.

Note: Esterases, including native enzymes from cells, can show activity in this assay. Avoid cell or lysed cell contamination in the sample.

Safety warnings

1. **Chemical Handling:** Always handle chemicals, including fluorescein diacetate and acetone, in a well-ventilated area while wearing appropriate personal protective equipment (PPE) such as gloves and safety goggles.
2. **Biological Material:** Treat algal cultures as potential biohazards. Follow your laboratory's biosafety procedures when handling and disposing of biological samples.
3. **Centrifuge Safety:** Ensure the centrifuge is balanced before use to avoid equipment damage or injury.
4. **Instrument Safety:** Operate the plate reader and other instruments according to the manufacturer's guidelines to prevent damage and ensure accurate readings.
5. **Waste Disposal:** Dispose of chemical and biological wastes according to your laboratory's waste disposal protocols.



Before start

1. **Algal Culture Sample:** For measuring enzyme activity.
2. **96-Well Clear Plate:** For sample measurement.
3. **Centrifuge:** To separate supernatant from cells.
4. **100 mM Phosphate Buffer (pH 8.0):** For assay preparation.
5. **Fluorescein Diacetate (FDA):** 2 mM solution in acetone.
6. **Plate Reader:** Capable of measuring absorbance at 750 nm, chlorophyll fluorescence at 440/680 nm, and FDA fluorescence at 490/yyy nm.
7. **Acetone:** For FDA solution preparation.
8. **Pipettes and Tips:** For accurate liquid transfer.
9. **Excel Software:** For data analysis.
10. **FITC Standard Curve:** For calculating activity per mol.



Checking cell density

- 1 Sample  160 μL of culture in a 96-well-clear plate.
- 2 Measure the absorbance at 750 nm and Chlorophyll fluorescence at 440/680nm. For example, in Infinite® M200 PRO plate reader, set the gain to 100, sensor distance 18000um, and top reading.

Sample preparation

13m

- 3 Spin down  1 mL of a sample at max speed (e.g.,  20000 g) for at least  00:03:00 .
- 4 Transfer the supernatant to a clean tube or use it quickly (e.g., less than a few minutes ~  00:10:00) to avoid any cells getting back into the solution. Also, handle samples with care to avoid mixing.

3m

10m

Assay preparation

10m

- 5 Add  89 μL of Phosphate buffer  100 millimolar (mM) ,  8 to a well. *The buffer can vary and depends on the enzyme being assayed. PHL7 has max activity at pH 8 and can be assayed even in highly concentrated buffer solutions such as  1 Mass Percent* Also, any buffering system works as long as the pH is buffered around pH 8, as Tris. Each enzyme has its own particularity.
- 6 Add  10 μL of sample to the well (e.g., supernatant)
- 7 Measure the absorbance at 750 nm and Chlorophyll fluorescence at 440/680nm. Step to check cell carried over, RFU - Relative Fluorescence units of >100 in Chlorophyll reading indicates cell or lysate present, which interferes with the activity assay. Values close to 100 can still be informative but need to be checked with care in the analysis step
- 8 Incubate plate @  37 °C for  00:10:00 .

10m

Assay performance

- 9 Set the plate reader to  37 °C prior to the reading activity. Enzymes display different activities in different temperatures, so we need to standardize one and keep it constant. PHL7 actually has an optimum temperature of  70 °C . Typically, plate readers can't handle that temperature.
- 10 Add  1 µL of FDA  2 millimolar (mM) in acetone to each well. Use multichannel for several wells. FDA concentration can be used in different values. For our assay, it just needs to be in excess so the substrate does not limit the reaction speed. We can be sure it was not by the shape of the kinetic curve. If a plateau is reached, it means the substrate was consumed, and at the steps before the plateau, the reaction was limited by the substrate reducing the apparent enzyme activity in that region of the graph. Acetone's final concentration shouldn't be higher than  1 % (v/v) , ideally.
- 11 Immediately after adding FDA, start reading the activity kinetics RFU FDA . In an Infinite  M200 PRO plate reader set ex/em to 490/520 nm, gain 40, sensor distance 18000um, top reading Kinetic cycle: 1 read every minute for 30 cycles

12

Data Analysis

- 13 Convert the FDA RFU to $\sqrt{\text{RFU}}$ to all data points. The square root is used to transform the data to make it linear. RFU increase in a reaction of order 2, meaning a quadratic function. Probably due to 2 bonds needing to be broken by the enzyme to restore FITC fluorescence.
- 14 Add the sample name to the Excel file for the specific well
- 15 Discard results in which a chlorophyll signal higher than 100 was detected in the activity experiment well. Cells containing esterase also generate signal in the experimental setup, interfering with the measurement.
- 16 Calculate "=Linest" to each sample in Excel
- 17 Calculate activity per mol by using the FITC standard curve



18 Save the result