



Apr 26, 2024

Version 1

Fluorescence assay for Enterovirus A71 3C protease activity measurement V.1

DOI

dx.doi.org/10.17504/protocols.io.14egn6zjql5d/v1

Charline Giroud¹, Oleg Fedorov¹

¹Center for Medicines Discovery, University of Oxford

ASAP Discovery



Charline Giroud

Center for Medicines Discovery

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.14egn6zjql5d/v1>

Protocol Citation: Charline Giroud, Oleg Fedorov 2024. Fluorescence assay for Enterovirus A71 3C protease activity measurement. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.14egn6zjql5d/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: April 26, 2024

Last Modified: April 26, 2024

Protocol Integer ID: 98851

Keywords: FRET assay, enzymatic assay, IC50 assay, protease assay, Picornaviridae , Enterovirus, enterovirus a71 3c protease activity measurement, activity of viral protease, enterovirus a71, viral protease, 3c protease activity measurement, detection of the fluorescence, protease activity measurement, cleavage of the peptide, sequence specific to the tested protease, fluorescence emission, fluorescence, 3c protease, tested protease, peptide, assay, labelled peptide, protease, proximity of dabcyI

Funders Acknowledgements:

National Institutes of Health/National Institute Of Allergy and Infectious Diseases (NIH/NIAID)

Grant ID: U19AI171399

Disclaimer

The content is solely the responsibility of the authors and does not necessarily represent the official views of the National Institutes of Health.

Abstract

This protocol details the fluorescence assay for Enterovirus A71 (EV-A71) 3C protease activity measurement. This method is intended to measure the activity of viral proteases by using a specific labelled-peptide that allows the detection of the cleaved product. The substrate contains the cleavage-sequence specific to the tested protease and is labeled in C-terminal by the fluorophore Edans (ex 336 nm; em: 455 nm) and in N-terminal by the quencher Dabcyl (ex 472 nm). In the case of a non-cleaved substrate, the proximity of Dabcyl to Edans prevents the emission and the detection of the fluorescence at 455 nm. The cleavage of the peptide by the protease allows Edans' fluorescence emission and detection.

Attachments



[nyjwb5id7.docx](#)


203KB


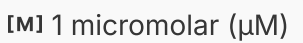
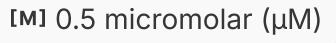
Materials




Reagents

■ Assay buffer:

	A	B
	Tris pH 7.0	50 mM
	NaCl	150 mM
	Glycerol	10%
	DTT (optional)	0.5 mM

■ **Incubation:**  01:00:00 at  Room temperature .

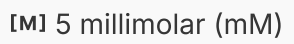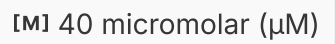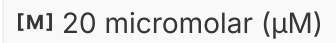■ **EV-D68 3C:** Protein stocks were stored at  -80 °C and used as 2x solution ( 1 micromolar (μM) ,  0.5 micromolar (μM) final assay concentration) in assay buffer.

■ **Substrate:** Dabcyl-KEALFQGPPQFE-Edans (LifeTein, USA) prepared as a stock solution at  5 millimolar (mM) in DMSO and used at 2x solution ( 40 micromolar (μM) ,  20 micromolar (μM) final concentration assay concentration) in assay buffer.

■ **Positive control:** GC376 (Pubchem CID 71481119),  50 micromolar (μM) top final assay concentration.

■ **Plates:** ProxiPlate-384 Plus, white, Greiner cat# 6008280.

■ **Liquid handler:** Echo® acoustic liquid handler (Beckman Coulter, USA).

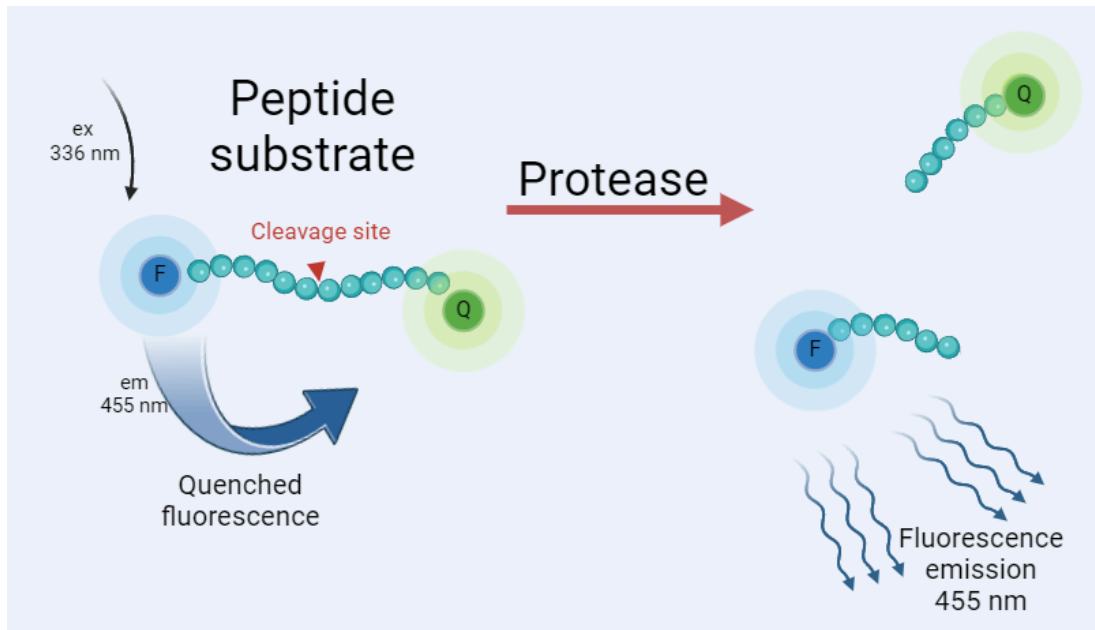
■ **Plate reader:** Pherastar FS, BMG Labtech (Germany), 350-460 FI optic module, the plate is read every 30 s for 2 hours and shacked during 5 s before the first reading.

Troubleshooting

EV-A71 3C Pro Assay

3h

- 1 This method is intended to measure the activity of viral proteases by using a specific labelled-peptide that allows the detection of the cleaved product. The substrate contains the cleavage-sequence specific to the tested protease and is labeled in C-terminal by the fluorophore Edans (ex 336 nm; em: 455 nm) and in N-terminal by the quencher Dabcyl (ex 472 nm). In the case of a non-cleaved substrate, the proximity of Dabcyl to Edans prevents the emission and the detection of the fluorescence at 455 nm. The cleavage of the peptide by the protease allows Edans' fluorescence emission and detection.



Asset URL:

EV-A71 3C Pro IC50 Measurement

3h

- 2 **Assay buffer:** 50 millimolar (mM) Tris pH 7.0, 150 millimolar (mM) NaCl, 10 % (v/v) glycerol and 1 millimolar (mM) TCEP (optional).
- Incubation:** at .
- EV-A71 3C:** protein stocks were stored at -80C and used as 2x solution (5 micromolar (μ M) , 2 2.5 micromolar (μ M) final assay concentration) in assay buffer.
- Positive control:** GC376 (Pubchem CID 71481119), 50 μ M top final assay concentration.

3h 0m 35s



Substrate: Dabcyl-KEALFQGPPQFE-Edans (LifeTein, USA) prepared as a stock solution at 10 mM in DMSO and used at 2x solution [M] 20 micromolar (μM) (

[M] 10 micromolar (μM) final concentration assay concentration) in assay buffer.

Plates: Non-binding, black 384-plate, Greiner for the assay.

Liquid handler: Echo® acoustic liquid handler (Beckman Coulter, USA).

Plate reader: Pherastar FS, BMG Labtech (Germany), 350-460 FI optic module, the plate is read every ⌚ 00:00:30 for ⌚ 02:00:00 and shaken ⌚ 00:00:05 before each reading.

3 50 μL of 2x protein solution were added to each well containing the compounds to be tested previously dispensed onto the plate.

4 Incubate the mix for ⌚ 01:00:00 at 🌡 Room temperature and initiate the enzymatic reaction by adding of 50 μL of 2x ([M] 10 micromolar (μM)) substrate solution using the plate reader injector.

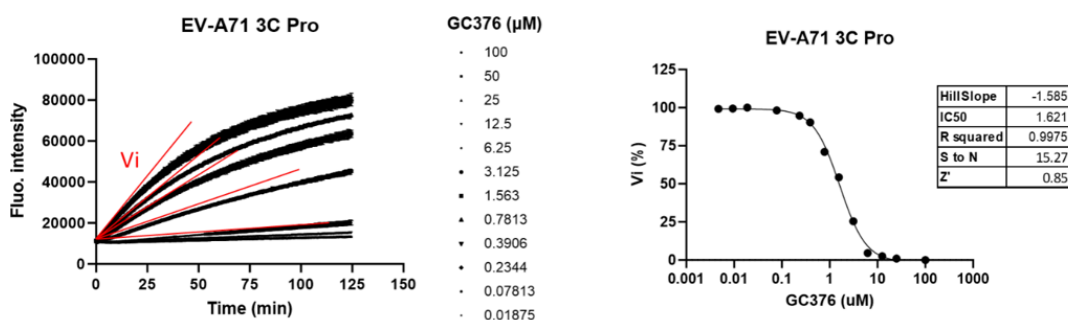
1h



5 Read the fluorescence intensity at 350/460 nm every 30 seconds for ⌚ 02:00:00 in kinetic mode, which includes a shaking step of the plate before each measurement.

2h

6 Calculate the IC₅₀ by plotting the initial velocity against various concentrations of tested inhibitors by using a four-parameter dose-response curve in Prism (v8.0) software.



Asset URL: