



May 23, 2023

Fluorescence in situ hybridization (FISH)

DOI

dx.doi.org/10.17504/protocols.io.kqdg39mmeg25/v1

Peter Sims¹

¹Columbia University



Ngan Kim Tran

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kqdg39mmeg25/v1>

Protocol Citation: Peter Sims 2023. Fluorescence in situ hybridization (FISH). **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.kqdg39mmeg25/v1>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 23, 2023

Last Modified: May 31, 2024

Protocol Integer ID: 82320

Keywords: ASAPCRN, fluorescence in situ hybridization, situ hybridization, fluorescence, fish

Funders Acknowledgements:

ASAP

Grant ID: ASAP-000375

Abstract

Fluorescence in situ hybridization (FISH) mRNA

Troubleshooting

- 1 Mouse brain sections were prepared, removed from cryoprotectant solution, and washed three times in tris-buffered saline (TBS) at room temperature.
Sections were incubated with hydrogen peroxide (ACD) for 15 min at room temperature, washed several times in TBS, and then mounted to Superfrost slides.
Sections were allowed to dry for 10 min and a hydrophobic barrier (PAP pen, Vector Labs) was created around the tissue.
Tissue was incubated in 50% EtOH, then 70% EtOH, then 100% EtOH for 5 min each.
Sections were rehydrated in TBS for several minutes, digested with **Protease**IV (ACD) for 25 min at room temperature, and rinsed twice with TBS before proceeding to the RNA Scope Multiplex Fluorescent v2 assay (ACD).
- 2 The RNA Scope Multiplex Fluorescent v2 assay was conducted according to the manufacturer's instructions, with all incubations taking place in a humidified chamber at 40 °C.
Two 5-minute washes in excess RNA Scope Wash Buffer (ACD) took place between each incubation in sequential order: probes (2-hours), AMP1 (30 min), AMP2 (30 min), AMP3 (15 min), HRP-C1/2/3 (15 min), TSA Cy3 (30 min), HRP blocker (30 min), HRP-C1/2/3 (15 min), and TSA Cy5 (30 min).
Samples were washed twice more in RNA Scope Wash Buffer, then twice more in TBS.
Samples were then blocked and immunostained for **tyrosine hydroxylase**.
After **immunostaining**, samples were mounted in Fluoromount G and stored at 4 °C for up to 1 week before imaging.