



Dec 02, 2021

Fluorescence microscope (Zeiss)

DOI

dx.doi.org/10.17504/protocols.io.b2iwqcfe

x.chu¹

¹Imperial College London

Behavioural Genomics



x.chu

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.b2iwqcfe>

Protocol Citation: x.chu 2021. Fluorescence microscope (Zeiss). **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.b2iwqcfe>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: December 02, 2021

Last Modified: December 02, 2021

Protocol Integer ID: 55606

Keywords: fluorescence microscope in the lab, fluorescence microscope, fluorescence, fluorescence mode, microscope, zeiss, time in fluorescence mode, operating lamp, led radiation, lab


Abstract

How to operate the fluorescence microscope in the lab. Beware of safety hazards: This instrument uses led radiation. Do not stare at the operating lamp for a prolonged period of time in fluorescence mode.

Troubleshooting

Microscope layout

1

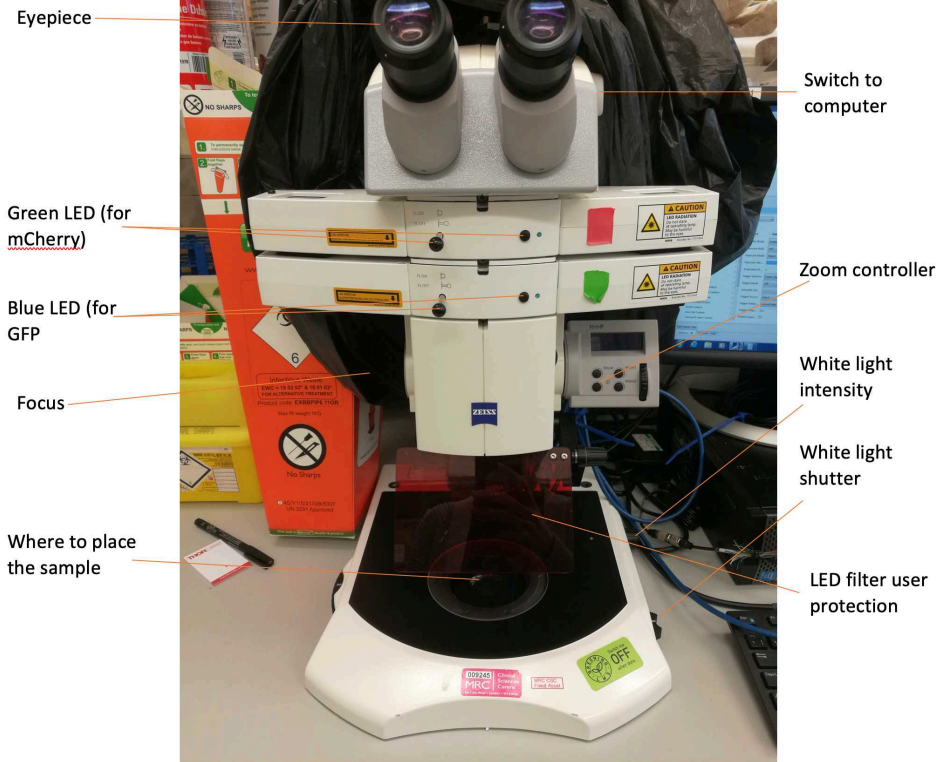

Schematics of microscope

Operating microscope

- 2 Turn on the white light intensity and shutter
- 3 Turn on the zoom controller
- 4 Turn on the LED, if using fluorescent mode
- 5 Turn on the software on the computer: Pylon viewer

- 6 Place sample in the microscope and adjust the eyepiece
- 7 Focus the sample using the wheel on the left-hand side whilst observing through the eyepiece
- 8 Using the white light shutter on the right to block out the white light, only fluorescence from the sample reaches the eyepiece.
- 9 To capture an image, switch to computer mode using the switch on the upper right-hand side.
- 10 Press the camera symbol in the software to capture an image. Parameters such as acquisition time can be set in the software.
- 11 The best results are obtained when the plastic cover is on and the background light in the room is kept at a minimum.
- 12 Turn off the white light, LED, and zoom controller when finished using the microscope.
- 13 Enclose with plastic cover.

Trouble-shooting

- 14 The software does not connect to the camera:
Make sure the USB cable is plugged in to a USB 3.0 port on the computer. Check that the camera is on. Make sure the correct driver is installed.
- 15 The zoom controller, white light or LED don't turn on:
Make sure that the microscope is plugged in.
- 16 Nothing can be seen through the eyepiece:
Make sure nothing is blocking the path between the sample and the objective lens. Make sure that the microscope is not switched to camera view.