

Jan 31, 2024

Fontana-Masson staining

DOI

<https://dx.doi.org/10.17504/protocols.io.n92ldmxoxl5b/v1>

Cristian González-Cabrera¹, Csilla Novák¹, Andrés auricio Jaramillo Flautero¹, Celine Winter¹, Matthias Prigge¹

¹Neuromodulatory Network Group, Leibniz Institute for Neurobiology, Magdeburg

TeamPrigge



Team Prigge

Leibniz Institute for Neurobiology

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n92ldmxoxl5b/v1>

Protocol Citation: Cristian González-Cabrera, Csilla Novák, Andrés auricio Jaramillo Flautero, Celine Winter, Matthias Prigge 2024. Fontana-Masson staining. protocols.io <https://dx.doi.org/10.17504/protocols.io.n92ldmxoxl5b/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: January 31, 2024

Last Modified: May 31, 2024

Protocol Integer ID: 94459

Keywords: Neuromelanin, Staining, pigments, ASAPCRN, masson staining protocol, other pigments in tissue sample, pigment analysis in tissue, pigment analysis, staining, other pigment, tissue sample, melanin, silver nitrate working solution, fontana, thorough washing steps between each incubation, staining protocol

Funders Acknowledgements:

ASAP

Grant ID: 020505

Abstract

The Fontana-Masson Staining Protocol is a detailed method for staining free-floating sections (20-50 microns) to detect melanin and other pigments in tissue samples. The procedure involves preparing a silver nitrate working solution, treating sections with triton, and sequential incubation in pre-heated silver nitrate solution, gold chloride, and thiosulfate solutions, with thorough washing steps between each incubation. The process concludes with sections being placed in PBS for mounting and counterstaining. This protocol is critical for histological studies in biology, particularly for pigment analysis in tissues.

Materials

Silver nitrate solution

Ammonium hydroxyde (concentrated)

PBS with 0.3% triton

Gold chloride solution

Thiosulfate solution

Distilled water

Tap water

Fontana-Masson staining

3h 14m 30s

- 1 Use free floating sections between 20-50 microns.
- 2 Silver Nitrate working solution:
 - Prepare a 1:3 solution of silver nitrate solution in distilled water. Prepare enough solution for the amount of sections you want to stain.
 - Add drop by drop concentrated ammonium hydroxide, while shaking, until the solution gradually turns transparent. Be careful and stop the addition of ammoniacal solution as soon as you get transparency.

10m





3 Place the sections to stain (up to 7-8 sections) in a  2 mL Eppendorf tube. Incubate the sections in 0.3% triton for two hours at room temperature while shaking

2h

4 Wash 3 times with distilled water for 5 min each.

15m

5 Pre-heat the silver nitrate working solution for 5 minutes at  60 °C .

5m

6 Incubate the sections in the preheated solution for 20 minutes at  60 °C .

20m

7 Wash the sections 3 times with abundant distilled water.

8 From now on, perform the following steps in individual sections:

8.1 Incubate the section with gold chloride solution. Apply enough solution to cover the section (approximately  100-150 μL). Manually shake the tube for 30 seconds.

30s

8.2 Wash the section 3 times with abundant distilled water.

5m

8.3 Incubate the section with thiosulfate solution. Apply enough solution to cover the section (approximately  100–150 μL). Manually shake the tube for 2 minutes.

2m

8.4 Wash with tap water for 2 minutes.

2m

8.5 Wash 3 times with distilled water for 5 min each.

15m

9 Place the sections in PBS for mounting, counterstaining, etc.

Results

10 Slices should now show an amplified pigmentation as show in figures

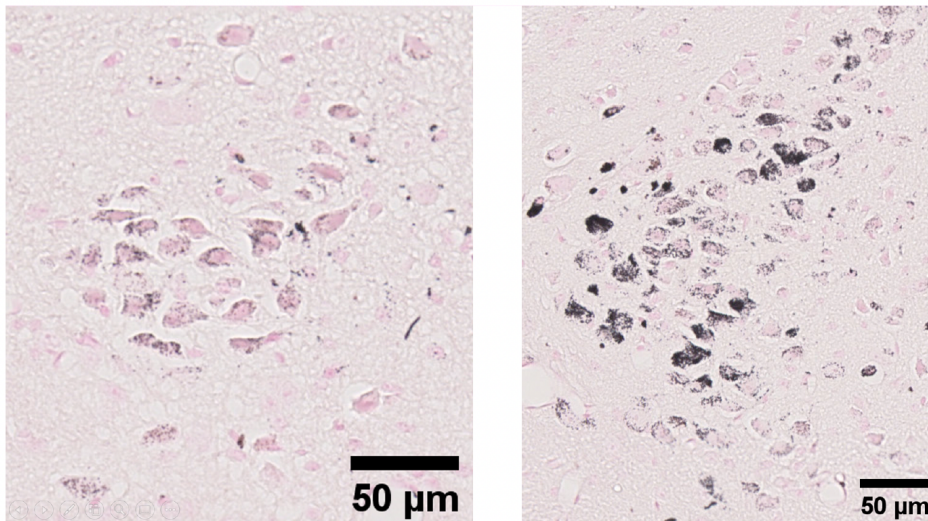


Fig. 6 Presence of neuromelanin in the left locus coeruleus of AAV-hTyr-injected mice 4449 (left) and 4448 (right).