

May 16, 2024

Version 2

Fragment Analyzer Operation for PCR Products V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.5jyl8mxwrg2w/v2>

Allen Institute for Brain Science¹

¹Allen Institute

Allen Institute for Brain Science
Tech. support
[Click here to message tech. support](#)



Allen Institute

Allen Institute

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.5jyl8mxwrg2w/v2>

Protocol Citation: Allen Institute for Brain Science 2024. Fragment Analyzer Operation for PCR Products. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.5jyl8mxwrg2w/v2> Version created by [Allen Institute](#)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: May 16, 2024

Last Modified: May 16, 2024

Protocol Integer ID: 99981

Keywords: MB0073, quality control, QC, genotyping, polymerase chain reaction, products, fragment analyzer operation for polymerase chain reaction, fragment analyzer operation for pcr product, fragment analyzer by advanced analytical, genotyping pcr product, fragment analyzer operation, fragment analyzer, pcr product, pcr, sample concentration, sample,

Abstract

This protocol describes fragment analyzer operation for polymerase chain reaction (PCR) products. The Fragment Analyzer by Advanced Analytical is used as a quality control step for genotyping PCR products. Peak size, shape, and sample concentration are recorded and used to evaluate samples on a pass/fail level.

Note: Research reported in this publication was supported by the National Institute Of Mental Health of the National Institutes of Health under Award Number U19MH114830. The content is solely the responsibility of the authors and does not necessarily represent the official views of the National Institutes of Health.

Attachments



MB0073_Fragment_Anal

...

30KB

