

Aug 23, 2024

Version 2

Freezing Adherent Cell Lines V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.5qpvokxq9l4o/v2>

Carolina Lopez¹

¹Washington University



Carolina Lopez

Washington University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.5qpvokxq9l4o/v2>

Protocol Citation: Carolina Lopez 2024. Freezing Adherent Cell Lines. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.5qpvokxq9l4o/v2> Version created by **Sydney Faber**

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 22, 2024

Last Modified: August 23, 2024

Protocol Integer ID: 106113

Keywords: freezing adherent cell line, adherent cell lines this protocol, adherent cell line, mdck cell, adherent cell, llcmk2 cell, examples of these cell, a549 cell, cell, freezing

Abstract

This protocol describes how to freeze Adherent cells. Examples of these cells are: A549 cells, LLCMK2 cells, and MDCK cells.

Materials

FREEZING MEDIUM:

90% of TCM

10% of DMSO

TISSUE CULTURE MEDIUM (TCM): Filter through 0.2 μ m filter

	Component	Amount	Conc. Supp.	Product information
	DMEM	500 mL		Gibco Cat. # 11965092 (#11965118-cs)
	Gentamicin	500 μ L	50 μ g/mL	Gibco Cat #15750060 (#15750078-pk) – 50 mg/mL
	Sodium Pyruvate	5.0 mL	1mM	Corning Cat #25-000-CI – 100mM
	L-Glutamine	5.5 mL	2 mM	Sigma Aldrich Cat #G7513 – 200mM
	FBS	50 mL	10%	

DMSO (Dimethyl sulfoxide): Sigma-Aldrich Cat# D2650

Freezing Adherent cells

1

*Try to freeze cells at the lowest passage possible and confirm the cells are negative for mycoplasma before freezing.

1. Wash flask 2x with sterile 1X PBS.
2. Add 2mL of trypsin/T75. Incubate at 37°C for 2-3 mins or until cells are detached from the flask.
3. Add 8mL of cell culture media and pipette up and down. Transfer all the media to a 15mL tube.
4. Centrifuge at 1200 rpm for 5 min.
5. Discard supernatant and resuspend cell pellet in 1mL of cell culture media. Add 9mL of media and pipette up and down to homogenize.
6. Count the cells. You will need your cell concentration/# for step 6.
7. Spin the tube containing the cells in media again at 1200 rpm for 5 min.
8. Meanwhile, prepare the freezing medium (see materials for recipe). Prepare the amount of freezing medium needed to dilute the cell pellet so that you have a final concentration of 1×10^6 cells/mL.
9. Discard the supernatant and resuspend the pellet in 1mL of freezing medium. Add the remaining volume of the freezing medium. Make sure you homogenize the solution well.
10. Add 1mL of freezing medium containing 1×10^6 cells/mL to cryotubes (caps screw from the outside). Label all the tubes with the following information:
 - **Cell Name**
 - **Generation/Passage**
 - **Date**
 - **Name**
 - **Cell concentration (if there is space)**
11. Transfer the cells to a Mr. Frosty Freezing container containing isopropanol and then transfer the container to the -80°C for 1 day. After 1 day, transfer the tubes to the liquid nitrogen. You can keep some vials in the -80°C for a couple of months, however, the -80°C is not for long-term storage.