

Nov 07, 2025

GCase activity assay

DOI

<https://dx.doi.org/10.17504/protocols.io.3byl46yb8go5/v1>

Sara Lucas Del Pozo^{1,2}, Giuseppe Uras^{1,2,3}, Federico Fierli^{1,2}, Veronica Lentini^{1,3}, Sofia Koletsi^{1,2}, Carlos Lazaro-Hernandez^{1,4}, Kai-Yin Chau^{1,2}, Derralynn A. Hughes⁵, Anthony H.V. Schapira⁵

¹Department of Clinical and Movement Neurosciences, UCL Queen Square Institute of Neurology, London, UK;

²Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD, USA;

³Department of Biomedical Science, University of Sassari, Viale San Pietro, Sassari 07100, Italy;

⁴Neurology Department, Vall d'Hebron University Hospital, Barcelona, Spain;

⁵Lysosomal Storage Disorders Unit, Royal Free Hospital NHS Foundation Trust and University College London, London, UK



Sofia Koletsi

University College London, University of London

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.3byl46yb8go5/v1>

Protocol Citation: Sara Lucas Del Pozo, Giuseppe Uras, Federico Fierli, Veronica Lentini, Sofia Koletsi, Carlos Lazaro-Hernandez, Kai-Yin Chau, Derralynn A. Hughes, Anthony H.V. Schapira 2025. GCase activity assay. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.3byl46yb8go5/v1>

Manuscript citation:

<https://doi.org/10.1016/j.nbd.2025.107053>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 02, 2023

Last Modified: November 07, 2025

Protocol Integer ID: 231665

Keywords: ASAPCRN, gcse protein, gcse activity, crucial part of the assay, acidic enzyme assay buffer, glucopyranoside, enzyme, lysosome, assay, enzyme in an active confirmation, utilised bile salt, bile salt, low ph, blue fluorogenic substrate, utilising blue fluorogenic substrate, taurocholate

Funders Acknowledgements:

Aligning Science Across Parkinson's Disease

Grant ID: ASAP-000420

Abstract

This protocol describes a commonly used approach to measure GCase activity by utilising blue fluorogenic substrate 4-methylumbelliferyl- β -D-glucopyranoside (4-MUG). In order to simulate the lysosome's low pH, this method uses GCase protein that has been isolated from a cellular model and diluted in an acidic enzyme assay buffer. The inclusion of a lipid or detergent to keep the enzyme in an active confirmation is a crucial part of the assay. Taurocholate, a regularly utilised bile salt, is used in this protocol.

Materials

Assay Buffer:

A	B
Citric acid	50 mM
K ₂ HPO ₄	176 mM
Sodium Taurocholate (Sigma Aldrich REF 86339-5G)	10 mM
Tween 20 (Sigma Aldrich REF P9416-50ML)	0.01%
dH ₂ O	

NOTE: The pH should adjusted down to make it closer to pH 5.8

Complete Lysis Buffer:

A	B
Cold Assay Buffer	10 ml
20% Triton X (Sigma Aldrich PCode: 1002733102)	100 µl
Protease Cocktail Inhibitor (Thermob Scientific REF A32965)	100 µl

STOP solution:

NaOH	0.5 M
Glycine	0.5 M

NOTE: The solution need to be adjusted to 1000ml with distilled H₂O (ie addition of 650ml) and the solution also needs to be adjusted to a pH of 10 approximately. The solution could be stored at room temperature.

Substrate:

[M] 1 millimolar (mM) 4-methylumbelliferyl β-D glucopyranosidase (4-MUG) (Sigma Aldrich [®] REF M3633-250MG)



Note

Dilute 4MUG in Assay Buffer.

⊗ Sodium taurocholate hydrate **Merck MilliporeSigma (Sigma-Aldrich) Catalog #86339**

⊗ TWEEN® 20 **Merck MilliporeSigma (Sigma-Aldrich) Catalog #P9416**

⊗ Pierce™ Protease Inhibitor Tablets, EDTA-free **Thermo Fisher Catalog #A32965**

⊗ 4-Methylumbelliferyl β -D-glucopyranoside **Merck MilliporeSigma (Sigma-Aldrich) Catalog #M3633**

⊗ Phosphate buffered saline (PBS) without Ca/Mg **Thermo Fisher Scientific Catalog #14190144**



GCase activity assay

1d 1h 20m

- 1 Seed the total of 3×10^4 cells in each well of a 96-well plate (Corning® Assay Plate REF 3340) and left to incubate for 24:00:00 . 1d
- 2 Remove the cell medium and wash each well 2 times with 1X PBS (Gibco® REF 14190-144) at Room temperature
- 3 Lyse the cells using 50 μL cold Complete Lysis Buffer for 00:10:00 at 4 °C in a shaker. 10m
- 4 Afterwards, wash the cells two times in 200 μL of sterile 1X PBS.
- 5 Subsequently, supplement 50 μL of Assay Buffer with 1 millimolar (mM) 4-MUG that are added to each well and left the samples to incubate Overnight at Room temperature and protected from light. 10m
- 6 Stop the reaction by adding 100 μL of STOP solution to each well, and then is left to incubate for 01:00:00 at Room temperature . 1h
- 7 Analyse the plates using CytationTM® plate reader using endpoint protocol with wavelength ex/em set up at 360/460nm.
- 8 Normalise the obtained data to control and expresses as a fraction.