

Jun 28, 2022

Generating a low-copy overexpression cell line using PhiC31 integration

DOI

<https://dx.doi.org/10.17504/protocols.io.36wgg71kxvk5/v1>

Goran Tomic¹

¹The Francis Crick Institute

GT_Protocols



Goran Tomic

Relation Therapeutics

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.36wgg71kxvk5/v1>

Protocol Citation: Goran Tomic 2022. Generating a low-copy overexpression cell line using PhiC31 integration. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.36wgg71kxvk5/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 28, 2022

Last Modified: June 28, 2022

Protocol Integer ID: 65542

Keywords: phiC31, overexpression, attb, phic31 integrase, phic31 integration, phic31 integrase vector system, pcmvint plasmid, using phic31 integration, px28 plasmid, addgene plasmid, pgk promoter, standard cl1 tissue culture room, copy overexpression cell line, plasmid, gene, attb construct, bgh pa sequence, clone

Abstract

This is an in-house adaptation of the commercially available Jump-In system by Thermo. It generates clones with a lower overexpression level than those generated by lentiviral delivery and can be carried out in a standard CL1 tissue culture room. The gene of interest in the attB construct is assembled in-house and is driven by the PGK promoter and contains the BGH pA sequence (assembled from pX28 plasmid (Addgene Plasmid #46850), which had the required sequences). pCMVInt plasmid that expresses the PhiC31 integrase was purchased from Addgene (Plasmid #18935).

This protocol is made combining the protocol for Jump-In system (ThermoFisher) and phiC31 Integrase Vector System (System Biosciences).

Attachments



PhiC31_schematic cop...

140KB

Materials

 pCMVInt **addgene Catalog #18935**

Before start

Perform a quick Blasticidin curve to determine the optimal concentration that kills the cells (0-20 ug/mL). Use 24 well plates and try, 0, 2.5, 5, 10, 15, 20 ug/mL Bsd.



- 1 Day 1: Use a 6-well plate (0.15×10^6 cells per well). Incubate for 24-48 hours. Transfect if the well is 70-90% confluent.
- 2 Day 2: Make DNA-Lipofectamine™ 3000 complexes in serum-free medium such as Opti-MEM™ Reduced Serum Medium.

For a 6 well plate: 2.5 ug of DNA + 125 uL Opti-MEM + 5 uL P3000 reagent (2 uL/ug DNA). Mix PhiC31 integrase plasmid in a 50:1 ratio (w/w) over the donor i.e. 2500 ng of PhiC31 + 50 ng of GOI/attB plasmid

125 uL OptiMEM + 3.75 uL Lipofectamine 3000

Mix 1:1 and incubate for 15 min at RT

Add 250 uL per well to cells in culture medium

- 3 Day 3: Split the cells into 10 cm plates at 1:10 and 1:20 ratio and add blasticidin. Observe daily and change blasticidin media every 3-4 days until colonies appear. Validate the expression of the GOI using qPCR, WB, etc.

N.B. To increase the probability of getting single-copy integration of the targeting expression construct, use the amount of vector that results in less than five drug-resistant colonies in the absence of integrase. Using that amount in the presence of integrase should result in >15 colonies.

It is good to have an empty vector control (to generate Bsd-resistant cells, but without GOI).

Important! We observed that in some clones the expression of the GOI can reduce over time. Best to freeze an early passage of the selected single clones after validation, and check the expression of the GOI before any important experiments.