

Jul 21, 2025

Generation of CRISPR/Cas9 edited mammalian cells with small epitope tags

DOI

<https://dx.doi.org/10.17504/protocols.io.j8nlkrxo5v5r/v1>

Hely Rodriguez Cruz¹, Michael Hanna¹, Pietro De Camilli¹

¹Yale University



Michael G Hanna

Yale University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.j8nlkrxo5v5r/v1>

Protocol Citation: Hely Rodriguez Cruz, Michael Hanna, Pietro De Camilli 2025. Generation of CRISPR/Cas9 edited mammalian cells with small epitope tags. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.j8nlkrxo5v5r/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 05, 2025



Last Modified: July 21, 2025

Protocol Integer ID: 219601

Keywords: ASAPCRN, generation of crispr, mammalian cells with small epitope tag, edited mammalian cell, immortalized mammalian cell culture line, small epitope tag, crispr, mammalian cell culture line, gene, cell, cas9

Abstract

This protocol explains how to endogenously edit immortalized mammalian cell culture lines to engineer an in-frame small epitope tag into a gene of interest.

Materials

- Alt-R™ CRISPR-Cas9 crRNA (synthesized by IDT)
- Alt-R™ CRISPR-Cas9 tracrRNA (IDT, Cat# 1075934)
- Alt-R™ S.p. HiFi Cas9 Nuclease V3 (IDT, Cat# 1081060)
- Nuclease-free duplex buffer (IDT, Cat# 11-01-03-01, included with Alt-R CRISPR-Cas9 tracrRNA)
- Single stranded repair oligo (IDT)
- PCR tube
- Thermal cycler
- 1xPBS
- Trypsin-EDTA (0.25%), phenol red
- 15 mL conical tube
- 1.5mL microcentrifuge tubes
- Microcentrifuge
- Lonza 4D Nucleofector System Core Unit with X Unit
- Lonza nucleofection solution
- Lonza nucleofector cuvette
- DMEM complete medium (10%FBS, 1% Pen/Strep, 0.2% Plasmocin)
- 6-well cell culture plate
- HDR enhancer (IDT, Cat# 10007910)



Assemble the gRNA

15m

- 1 Resuspend Alt-R CRISPR-Cas9 crRNA and Alt-R CRISPR-Cas9 tracrRNA in nuclease-free duplex buffer at 200 μ M
- 2 Mix 5 μ l of 200 μ M crRNA and 5 μ l of 200 μ M tracrRNA together in a PCR tube
- 2.1 Heat to 95°C for 5 minutes in a thermal cycler
- 2.2 Let cool to RT

5m

10m

Assemble the RNP

30m

- 3 In a microcentrifuge tube, mix 1.7 μ l (104 pmol) of Alt-R S.p. Cas9 nuclease together with 1.2 μ l (120 pmol) Alt-R CRISPR-Cas9 crRNA/tracrRNA duplex and 2.1 μ l of 1xPBS solution.
- 3.1 Incubate for 30 minutes at RT

30m

Prepare repair oligo

- 4 Resuspend repair oligo in 1xPBS at a concentration of 100 μ M (pmol/ μ l).

Prepare cell & electroporation

10m

- 5 Lift A549 cells from dish using trypsin digest and collect in a 15 mL conical
- 6 Count cells and pellet ~800,000 cells in a microcentrifuge tube at 300g for 3 minutes.
- 7 Resuspend cell pellet in 100 μ L desired Lonza nucleofection solution

5m

5m



- 8 Add 2 μ L of repair oligo (100 μ M) to 5 μ L of pre-assembled RNP complex from Step 3
- 9 Add RNP complex and repair oligo to resuspended cells in nucleofector cuvette
- 10 Electroporate cells according to Lonza nucleofection kit specifications

Post electroporation

- 11 Remove cells from nucleofector cuvette and directly add to 2mL warmed complete culture media in one well of a 6-well plate
- 12 Add 3.4 μ L of HDR enhancer to the 2 mL culture (optional)