



Nov 20, 2025

Genome Edited (KO) Cell Line Generation – PCR Validation



Forked from [Genome Edited \(KO\) Cell Line Generation –](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.5jyl88wj6l2w/v1>

Amanda Bentley-DeSousa^{1,2,3,4,5}, Devin Clegg^{1,2,3,4,5}, Shawn Ferguson^{1,2,3,4,5,6}

¹Department of Cell Biology, Yale University School of Medicine, New Haven, Connecticut 06510, USA;

²Neuroscience, Yale University School of Medicine, New Haven, Connecticut 06510, USA;

³Program in Cellular Neuroscience, Neurodegeneration and Repair;

⁴Wu Tsai Institute Yale University School of Medicine, New Haven, Connecticut 06510, USA;

⁵Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD, 20815, USA;

⁶Kavli Institute for Neuroscience, Yale University School of Medicine, New Haven, Connecticut 06510, USA



Devin Clegg

Yale University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.5jyl88wj6l2w/v1>



Protocol Citation: Amanda Bentley-DeSousa, Devin Clegg, Shawn Ferguson 2025. Genome Edited (KO) Cell Line Generation – PCR Validation. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.5jyl88wj6l2w/v1>

Manuscript citation:

<https://doi.org/10.1101/2023.10.31.564602>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 01, 2025

Last Modified: November 20, 2025

Protocol Integer ID: 228804

Keywords: RAW 264.7, KO, Synthego, ko cell lines in raw, ko cell line, cell line generation, cell lines in raw, genome edited, genome, cell, line, ko, raw, pcr validate ko cell line, pcr validate ko cell lines in raw, pcr validation, pcr validation this protocol details step

Funders Acknowledgements:

Aligning Science Across Parkinson's Disease

Grant ID: ASAP-025173

Abstract

This protocol details steps taken to create and PCR validate KO cell lines in RAW 264.7 cells.

Materials

DMEM (Thermo Fisher Scientific, 11965-092)

FBS (Thermo Fisher Scientific, 16140-071)

Penicillin/Streptomycin (Thermo Fisher Scientific, 15140122)

CellStripper (Corning, 356230)

Synthego CRISPR Gene Knockout V2 Mouse Kits

Lipofectamine CRISPRiMAX (ThermoFisher Scientific, CMAX00003)

Cas9 (Synthego CRISPR Gene Knockout V2)

Opti-Mem (Thermo Fisher Scientific, 31985062)



Day 0

- 1 Plate 2.5×10^5 RAW 264.7 cells per well in a 6-well dish.

Day 1

- 2 Transfect cells using Lipofectamine CRISPRiMAX, gene specific sgRNAs and recombinant Cas9 (Synthego CRISPR Gene Knockout V2).

2.1 CRISPRiMAX protocol

In Tube 1 combine  62.5 μL Opti-MEM,  3.25 μL sgRNA (from 3 uM stock),  2.5 μL Cas9 (from 3 uM stock), and  2.5 μL Cas9+ then incubate for  00:05:00 . In Tube 2 combine  62.5 μL Opti-MEM and  3.75 μL CRISPRMAX then incubate for  00:05:00 . Combine tube 1 and tube 2. Incubate for  00:20:00 . Add drop-wise to 1 well containing  1.8 mL of media.

Day 3

- 3 Change the media.

Day 4

- 4 Plate single cells into 96-well dishes to obtain clonal cell lines. After subculturing, confirm cell lines via immunoblotting

- 4.1 Expand clonal cell lines to 24-well dish
(Design primers flanking Cas9 cut sites)

Harvesting DNA using Quick Extract

8m 20s

- 5 Dissociate cells from 24-well dish

- 5.1 Centrifuge Cells  1000 rpm, 00:03:00



5.2 Discard supernatant

6 Add  200 μL of Quick Extract to cells

7 Vortex sample for  00:00:10

10s

8 Incubate sample at  65 °C for  00:06:00

6m

9 Vortex sample for  00:00:10

10s

10 Incubate sample at  98 °C for  00:02:00

2m

PCR Validation

33m

11 Run PCR using designed primers

11.1 PCR Reaction Setup

GoTaq® Green Master Mix, 2X  12.5 μL

upstream primer, 10 μM  1 μL

downstream primer, 10 μM  1 μL

DNA template  3 μL

Nuclease-free water  7.5 μL

Total  25 μL

12 Thermocycler:

3m

Initial denaturation  98 °C  00:02:00

denaturation  98 °C  00:00:30

Annealing  00:00:30 Temp. is determined by primers



Extension  72 °C 1 minute per kilobase

30 Cycles

Hold  4 °C

13 Run PCR reactions on 1% agarose gel

100v for  00:30:00

30m