

Dec 18, 2019

GENOTYPING OF iPSCS WITH GENE INSERTIONS (Support Protocol 1)

 In 1 collection

DOI

dx.doi.org/10.17504/protocols.io.5u9g6z6

Michael S. Fernandopulle¹, Ryan Prestil¹, Christopher Grunseich¹, Chao Wang², Li Gan², Michael E. Ward¹

¹National Institute of Neurological Disorders and Stroke, National Institutes of Health, Bethesda, Maryland;

²Gladstone Institute of Neurological Disease, Gladstone Institutes, San Francisco, California

Neurodegeneration Method Development Community
Tech. support email: ndcn-help@chanzuckerberg.com



Julia Rossmanith

protocols.io

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.5u9g6z6

External link: <https://doi.org/10.1002/cpcb.51>

Protocol Citation: Michael S. Fernandopulle, Ryan Prestil, Christopher Grunseich, Chao Wang, Li Gan, Michael E. Ward 2019. GENOTYPING OF iPSCS WITH GENE INSERTIONS (Support Protocol 1). **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.5u9g6z6>

Manuscript citation:

Fernandopulle, M. S., Prestil, R., Grunseich, C., Wang, C., Gan, L., & Ward, M. E. (2018). Transcription-factor mediated differentiation of human iPSCs into neurons. Current Protocols in Cell Biology, e51. doi:<https://doi.org/10.1002/cpcb.51>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: July 25, 2019

iPSCs

Last Modified: December 18, 2019

Protocol Integer ID: 26241

Keywords: i3LMN, i3Neurons, iPSC, iPSC-derived neurons, transcription factor-mediated differentiation

Abstract

This protocol will use genomic DNA isolated from the purified iPSC clones with CLYBL or AAVS1 gene insertion to determine if integration of the transgene has occurred correctly, in a heterozygous or homozygous fashion, and if the floxed selection genes are present. Primer sequences, amplicon sizes, PCR mix composition, and thermal conditions are included in the steps.

Attachments



[fernandopulle2018.pdf...](#)

1.7MB

Materials

- Primers (see recipe in Reagents and Solutions)
- PCR reagents (e.g., Platinum SuperFi PCR Master Mix, Invitrogen, cat. no. 12358250)

 Platinum™ SuperFi™ PCR Master Mix **Thermo Fisher Scientific Catalog #12358010**

- Genomic DNA
- Additional reagents and equipment for PCR (Kramer & Coen, 2001) and agarose gel electrophoresis (Voytas, 2001)

Safety warnings

 Please see SDS (Safety Data Sheet) for hazards and safety warnings.

- 1 Set up PCR reaction (also see Kramer & Coen, 2001).
For each 10 µl reaction, use the primer pairs recommended in **Reagents and Solutions** for the given safe-harbor site and donor construct chosen:

1 µL	10 micromolar (µM)	Primer 1
1 µL	10 micromolar (µM)	Primer 2
1 µL	purified genomic DNA from iPSC clone	
2 µL	H ₂ O	
5 µL	2 × PCR Master Mix.	

- 2 Run PCR reactions

95 °C	3min
95 °C	30 sec
64 °C for CLYBL primers, 58 °C for AAVS1 primers	30 sec
72 °C	1min
Repeat (to step 2) 34 times	
72 °C	5min
Hold at 12 °C	Indefinitely

- 3 Run PCR products on 1 % agarose gel (Voytas, 2001) with interpretation given above for each primer pair.

Note

If performing the PCR on samples following Cre excision, increasing the amount of time during the extension step may be necessary to detect the larger un-excised template if a heterogeneous culture is obtained.