

Nov 12, 2022

Version 3

Germination and seedling establishment for deep-flow hydroponics: The benefit of slant boards V.3

DOI

<https://dx.doi.org/10.17504/protocols.io.5jyl89b86v2w/v3>

noah.langenfeld¹, Bruce Bugbee¹

¹Utah State University

USU Crop Physiology La...


PLOS ONE Lab Protocols

1 more workspace



Noah Langenfeld

Utah State University

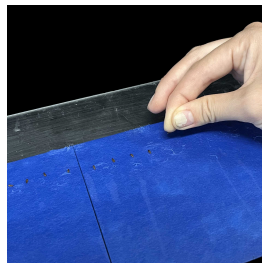

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.5jyl89b86v2w/v3>

External link: <https://doi.org/10.1371/journal.pone.0275710>



Protocol Citation: noah.langenfeld , Bruce Bugbee 2022. Germination and seedling establishment for deep-flow hydroponics: The benefit of slant boards. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.5jyl89b86v2w/v3> Version created by **Noah Langenfeld**

Manuscript citation:

Langenfeld NJ, Bugbee B (2022) Germination and seedling establishment for hydroponics: The benefit of slant boards. PLOS ONE 17(10): e0275710.

<https://doi.org/10.1371/journal.pone.0275710>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 03, 2022

Last Modified: November 12, 2022

Protocol Integer ID: 72221

Keywords: slant boards germination, benefit of slant boards germination, germination paper, seedlings with long root, polycarbonate sheets with germination paper, procedures for rapid germination, germination, rapid germination, seedling, seedling establishment, hydroponic, flow hydroponic, germination of seed, hydroponic system, slant board construction, slant board, seed, nutrient solution, board, transplanting, thin paper before vertical placement, damaging root, long root, polycarbonate sheet, exposure of the primary root

Abstract

Germination and seedling establishment for transplanting into hydroponics often uses porous substrates, but fine roots grow into these substrates, and they cannot be removed without damaging roots. Seedlings transplanted without removal of substrates can cause interactions with solution chemistry or addition of particulates to the nutrient solution. Germination of seeds on slant boards is clean, uniform, and reduces the time to transplanting. Slant boards facilitate long roots, which maximize exposure of the primary root to the nutrient solution after transplanting. The "boards" are made from thin acrylic or polycarbonate sheets with germination paper on top. Seeds are held in place by covering with thin paper before vertical placement of the boards in a container. Four to twelve days later, the seedlings with long roots can be removed from the paper without damage and transplanted into the hydroponic system. Here we describe slant board construction and procedures for rapid germination and transplanting.

(The last step contains a supplemental video with extra context and tips, as part of the protocols.io Spotlight series, featuring conversations with protocol authors.)



Materials

- Seeds
- 2 mm thick acrylic board
- Germination blotter paper
- Single-ply cellulose paper (Kimwipes)
- Acrylic holder
- Exterior container (plastic storage tote)
- Water

Seeding

- 1 Clean a rigid plastic slant board and lay flat on a clean surface.
- 2 Cut a piece of germination paper 10 cm in height and the same length as the slant board.
- 3 Place germination paper on top of slant board.
- 4 Place seeds on top of germination paper (about 9 cm from the bottom of the board and 1 cm from the top of the germination paper) in a line parallel to the top and bottom of the slant board. Seeds should be about 5 mm apart.
- 5 Place a single layer of thin absorbent paper (Kimwipe) over the seeds.
- 6 Use a squirt bottle to gently wet the entire germination board and thin absorbent paper with water to lock seeds and paper in place on the board. Do not use nutrient solution.
- 7 Transfer the slant board into the exterior container. The angle of the board should be between 70° and 80°.
- 8 Fill the exterior container with purified water with calcium until the water level is about 2 cm above the bottom of the germination paper.

Monitoring and transplanting

- 9 After at least one cotyledon leaf has appeared, gently peel back the thin absorbent paper to expose the leaves. The thin absorbent paper should remain covering the roots.
- 10 After the roots have grown long enough to reach the solution, remove the seedlings from the paper. Fingers or a tweezers can be used to remove the seedlings. Avoid excessive pressure to prevent root damage.
- 11 Place the seedling in a neoprene collar. Ensure the roots are not pinched in the collar. The cotyledons can be used to support the seedling and stop it from falling into the solution.



- 12 Place the neoprene collar with the seedling into the main culture tank. Ensure to roots can reach the nutrient solution to avoid desiccation.
- 13 (The following video contains extra context and tips, as part of the protocols.io Spotlight series, featuring conversations with protocol authors.)

https://www.youtube.com/embed/A4P_Pjg-Hvs