

Sep 20, 2019

Gibson Assembly

DOI

<https://dx.doi.org/10.17504/protocols.io.7h5hj86>

宏亮¹

¹Northeast Forest University

2019 iGEM NEFU_China

Tech. support email: shengyiyanwork@gmail.com



宏亮 董

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.7h5hj86>

Protocol Citation: 宏亮 2019. Gibson Assembly. protocols.io <https://dx.doi.org/10.17504/protocols.io.7h5hj86>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: September 19, 2019

Last Modified: September 20, 2019

Protocol Integer ID: 27933

Keywords: Gibson Assembly, gibson assembly this protocol, gibson assembly, different kit, assembly

Abstract

This protocol includes two different kits. Choose one to use.

Materials

The old kit: Trelief™ SoSoo Cloning Kit Ver.2

The new kit: ClonExpress® II One Step Cloning Kit

Safety warnings

 Please wear gloves for the experiment.



The old kit

- 1 Set up a small box with ice, put DNA and Gibson Mix into it before going into the metal bath or the Bio-rad S1000TM thermocycler cycler .

- 2 Add the following reagent to a PCR tube (4 μ l).

Gibson Mix	2 μ l
insert	X μ l
vector	Y μ l

$X+Y=2$ ul , $X : Y=3 : 1\sim 5 : 1$.

Temperature : Time= 50°C : 15-30 min

The new kit

- 3 Set up a small box with ice, put DNA and Gibson enzyme and Gibson buffer into it before going into the metal bath or the Bio-rad S1000TM thermocycler cycler .

- 4 Add the following reagent to a PCR tube (10 μ l).

Gibson enzyme	1 μ l
Gibson buffer	2 μ l
insert	X μ l
vector	Y ul

$X=0.02\times$ vector length /vector concentration

$Y=0.02\times$ insert length /insert concentration

Temperature : Time= 37°C : 30-60 min