

Jun 02, 2020

## Glycine-HCl Buffer

 In 1 collection

DOI

[dx.doi.org/10.17504/protocols.io.bfybjpsn](https://dx.doi.org/10.17504/protocols.io.bfybjpsn)

Neillier Junior<sup>1</sup>

<sup>1</sup>University of Manitoba



**Neillier Junior**

University of Manitoba, Universidade Federal de Viçosa



OPEN  ACCESS



DOI: [dx.doi.org/10.17504/protocols.io.bfybjpsn](https://dx.doi.org/10.17504/protocols.io.bfybjpsn)

**Protocol Citation:** Neillier Junior 2020. Glycine-HCl Buffer. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.bfybjpsn>

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** May 04, 2020

**Last Modified:** June 02, 2020

**Protocol Integer ID:** 36579

**Keywords:** pH, Henderson-Hasselback, Biochemistry, Molecular Biology,



## Abstract

A buffer solution has the function of resisting changes in pH even when adding powerful acids or bases. However, in the physiological environment the buffered system also provides cofactors for enzymatic reactions, critical salts and even essential nutrients for cells and tissues. Therefore, when trying to reproduce biological conditions in vitro, we must make the appropriate choice of the buffer. After all, it will provide the appropriate medium in which reactions will occur.

## Materials

- Deionized Water
- pH Meter (sensitive)
- Glycine
- Hydrochloric Acid

## Safety warnings

 Wear personal protective equipment: gloves, lab coat and mask.

## Before start

Organize your workspace.


Make sure all solutions and equipment are available.



## Glycine-HCl Buffer

1

### Note

pH range:  2.2 to  3.6

(a) 0.1 M Glycine: 7.5 g L<sup>-1</sup> (M.W.: 75.0 g mol<sup>-1</sup>)

(b) 0.1 M Hydrochloric acid

Mix  50 mL glycine and indicated volume of hydrochloric acid.

<b>mL of HCl</b>	44.0	32.4	24.2	16.8	11.4	8.2	6.4	3.6
<b>pH</b>	2.2	2.4	2.6	2.8	3.0	3.2	3.4	3.6

2

Mix and adjust the final volume to  100 mL with deionized water.

3

Adjust the final pH using a sensitive pH meter