


Aug 07, 2023

Version 1

🌐 Grid patterning protocol for Cryo-FIB/ET workflow V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.e6nvwdjewlmk/v3>

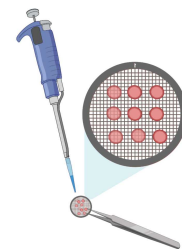

Farhaz Shaikh¹, Josh Hutchings¹, Tamar Basiashvili², Elizabeth Villa^{1,3}

¹Biological Sciences, UCSD; ²Chemistry and Biochemistry, UCSD; ³HHMI



Tamar Basiashvili

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.e6nvwdjewlmk/v3>

Protocol Citation: Farhaz Shaikh, Josh Hutchings, Tamar Basiashvili, Elizabeth Villa 2023. Grid patterning protocol for Cryo-FIB/ET workflow. [protocols.io https://dx.doi.org/10.17504/protocols.io.e6nvwdjewlmk/v3](https://dx.doi.org/10.17504/protocols.io.e6nvwdjewlmk/v3)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](https://creativecommons.org/licenses/by/4.0/), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: August 07, 2023

Last Modified: August 07, 2023

Protocol Integer ID: 86062

Keywords: Cryo-FIB/ET, Cryo grid preparation, Cryo-FIB milling, In situ structural biology, Grid, Photo-Micropatterning , electron microscopy grid preparation, microscopy grid preparation, architecture of organelle, focused ion beam milling, electron tomography, protocol for cryo, organelle, cellular deposition, et workflow cryo, fib milling, micropatterns on the grid, cryo, accessibility for cryo, successful cryo

Funders Acknowledgements:

Aligning Science Across Parkinson's

Grant ID: ASAP-000519

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to protocols.io is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with protocols.io, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

Cryo-Electron Tomography combined with Cryo-Focused Ion Beam milling provides a novel way to study the structure of proteins and the architecture of organelles in situ. Electron microscopy grid preparation is a key step toward a successful Cryo-FIB/ET workflow execution. Here we provide an optimized protocol to generate photo-micropatterns on the grids to enhance cellular deposition and accessibility for Cryo-FIB milling.



Materials

- Use Gold AU 200 R1/4 Grids for mammalian cells.
- Pelco Plasma cleaner.
- A 10ul PLPP aliquot in ice
- Tweezers
- 1.5 ml centrifuge tubes that each hold 1 mL of DPBS buffer
- Two P10 pipette with tips
- Blotting paper



1. Passivation

- 1 Place a PDMS stencil (14 mm with 2×2 4 mm wells) into a 35 mm confocal dish (MatTek P35G-1.5-20-C).
- 2 Glow-discharge both sides of the grids—standard settings suffice (20 mA, 1 min hold, 3-sec hold). Discharge gold side up first. Then discharge carbon film side up.
- 3 While grids are discharging you can place 1 uL of water in the center of each of the PDMS stencil sections. This is intended to keep grids in place.
- 4 Get PLL-PEG (100 ul of 1ug/ml) from the -20 freezer. Dilute two fold to working concentration with 100 ul of HEPES buffer (10 mM, pH 7.4).
- 5 After grids are discharged, add the grid to the center of each of the PDMS wells. Make sure the carbon film side is up.
- 6 Apply 10 uL of the PLL-PEG solution onto each grid. Start a timer for 1 hour. After 1 hour, you can begin patterning.  01:00:00


1h

2. Preparing Microscope

- 7 Switch on a light source, microscope, microscope light, Primo (button on the back and key to the right), PC, then PCO camera – in that order.
- 8 Switch objective to 20x lens with correction collar (S plan Fluor20x), set filter cube #6, and set condenser turret to #7.
- 9 Open Micromanager. Select Leonardo from plug-ins menu.
- 10 We recommend using a glass slide with yellow highlighter to verify and focus the Primo laser. The laser can be toggled on/off in Leonardo when autoscale is on.

3. Patterning

- 11 Remove PLL-PEG solution from the grid.
- 12 Wash with 10 uL of DPBS four times, try not to touch the grid. During each wash, pipette up and down four times.
- 13 Blot extra liquid around grids.
- 14 Add 1 uL of PLPP onto the grid.
- 15 Place the dish onto the microscope stage, drive stage to grid center and focus on the foil giving the sharpest view of holes in the foil mesh.
- 16 Set contrast by pressing and unpressing autoscale.
- 17 Automatic Patterning: Load TEM wizard and specify settings.
 - Select "Add Grid"
 - Select "Automatic execution"
 - Select "Lock"
 - Select "Rotate"
 - Select "Finish"

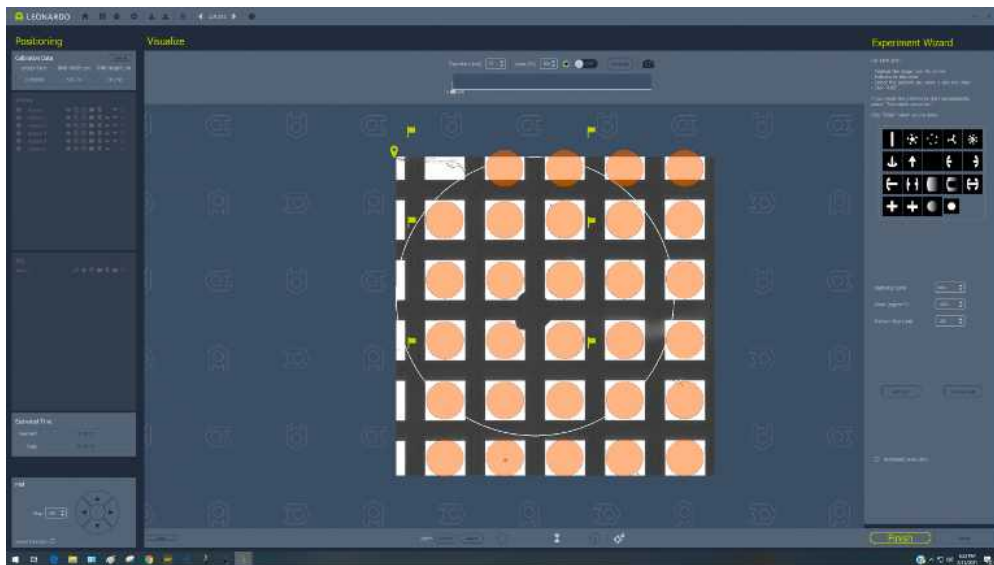



- 18 Manual Patterning:
 - Select "MicroPatterning"
 - Select desired pattern
 - Set desired ration and angle
 - Set dose to 1000
 - Select "Lock"
 - Select "Play"

4. Finishing up

- 19 Wash the grids with 10ul of DPBS four times, try not to touch the grid. During each wash, pipette up and down four times.
- 20 Leave 10ul of DPBS on the grids.
- 21 Repeat the patterning steps for each grid you have.

Protocol references

Toro-Nahuelpan, M., Zagoriy, I., Senger, F. *et al.* Tailoring cryo-electron microscopy grids by photo-micropatterning for in-cell structural studies. *Nat Methods* **17**, 50–54 (2020). <https://doi.org/10.1038/s41592-019-0630-5>