



Oct 26, 2018

Growth media for *V. natriegens*

DOI

dx.doi.org/10.17504/protocols.io.u3ceyiw

Josef Hoff¹

¹Phillips University



Josef Hoff

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.u3ceyiw>

Protocol Citation: Josef Hoff 2018. Growth media for *V. natriegens*. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.u3ceyiw>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 26, 2018



Last Modified: October 26, 2018

Protocol Integer ID: 17220

Keywords: additional v2 salt, additional with additional v2 salt, v2 salt, natriegens media, brain heart infusion broth, growth media, vibrio natriegen, nature methods volume, salt, host for molecular biology, bhiv2, mgcl2, mm mgcl2, mm nacl



Abstract

Media for *V. natriegens*

LB2:

Standard LB media (10.0 g/L Tryptone, 5.0 g/L Yeast Extract, 10.0 g/L NaCl) with additional 1.5% NaCl

LBv2 Salts:

Standard LB media (10.0 g/L Tryptone, 5.0 g/L Yeast Extract, 10.0 g/L NaCl) with additional with additional v2 salts (204 mM NaCl, 4.2 mM KCl, and 23.14 mM MgCl₂)

BHlv2:

Brain heart infusion Broth (37 g/L Teknova Brain Heart Infusion Broth Dry Media (Cat. No. B9500) with additional v2 salts (204 mM NaCl, 4.2 mM KCl, and 23.14 mM MgCl₂)

note: BHI can not be autoclaved together with the salts, they have to be autoclave separate and later on added.



Matthew T Weinstock, Eric D Heseck, Christopher M Wilson, Daniel G Gibson

Vibrio natriegens as a fast-growing host for molecular biology

Nature Methods volume 13, pages 849–851 (2016)

Attachments



Media for V.pdf

35KB

Troubleshooting

