

Nov 12, 2023

Version 1

GST fusion protein production V.1

DOI

dx.doi.org/10.17504/protocols.io.4r3l22y14l1y/v1

Leonardo A Parra-Rivas¹

¹University of California, San Diego



Leonardo A Parra-Rivas

University of California, San Diego

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l22y14l1y/v1>

Protocol Citation: Leonardo A Parra-Rivas 2023. GST fusion protein production. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.4r3l22y14l1y/v1>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 12, 2023



Last Modified: November 12, 2023

Protocol Integer ID: 90819

Keywords: protein

Abstract

GST fusion protein production

Troubleshooting

- 1 Full-length recombinant human WT α -syn, α -syn S129A, and S129D were expressed in *Escherichia coli* BL21 (DE3) (New England Biolabs, Cat # C2530H) using the bacterial expression vector pGEX-KGmyc. Following transformation protein expression was induced with 0.05 mM IPTG (isopropyl- β -D-thiogalactopyranoside), and either incubated at 37 °C for 2 hours or at room temperature for 6 hours, with shaking.
- 2 The cells grown on Terrific Broth (Thermo Scientific Cat# BP9728-2) were harvested by centrifugation at 4500× *g* at 4 °C for 20 min, and pellets were stored at -80 °C until use.
- 3 For protein purification, protein pellets were resuspended in 30 ml Lysis Buffer containing 1X PBS, 0.5 mg/ml lysozyme, 1 mM PMSF, DNase, and EDTA-free protease cocktail inhibitor (Roche Cat# 11836170001) for 15 min on ice, briefly sonicated (3 sets with 33 strikes and 30-second breaks on ice between sets), and removed the insoluble material by centrifugation at 15,000 *g* at 4 °C for 30 min.
- 4 The clarified lysate was incubated with 500 ml of glutathione-Sepharose 4B (Sigma Cat# 17-0756-01), preequilibrated with 1X PBS containing 0.1% Tween 20 and 5% glycerol (binding buffer), on a tumbler at 4 °C overnight.
- 5 The GST-bound proteins were washed four times with 30 ml binding buffer and maintained at 4 °C for pull-down assays, *in vitro* phosphorylation, and *in vitro* dephosphorylation experiments.
- 6 GST-bound proteins were occasionally eluted by TEV cleavage. In brief, 15 μ g of the fusion protein was mixed with 5 μ l of TEV protease reaction buffer (10X) and 1 μ l of TEV protease (New England Biolabs, Cat#P8112), followed by overnight incubation at 4 °C.