

Jul 15, 2024

High Molecular Weight (HMW) DNA Extraction Protocol for Tissue Sample

DOI

<https://dx.doi.org/10.17504/protocols.io.kqdg3xdrpg25/v1>

Aswini Leela Loganathan¹

¹Monash University Malaysia



Aswini Leela Loganathan

Monash University Malaysia

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kqdg3xdrpg25/v1>

Protocol Citation: Aswini Leela Loganathan 2024. High Molecular Weight (HMW) DNA Extraction Protocol for Tissue Sample. protocols.io <https://dx.doi.org/10.17504/protocols.io.kqdg3xdrpg25/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: January 22, 2024



Last Modified: July 15, 2024

Protocol Integer ID: 93868

Keywords: dna extraction protocol, dna extraction protocol for tissue sample, high molecular dna extraction for tissue sample, high molecular dna extraction, dna extraction, gdna for pacbio sequel iie library preparation, obtaining gdna, organic extraction protocol, gdna, high molecular weight, tissue sample, dna, extraction, pacbio sequel iie library preparation

Abstract

This is an organic extraction protocol used for high molecular DNA extraction for tissue samples. This protocol is suitable for obtaining gDNA for PacBio Sequel Iie library preparation and sequencing.



HMW DNA Extraction Method

9h 12m 10s

- 1 **Lysis the tissue sample.**
Add  500 μL of lysis buffer to 50 - 150 mg of Sample of sample. Make sure the tissue sample has been finely cut.
- 2 **Denatures and digest proteins that are subsequently hydrolyzed with Proteinase K.** 10s
Add  10 μL of Proteinase K and vortex for  00:00:10 .
- 3 Incubate on a shaking incubator at  55 $^{\circ}\text{C}$ for  02:00:00 (or till dissolved). 2h
- 4 **Remove RNA with RNase .**
Add  5 μL of RNase and vortex briefly.
- 5 Incubate in a shaking incubator for at  37 $^{\circ}\text{C}$.
- 6 **Partitioning of lipids and debris into an organic phase using P:C:IA.**
Add  500 μL of P:C:IA (25:24:1). Vortex until an emulsion is formed.
- 7 Centrifuge at  Room temperature for  00:10:00 at  10000 x g . 10m
- 8 Pipette  200 μL of the aqueous layer into a new tube.
- 9 **Neutralize the charges on the sugar-phosphate backbone of the DNA with Sodium Acetate.**
Add 1/10th :  20 μL 3M Sodium Acetate. Vortex gently (avoid creating bubbles).
- 10 **Precipitation Step.**
Add 2 Vol :  440 μL of ice-cold 100% Ethanol. Mix gently and slowly by inverting the tube.



- 11 Incubate at -20 °C for 02:00:00 . 2h
- 12 Centrifuge at 10000 x g, Room temperature, 00:02:00 . 2m
- 13 **Wash Step.**
Wash with ice-cold 70% ethanol.
- 14 Centrifuge for 00:02:00 at 10000 x g . 2m
- 15 Repeat **Wash Step** (Step 13 and 14) **2** more times.
- 16 Air dry for at least 00:30:00 . 30m
- 17 **Final elution.**
Add 35 µL of elution buffer (EB) and mix gently using a wide bore tips or by tapping.
- 18 **Storage.**
Store the gDNA at 4 °C